Interface Kit
RS232 – USB
Instruction Manual

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DOMINIQUE DUTSCHER SAS

INTRODUCTION

The OHAUS RS232-USB Interface Kit is a unique solution that allows users to connect a balance to a computer using a Universal Serial Bus (USB). This Interface kit needs to download the software drivers from OHAUS website to create the required virtual port on the computer.

The RS232-USB Interface Kit can be used with OHAUS products that have the RS232 interface.

SYSTEM REQUIREMENTS

- PC running Windows 98®, Windows 98SE®, Windows ME®, Windows 2000®, Windows XP® or Windows 7®
- Be forward compatible with iMac 10.7 (contact OHAUS)
- Available USB port (Type A, 4-pin, female)

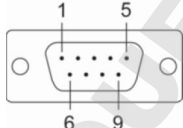
PRODUCT SPECIFICATION

◆ USB Interface

- Compliance: USB1.1 compliant, USB 2.0 compatible
- Connector: USB type A

◆ RS232 Interface

This Kit uses male DB9 connectors. Pin assignments are shown in the following diagram:

DB9 (Male)	Pin	RS232
	1	N/A
	2	RxD (in)
	3	TxD (out)
	4	N/A
	5	GND
	6	N/A
	7	N/A
	8	N/A
	9	N/A

◆ Technical Data

- Indoor use only
- Operating Temperature: 0 to 40°C
- Storage Temperature: -10 to 55°C
- Ambient Relative Humidity: maximum relative humidity 80 % for temperatures up to 30°C, decreasing linearly to 50% relative humidity at 40°C, noncondensing
- Altitude: Up to 2000 m

INTERFACE INSTALLATION

Plug the USB connector into any upstream USB port, such as a USB port on your computer. Windows will automatically detect the kit and begin installing the driver. When Windows finished installing the driver for the interface kit, it will detect a new COM port. Another RS232 connector of the kit can be plugged into the female RS232 connector on the balance.

1.

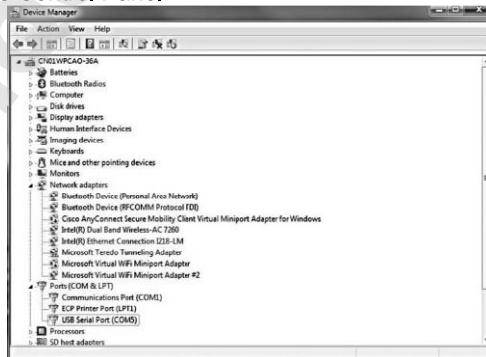
Example of Windows XP Hardware Wizard

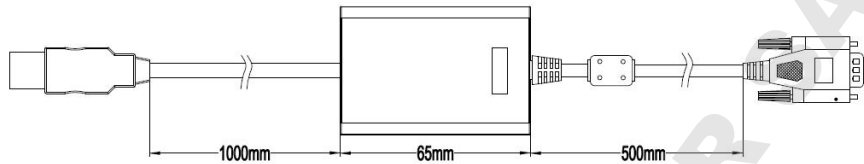
2. After clicking Finish, the virtual port should be ready for use.

Windows® typically adds the virtual port in sequence after the highest number COM port. For example, on PC's equipped with up to 4 COM ports, the virtual port will be COM5.

When using the USB interface with programs that limit the number of COM port designations, it may be necessary to assign one of these port numbers to the new virtual port. This can be done in the Port Settings of the Device Manager utility, found in the Windows Control Panel.

DIMENSIONS

COMPLIANCE

Compliance to the following standards is indicated by the corresponding mark on the product.

Marking	Standard
	EN 55032 (Class B), EN 55024
	EN 55032 (Class B)
	FCC Part 15 (Class B)

FCC Note:

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions: (1) this device may not cause harmful interference, and (2) this device must accept any interference received, including interference that may cause undesired operation.

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.

- Consult the dealer or an experienced radio/TV technician for help.
Please note that changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

Industry Canada Note

This product complies with Canadian ICES-003.

Disposal

In conformance with the European Directive 2002/96/EC on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.

Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.

Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.

For disposal instructions in Europe, refer to www.ohaus.com/weee.
Thank you for your contribution to environmental protection.

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P/N 30304103 B © 2016 Ohaus Corporation, all rights reserved
OHAUS Corporation