

# Instructions for Water Samples



1. Remove screen from faucet.
2. Run Water for 3 to 6 minutes.
3. Open container by pushing up on cap where indicated (Fig. 1).  
**NOTE:** Do not use if arrow was broken before opening.  
**NOTE:** Do not take pill out of container.
4. Fill container to the E.P.A. 100 mL fill line.
5. Close the container by pressing cap from the hinge location and pushing forward, it should securely snap shut (Fig. 2).
6. Close tab (Fig. 3) then pull the tie through round hole on post and lock vial (Fig. 4).
7. Complete information on sample label.



Figure 1



Figure 2



Figure 3



Figure 4

Note: The color of the tablet can range from white to dark gray or purple and it can be speckled in appearance. The color of the tablet may be seen to change while it remains in the sealed container when the container is exposed to light. Some tablets may fracture within the sealed container prior to use. Tablet color changes and the potential for tablet fracturing are known and expected conditions resulting from the sterilization processes used to manufacture this product. Tablets of different color or tablets that have fractured within the sealed container continue to provide the 10 mg of Sodium Thiosulfate needed to neutralize free chlorine in water samples collected.

## Corning Incorporated Life Sciences

836 North St.  
Building 300, Suite 3401  
Tewksbury, MA 01876  
t 800.492.1110  
t 978.442.2200  
f 978.442.2476

[www.corning.com/lifesciences](http://www.corning.com/lifesciences)

## Worldwide Support Offices

**ASIA / PACIFIC**  
Australia/New Zealand  
t 0402-794-347

**China**  
t 86 21 2215 2888  
f 86 21 6215 2988

**India**  
t 91 124 4604000  
f 91 124 4604099

**Japan**  
t 81 3-3586 1996  
f 81 3-3586 1291

**Korea**  
t 82 2-796-9500  
f 82 2-796-9300

**Singapore**  
t 65 6733-6511  
f 65 6861-2913

**Taiwan**  
t 886 2-2716-0338  
f 886 2-2516-7500

## EUROPE

**France**  
t 0800 916 882  
f 0800 918 636

**Germany**  
t 0800 101 1153  
f 0800 101 2427

**The Netherlands**  
t 31 20 655 79 28  
f 31 20 659 76 73

**United Kingdom**  
t 0800 376 8660  
f 0800 379 1117

## All Other European Countries

t 31 (0) 20 659 60 51  
f 31 (0) 20 659 76 73

## LATIN AMERICA

**Brasil**  
t (55-11) 3089-7419  
f (55-11) 3167-0700

**Mexico**  
t (52-81) 8158-8400  
f (52-81) 8313-8589

The Corning Family of Brands

CORNING costar cellgro PYREX AXYGEN GOSSELIN™