

### Nalgene Color-Coded "Through the Closure Style" Wash Bottles

Catalog No.	Color	Capacity (ml / oz)	Closure Size, mm	No. per Pkg	No. per Case
2422-2500	Yellow	500 / 16	28	6	24
2422-3500	Blue	500 / 16	28	6	24
2422-4500	Green	500 / 16	28	6	24
2422-5500	Red	500 / 16	28	6	24

### Nalgene Color-Coded "Unitary Style" Wash Bottles

Catalog No.	Color	Nom. Capacity (ml / oz)	Fill Line Capacity, ml	Closure Size, mm	No. per Pkg	No. per Case
2423-0500	Red, Yellow, Blue, White	500 / 16	480	38-415	4	16

### Cleaning Guidelines:

Clean Nalgene Wash Bottles with a mild, non-abrasive detergent, such as Nalgene L900 Liquid Detergent (Cat. No. 900-4000). Follow with a rinse in tap water and final rinse in distilled water. To avoid scratching the plastic, do not use abrasive cleaners or scouring pads.

### Warnings:

1. Bottles are not intended for containment of volatile solvents since vapor pressure may build inside bottle and cause fluid to siphon through spout. Please refer to our Nalgene RTK Safety Wash Bottles (Cat. No. 2425-XXXX and 2436-XXXX), and 2421-XXXX Solvent Wash Bottles for self-venting systems.
2. DO NOT fill wash bottles above the "FILL LINE".
3. DO NOT autoclave bottles.
4. Inspect bottles periodically for signs of stress (cracking or crazing/whitening of the plastic). When signs of stress are detected, please discontinue use.

### Additional Nalgene Right To Know Products include:

- Vented Unitary Safety Wash Bottles (Cat. No. 2436-XXXX)
- RTK Safety Wash Bottles (Cat No. 2425-XXXX)