
Nalgene™ Desiccator Cabinets Cat. No. 5317

Nalgene Desiccator
Cabinets provide dustand moisture-free
storage of anhydrous or
hygroscopic, biological
and chemical compounds, as well as
electronic assemblies
and any components
requiring controlled
storage conditions.

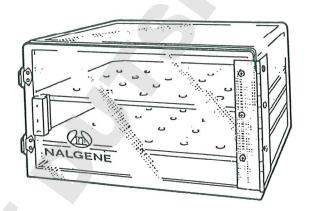

- Constructed of durable 3/8-in. thick acrylic.
- Completely transparent to permit an undistorted view of the contents.
- Cabinet door has a neoprene gasket to ensure a tight seal. Door is secured by two stainless steel latches and a full-height hinge.
- The 1/4-in. thick vented shelves are removable and adjustable.
- Desiccant tray is 1-in. deep to contain all common dry desiccants.
- Cabinet dimensions:

Cat. No. 5317 No. of shelves No. of grooves H x W x D, cm	-0070 2 3 17.8 x 30.5 x 30.5 7 x 12 x 12	-0120 3 5 30.5 x 30.5 x 30.5 12 x12 x 12	-0180 4 7 45.7 x 30.5 x 30.5 18 x 12 x12
--	--	--	--

Care and Use Instructions

CAUTION:

- Do not use concentrated sulfuric acid as a desiccant.
- Do not expose cabinets to temperatures above 50°C.
- Acrylic is strong, but it can crack or shatter if struck against a hard surface.

Cleaning

- Clean with a mild soap or detergent solution. Do not use brushes or abrasive cleaners, which will scratch the plastic.
- Acrylic labware has good chemical resistance to dilute or weak acids, aldehydes and aliphatic hydrocarbons. Do not use with acid cleaning solutions or organic solvents, including alcohols.
- Sterilize with ethylene oxide gas or gamma radiation. Do not autoclave.

For further information on this and other NalgeneTM Labware products, see the current Nalgene Labware Catalog or contact Marketing Communications, Nalge Company, Tel. (716) 586-8800, Telex 97-8242, Fax (716) 586-8431. ✓

Nalge Company

Nalgené Brand Products

Nalge Company A Subsidiary of Sybron Corporation Box 20365 Rochester, New York 14602-0365 U.S.A.

8-0404-32 1188

© Nalge Company, 1988


Printed in U.S.A.

Z (INTERNAL HEIGHT): 6.25 ±.06	Y (INTERNAL DEPTH): 11.25 ±.06 285.8 ±1.5 11.25 ±.06 285.8 ±1.5 11.25 ±.06 285.8 ±1.5	X (INTERNAL WIOTH): 11.25 ±.06 285.8 ±1.5 11.25 ±.06 285.8 ±1.5 11.25 ±.06 285.8 ±1.5	H (HEIGHT)	B (DEPTH WITH HANDLE):	A (WIDTH)	*	NOM. WEIGHT (16s):	NO. OF GROOVES	NO. OF SHELVES	CATALOG NO	
	11.25 ±.06	11.25 ±.06	7	Ξ	12	IJ.	_	E L	2	5317-0070	
158.8 ±1.5	285.8 ±1.5	285.8 ±1.5	177.8	330.2	304.8	MM	ΙΞ			0070	
158.8 ±1.5 11.25 ±.06 285.8 ±1.5 17.25 ±.06 438.1 ±1.5	11.25 ±.06	11.25 ±.06	12	13	12	IN.	2	10	E	5317-0120	
285.8 ±1.5	285.8 ±1.5	285.8 ±1.5	304.8	330.2	304.8	MM	22			-0120	
17.25 ±.06	11.25 ±.06	11.25 ±.06	18	13	12	IN.	<u> </u>	7	4	5317-0180	
438.1 ±1.5	285.8 ±1.5	285.8 ±1.5	457.2	330.2	304.8	MM	ĩ			0810	
 _	_									_	_

NEDPRENE	GASKET:
STAINLESS STEEL	LATCHES & HINGES
ACRYLIC	SHELVES & TRAY
ACRYLIC	BODY:
	- TA

NOTES:

- ALL DIMENSIONS ARE NOMINAL AND IN INCHES AND MILLIMETERS UNLESS OTHERWISE NOTED.
- 2. THIS IS NOT A PRODUCTION DRAWING, BUT IS FOR INFORMATIONAL PURPOSES ONLY, AND IS SUBJECT TO CHANGE WITHOUT NOTICE.

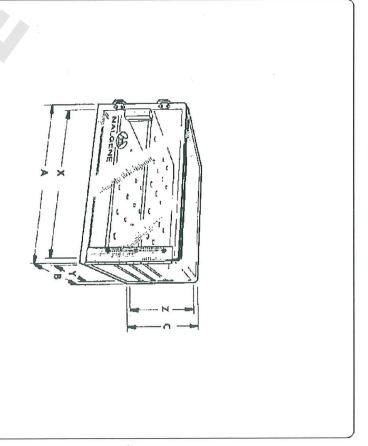

SPECS ON BACK

THIS PRODUCT SPECIFICATION
SHEET CONTAINS CONFIDENTIAL
AND/OR PROPRIETIORY INFORMATION
AND IS TO BE USED DNLY FOR THE
EXPRESSED PURPOSES AGREED UPON
BY AN AUTHORIZED DEFICER OF
THERMO FISHER SCIENTIFIC. CONFIDENTIAL

00	
0	2
_	4
ш	3
z	ā
_	77
_	65
_	7
0	0

Therno Fisher Scientific 75 Ponorono Creek Drive Rochester, NY 14625-2385, LISA

NOT TO SCALE

界 PSS -00-5317-XXXX-S 1108

_00L

DESICCATOR CABINETS