


Nalgene™ Desiccator Cabinets Cat. No. 5317

Nalgene Desiccator Cabinets provide dust- and moisture-free storage of anhydrous or hygroscopic, biological and chemical compounds, as well as electronic assemblies and any components requiring controlled storage conditions.

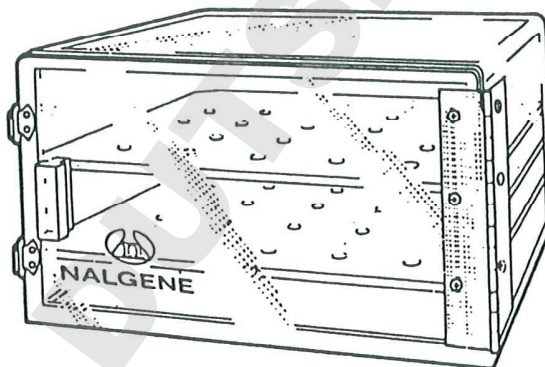

- Constructed of durable 3/8-in. thick acrylic.
- Completely transparent to permit an undistorted view of the contents.
- Cabinet door has a neoprene gasket to ensure a tight seal. Door is secured by two stainless steel latches and a full-height hinge.
- The 1/4-in. thick vented shelves are removable and adjustable.
- Desiccant tray is 1-in. deep to contain all common dry desiccants.
- Cabinet dimensions:

Cat. No. 5317	-0070	-0120	-0180
No. of shelves	2	3	4
No. of grooves	3	5	7
H x W x D, cm	17.8 x 30.5 x 30.5	30.5 x 30.5 x 30.5	45.7 x 30.5 x 30.5
in.	7 x 12 x 12	12 x 12 x 12	18 x 12 x 12

Care and Use Instructions

CAUTION

- Do not use concentrated sulfuric acid as a desiccant.
- Do not expose cabinets to temperatures above 50°C.
- Acrylic is strong, but it can crack or shatter if struck against a hard surface.

Cleaning

- Clean with a mild soap or detergent solution. Do not use brushes or abrasive cleaners, which will scratch the plastic.
- Acrylic labware has good chemical resistance to dilute or weak acids, aldehydes and aliphatic hydrocarbons. Do not use with acid cleaning solutions or organic solvents, including alcohols.
- Sterilize with ethylene oxide gas or gamma radiation. Do not autoclave.

For further information on this and other Nalgene™ Labware products, see the current Nalgene Labware Catalog or contact Marketing Communications, Nalge Company, Tel. (716) 586-8800, Telex 97-8242, Fax (716) 586-8431.

Nalge Company

Nalgene® Brand Products

Nalge Company
A Subsidiary of Sybron Corporation
Box 20365
Rochester, New York 14602-0365
U.S.A.

8-0404-32 1188


© Nalge Company, 1988

Printed in U.S.A.

CATALOG NO.	5317-0070	5317-0120	5317-0180
NO. OF SHELVES.....	2	3	4
NO. OF GROOVES.....	3	5	7
NOM. WEIGHT (LBS)....	13	22	31
	IN.	MM	IN.
A (WIDTH).....	12	304.8	12
B (DEPTH WITH HANDLE):	13	330.2	13
H (HEIGHT).....	7	177.8	12
X (INTERNAL WIDTH)....	11.25 ±.06	285.8 ±1.5	11.25 ±.06
Y (INTERNAL DEPTH)....	11.25 ±.06	285.8 ±1.5	11.25 ±.06
Z (INTERNAL HEIGHT)...	6.25 ±.06	158.8 ±1.5	11.25 ±.06

MATERIALS	
BODY.....	ACRYLIC
SHELVES & TRAY.....	ACRYLIC
LATCHES & HINGES.....	STAINLESS STEEL
GASKET.....	NEOPRENE

- NOTES:
1. ALL DIMENSIONS ARE NOMINAL AND IN INCHES AND MILLIMETERS UNLESS OTHERWISE NOTED.
 2. THIS IS NOT A PRODUCTION DRAWING, BUT IS FOR INFORMATIONAL PURPOSES ONLY, AND IS SUBJECT TO CHANGE WITHOUT NOTICE.

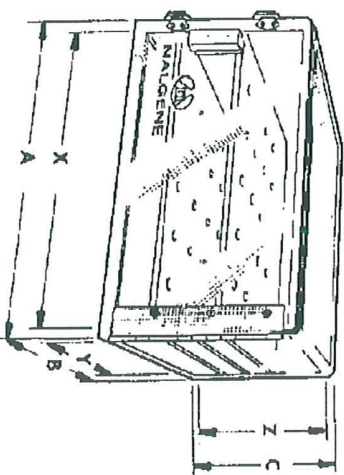

SPECS ON BACK

NOT TO SCALE

CONFIDENTIAL
 THIS PRODUCT SPECIFICATION SHEET CONTAINS CONFIDENTIAL AND/OR PROPRIETARY INFORMATION AND IS TO BE USED ONLY FOR THE EXPRESSED PURPOSES AGREED UPON BY AN AUTHORIZED OFFICER OF THERMO FISHER SCIENTIFIC.

DESICCATOR CABINETS

ThermoFisher
SCIENTIFIC
 Thermo Fisher Scientific
 75 Pomeroy Creek Drive
 Rochester, NY 14625-2385, USA

Part No.	J00	Rev	BR	Description	PSS-00-5317-XXXX-S	Rev	1108
----------	-----	-----	----	-------------	--------------------	-----	------