

# **NALGENE® PIPET WASHER-RINSERS**

## **(High-Density Polyethylene)**

**Cat. No. 5245**

The Nalgene Pipet Washer-Rinser can be used in a sink or on a drainboard. It operates as an automatic siphon. When the water level reaches the top of the siphon tube, the system drains, and the cycle repeats as long as water is supplied.

### **TO ASSEMBLE**

Connect upper tubulation to a water source with the 1/2-in. I.D. tubing provided. Connect 1-in. I.D. tubing to lower tubulation. Leave end of tubing free in sink or connect it directly in the drainline.

### **TO WASH AND RINSE PIPETS**

For best results, place pipets in a Nalgene Pipet Basket of the appropriate size. (See current Nalgene Labware Catalog, Cat. No. 5241.) Make sure basket handle does not interfere with incoming water.

Add detergent (powder, tablet, or liquid) to the pipet basket. Add pipets and place basket in the washer-rinser. Adjust water input to a low rate (about 2 liters per minute). A detergent tablet will dissolve slowly over several cycles. Rinsing will continue after the tablet dissolves.

To rinse pipets, use a higher water input rate—up to 12 liters per minute.

For more information, write or call Technical Service, Nalge Company. Tel. (716) 586-8800. Telex 97-8242. Fax (716) 586-8431

© Nalge Company, 1988

## **Nalge Company**

NALGENE® Brand products

Nalge Company  
A Subsidiary of Sybron Corporation  
Box 20365  
Rochester, New York 14602-0365 U.S.A.

8-0400-04 1088

Printed in U.S.A.