

GUIDANCE FOR USE

PROTECT Cryopreservation System

All samples should be kept cool until testing is performed.

SETUP FOR FREEZING	
<p>STEP 1 Make a suspension of the organism in the vial liquid.</p> 	<p>STEP 2 Shake vial to coat the beads. Stand for 30 seconds.</p> 
<p>STEP 3 Remove vial liquid using pastette (TS/126-H).</p> 	<p>STEP 4 Label, place in freezer. Done!</p> 
RECOVERY FROM FREEZING	
<p>STEP 5 Remove one bead using a sterile needle (TS/3-C).</p> 	<p>STEP 6 Roll bead directly onto medium or place in broth and incubate.</p> 

Have you considered? Fine Tip Pastettes - TS/126-H