

## Safety at high speed

Improve the safety of your membrane filtration workflow

You don't have to be exposed to obvious risks every time you execute your microbiological analyses. Discover a safe and fast way of performing your membrane filtration with single-use EZ-Fit® Filtration Units.

### Filtration set-up with reusable funnels

Flaming for disinfection can lead to burns or blistering.

Be careful when installing the funnel on the manifold, you might clip your fingers.




Flaming manifold head

5 sec



Flaming funnel base and inside

4 sec

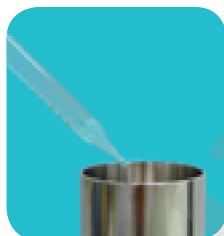

Cooling time with sterile water

3 to 5 sec



Placing filter on manifold head

5 sec



Installing filter funnel on manifold

3 sec




### Filtration set-up with sterile single-use EZ-Fit® Filtration Units



EZ-Fit® Filtration Unit set-up

1 sec

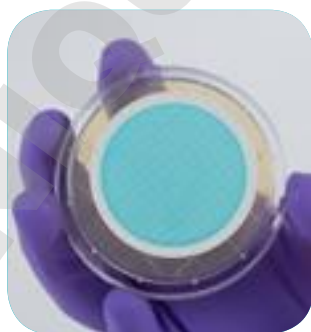

EZ-Fit® Filtration Units save you precious time while ensuring accurate results. The sterilized, disposable filtration devices for bioburden testing streamline your workflow.

Learn more:


[MerckMillipore.com/EZ-Fit](https://www.MerckMillipore.com/EZ-Fit)



For optimal microbial recovery, use EZ-Fit® Filtration Units with our Readyplate™ 55 plates. These 55 mm agar plates are small but mighty: the high filling volume ensures reproducible membrane transfer without bubbles and folds.

Learn more:

[MerckMillipore.com/ReadyPlate55](https://www.MerckMillipore.com/ReadyPlate55)

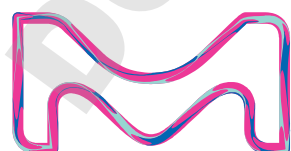

## EZ-Fit® Filtration Unit, blue base with pad

| Product Description  | Pack Size | Cat. No.  |
|--|-----------|-----------|
| EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 100 mL funnel, bulk packaging       | 48        | EFHAW100B |
| EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 100 mL funnel, individual packaging | 48        | EFHAW100I |
| EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 250 mL funnel, bulk packaging       | 48        | EFHAW250B |
| EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 250 mL funnel, individual packaging | 48        | EFHAW250I |
| EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 100 mL funnel, bulk packaging       | 48        | EFHAB100B |
| EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 100 mL funnel, individual packaging | 48        | EFHAB100I |
| EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 100mL funnel, bulk packaging        | 48        | EFHAB250B |
| EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 250 mL funnel, individual packaging | 48        | EFHAB250I |



## EZ-Fit® Filtration Unit, pink base

| Product Description  | Pack Size | Cat. No.  |
|--|-----------|-----------|
| EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 100 mL funnel, individual packaging, no pad | 48        | EFHVW10IS |
| EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 100 mL funnel, multi packaging, no pad      | 48        | EFHAW10MS |
| EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 250 mL funnel, bulk packaging, no pad       | 48        | EFHAW25BS |
| EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 100 mL funnel, multi packaging, no pad      | 48        | EFHAB10MS |
| EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 250 mL funnel, bulk packaging, no pad       | 48        | EFHAB25BS |
| EZ-Fit® Filtration Unit, white membrane, 0.22µm pore size, 100 mL funnel, multi packaging, no pad      | 48        | EFGSW10MS |
| EZ-Fit® Filtration Unit, white membrane, 0.22µm pore size, 100 mL funnel, individual packaging, no pad | 48        | EFGSW10IS |
| EZ-Fit® Filtration Unit, white membrane, 0.8µm pore size, 100 mL funnel, bulk packaging, no pad        | 48        | EFAAW10BS |
| EZ-Fit® Filtration Unit, white membrane, 0.8µm pore size, 250 mL funnel, bulk packaging, no pad        | 48        | EFAAW25BS |
| EZ-Fit® Filtration Unit, black membrane, 0.8µm pore size, 100 mL funnel, bulk packaging, no pad        | 48        | EFAAB10BS |
| EZ-Fit® Filtration Unit, black membrane, 0.8µm pore size, 250 mL funnel, bulk packaging, no pad        | 48        | EFAAB25BS |



## EZ Family

| Product Description   | Pack Size | Cat. No.   |
|---|-----------|------------|
| EZ-Stream® Vacuum Pump for liquid transfer                          | 1         | EZSTREAM1  |
| 5 meters autoclavable silicone tubing with 9.5 mm internal diameter | 1         | STREAMTUB  |
| EZ-Fit® Manifold, 1-place for EZ-Fit® Filtration Units              | 1         | EZFITEFUN1 |
| EZ-Fit® Manifold, 3-place for EZ-Fit® Filtration Units              | 1         | EZFITEFUN3 |
| EZ-Fit® Manifold, 6-place for EZ-Fit® Filtration Units              | 1         | EZFITEFUN6 |
| EZ-Fit® filtration head for EZ-Fit® Filtration Units                | 1         | EZFITMVHE1 |
| EZ-Fit® filtration head for EZ-Fit® Filtration Units                | 3         | EZFITMVHE3 |
| EZ-Fit® filtration head gasket                                      | 1         | EZFITMVG3  |



## ReadyPlate™ 55 agar plates

| Product Description                              | Pack Size | Cat. No.   |
|--|-----------|------------|
| ReadyPlate™ 55 CCA ISO 9308                      | 20x55mm   | 1467570020 |
| ReadyPlate™ 55 CCA ISO 9308                      | 20x55mm   | 1467570200 |
| ReadyPlate™ 55 Lactose TTC Agar                  | 20x55mm   | 1467590020 |
| ReadyPlate™ 55 TGE Agar                          | 20x55mm   | 1467610020 |
| ReadyPlate™ 55 Plate Count Agar                  | 20x55mm   | 1467630020 |
| ReadyPlate™ 55 Slanetz and Bartley Agar ISO 7899 | 20x55mm   | 1467650020 |
| ReadyPlate™ 55 CN Agar ISO 16266                 | 20x55mm   | 1467670020 |
| ReadyPlate™ 55 m-Green Agar                      | 20x55mm   | 1467690020 |



To place an order or receive technical assistance in Europe, please call Customer Service:  
 France: 0825 045 645 Spain: 901 516 645 Option 1  
 Germany: 069 86798021 Switzerland: 0848 645 645  
 Italy: 848 845 645 United Kingdom: 0870 900 4645

For other countries across Europe, please call: +44 (0) 115 943 0840  
 Or visit: [MerckMillipore.com/offices](http://MerckMillipore.com/offices)  
 For Technical Service visit: [MerckMillipore.com/techservice](http://MerckMillipore.com/techservice)

© 2018 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. Merck, EZ-Fit, Millipore and the Vibrant M are trademarks of Merck KGaA, Darmstadt, Germany. All other trademarks are the property of their respective owners.

Lit. No. MK\_FL2453EN 2018 - 13706 07/2018

