

## Safety at high speed

Improve the safety of your membrane filtration workflow

You don't have to be exposed to obvious risks every time you execute your microbiological analyses. Discover a safe and fast way of performing your membrane filtration with single-use EZ-Fit® Filtration Units.

### Filtration set-up with reusable funnels

Flaming for disinfection can lead to burns or blistering.

Be careful when installing the funnel on the manifold, you might clip your fingers.

					
Flaming manifold head	Flaming funnel base and inside	Cooling time with sterile water	Placing filter on manifold head	Installing filter funnel on manifold	
5 sec	4 sec	3 to 5 sec	5 sec	3 sec	

### Filtration set-up with sterile single-use EZ-Fit® Filtration Units



EZ-Fit® Filtration Unit set-up

1 sec



EZ-Fit® Filtration Units save you precious time while ensuring accurate results. The sterilized, disposable filtration devices for bioburden testing streamline your workflow.

Learn more:

[MerckMillipore.com/EZ-Fit](https://www.MerckMillipore.com/EZ-Fit)



For optimal microbial recovery, use EZ-Fit® Filtration Units with our Readyplate™ 55 plates. These 55 mm agar plates are small but mighty: the high filling volume ensures reproducible membrane transfer without bubbles and folds.

Learn more:

[MerckMillipore.com/ReadyPlate55](https://www.MerckMillipore.com/ReadyPlate55)



## EZ-Fit® Filtration Unit, blue base with pad

Product Description	Pack Size	Cat. No.
EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 100 mL funnel, bulk packaging	48	EFHAW100B
EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 100 mL funnel, individual packaging	48	EFHAW100I
EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 250 mL funnel, bulk packaging	48	EFHAW250B
EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 250 mL funnel, individual packaging	48	EFHAW250I
EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 100 mL funnel, bulk packaging	48	EFHAB100B
EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 100 mL funnel, individual packaging	48	EFHAB100I
EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 100mL funnel, bulk packaging	48	EFHAB250B
EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 250 mL funnel, individual packaging	48	EFHAB250I



## EZ-Fit® Filtration Unit, pink base

Product Description	Pack Size	Cat. No.
EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 100 mL funnel, individual packaging, no pad	48	EFHVW10IS
EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 100 mL funnel, multi packaging, no pad	48	EFHAW10MS
EZ-Fit® Filtration Unit, white membrane, 0.45µm pore size, 250 mL funnel, bulk packaging, no pad	48	EFHAW25BS
EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 100 mL funnel, multi packaging, no pad	48	EFHAB10MS
EZ-Fit® Filtration Unit, black membrane, 0.45µm pore size, 250 mL funnel, bulk packaging, no pad	48	EFHAB25BS
EZ-Fit® Filtration Unit, white membrane, 0.22µm pore size, 100 mL funnel, multi packaging, no pad	48	EFGSW10MS
EZ-Fit® Filtration Unit, white membrane, 0.22µm pore size, 100 mL funnel, individual packaging, no pad	48	EFGSW10IS
EZ-Fit® Filtration Unit, white membrane, 0.8µm pore size, 100 mL funnel, bulk packaging, no pad	48	EFAAW10BS
EZ-Fit® Filtration Unit, white membrane, 0.8µm pore size, 250 mL funnel, bulk packaging, no pad	48	EFAAW25BS
EZ-Fit® Filtration Unit, black membrane, 0.8µm pore size, 100 mL funnel, bulk packaging, no pad	48	EFAAB10BS
EZ-Fit® Filtration Unit, black membrane, 0.8µm pore size, 250 mL funnel, bulk packaging, no pad	48	EFAAB25BS



## EZ Family

Product Description	Pack Size	Cat. No.
EZ-Stream® Vacuum Pump for liquid transfer	1	EZSTREAM1
5 meters autoclavable silicone tubing with 9.5 mm internal diameter	1	STREAMTUB
EZ-Fit® Manifold, 1-place for EZ-Fit® Filtration Units	1	EZFITEFUN1
EZ-Fit® Manifold, 3-place for EZ-Fit® Filtration Units	1	EZFITEFUN3
EZ-Fit® Manifold, 6-place for EZ-Fit® Filtration Units	1	EZFITEFUN6
EZ-Fit® filtration head for EZ-Fit® Filtration Units	1	EZFITMVHE1
EZ-Fit® filtration head for EZ-Fit® Filtration Units	3	EZFITMVHE3
EZ-Fit® filtration head gasket	1	EZFITMVG3



## ReadyPlate™ 55 agar plates

Product Description	Pack Size	Cat. No.
ReadyPlate™ 55 CCA ISO 9308	20x55mm	1467570020
ReadyPlate™ 55 CCA ISO 9308	20x55mm	1467570200
ReadyPlate™ 55 Lactose TTC Agar	20x55mm	1467590020
ReadyPlate™ 55 TGE Agar	20x55mm	1467610020
ReadyPlate™ 55 Plate Count Agar	20x55mm	1467630020
ReadyPlate™ 55 Slanetz and Bartley Agar ISO 7899	20x55mm	1467650020
ReadyPlate™ 55 CN Agar ISO 16266	20x55mm	1467670020
ReadyPlate™ 55 m-Green Agar	20x55mm	1467690020



To place an order or receive technical assistance in Europe, please call Customer Service:  
 France: 0825 045 645 Spain: 901 516 645 Option 1  
 Germany: 069 86798021 Switzerland: 0848 645 645  
 Italy: 848 845 645 United Kingdom: 0870 900 4645

For other countries across Europe, please call: +44 (0) 115 943 0840  
 Or visit: [MerckMillipore.com/offices](http://MerckMillipore.com/offices)  
 For Technical Service visit: [MerckMillipore.com/techservice](http://MerckMillipore.com/techservice)

© 2018 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. Merck, EZ-Fit, Millipore and the Vibrant M are trademarks of Merck KGaA, Darmstadt, Germany. All other trademarks are the property of their respective owners.

Lit. No. MK\_FL2453EN 2018 - 13706 07/2018

