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Manual instruction : Up/Down timer 437/2E

3 countdowns 100h
3 memories / 3 rings of 1 min at the end of time
Clock – 1/100th sec Stopwatch on 100h

A- Programming 3 timers T1-T2 and T3 (3 different sounds)

Select «Clock/Timer»
a. Programming T1 : Press on for 4 seconds T1 (the digits flash) → program hours / minutes and seconds by pressing the HR, MIN and SEC buttons
→ memory by pressing T1 → Run or pause by pressing «Start/Stop» or T1
Ring 1 minute at the end of cycle (timer recount in ascending order) → Stop by pressing «Start/Stop» button or T1.
Reset by «Clear» button, redisplay time programmed by pressing «Start/Stop» button
b-Programming T2 and T3, same procedure as T1

B- Using 3 countdowns at the same time (T1-T2-T3)

For programming follow the instructions above
Start by «Start/Stop» button → Run or pause time during the count in ascending order by pressing T1, T2 and T3
Reset timers by pressing «Clear» button
Redisplay memorized time by pressing «Start/Stop» button

C- Programming clock

Call display by pressing «Clock/Timer»
Program by pressing for 4 seconds on «Clock/Timer» (the digits flash), program hours / minutes and seconds by pressing the HR, MIN and SEC buttons
Press again on «Clock/Timer» for storage

D- Mode Stopwatch

Same display as the clock : Run or pause by pressing «Start/Stop» button, reset by pressing «Clear» button

E- Replaceable battery

2 x LR44 batteries, observe the correct polarity.

Note : Be careful to remove the material and used batteries in an environmentally suitable location.

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