WUC-D

Ultraschallreiniger Bedienungsanleitung Ultrasonic Cleaner Operating Manual

gilt für / written for
WUC- D03H/ D06H/ D10H/ D22H

witeg Labortechnik GmbH

(English manual at page 9)

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Contents

1 Introduction	10
2 Safety Instructions	10
2.1 Description of symbols.....	10
2.2 Instructions	10
3 Product description	12
3.1 Key features	12
3.2 Operating panel	12
4 Packing content	13
5 Positioning the device and commissioning	13
6 Job description	13
6.1 Menu structure diagram	13
6.2 Main menu	13
6.3 Menu for timer setting.....	14
6.4 Menu for temperature setting	14
6.5 Ultrasonic setting	14
7 Maintenance and cleaning	15
8 Technical data	15
9 Certificate	16

DOMINIQUE DUTSCHER SAS

1 Introduction

Thank you for purchasing our **WUC-D Ultrasonic Cleaner**. You did a good choice buying a **Wisdom** product.

Before use, please read this manual and use it properly.

To cover the user's health and avoid the product's destruction you have to follow the safety instructions of the second chapter.

Please keep this operating manual in a save place for future reference and hand it out if you consign the device to a third person.

Witeg reserves to modify the capsule or make amendments for the improvement of quality and efficiency without prior notice.

2 Safety Instructions

2.1 Description of symbols

WARNING

Information for avoiding injuries or fatal accidents.

ATTENTION

Safety instruction to avoid a damage of the product.

ATTENTION

Information about technical parameters.

Environment

Information about the operating environment.

WARNING of explosions

Safety instruction to avoid explosions.

ATTENTION to high temperatures

Instructions to avoid injuries caused by high temperatures.

2.2 Instructions

For a safety and faultless using of the device it is indispensable that you read the following instructions purposely and follow them by the use.

- Connect the device only to the suitable power supply. Please have a look to the identification plate.

- Use the device only in electric circuits with ground fault circuit interrupter.
- Check the power supply cord before you stick it into the electrical outlet.
- Do not use power bars to extend the power supply cord.
- Only use extension cables with sufficient cable cross section.
- Do not touch the device with wet hands during the usage because this might cause an electrical shock.
- If unexpected sound, smell or smog is generated by the device, pull out the main plug and contact the manufacturer or your supplier immediately.
- Disconnect the power supply plug before cleaning or maintenance.
- If you use radiation or contaminated samples the responsibility is only up to you.
- Switch the device off during when you load the device.
- Avoid injuries: never put your hands or other parts of your body in the working device.
- Do not clean living creatures in this device.

- It is important that the device stands up on a horizontal, stable and firm subsurface to avoid vibrations.
- Make sure that the device can not slip away.
- To avoid heat congestion or fires make sure that there is enough space for air circulations.
- Avoid direct solar radiation.
- Do not use the device near highly combustible materials.
- Do not set the device outdoors.
- It is superiorly to set the device in a good illuminated place to avoid accidents engendered by handling errors.
- To set the device in wet or dusty places could cause overheating, short circuits or fires.

- Do not use this device in area with potentially explosive atmosphere.

- Repairs and amendments should only made by **witeg** authorized people. Contraventions exclude the warranty.

- Do not put heat-sensitive items at the device.
- Warn all attendees if you leave the device unattended.
- Pay attention when you lift up the lid, hot condensation water may drop down or hot steam could come out.
- Do not clean the hot device.


- Use this device only if you read the whole safety instructions.
- This device should only be used by persons who are familiarized with safety instructions made for laboratories.
- Pay attention that no liquid reaches the controller or inward of the capsule.
- You have to keep the device clean if you want to get a longer life of it.
- Wear corresponding protective work clothing during the usage.
- Pull out the power cable if you didn't want to use the device for a longer time.

3 Product description

3.1 Key features

- Digital timer and temperature adjustment.
- Frequency 40 kHz.
- Temperature up to 80 °C adjustable.
- Stainless steel bath.
- Drain valve for devices up to 10 liter capacity.
- Stainless Steel lid inclusive.
- Over temperature protection by safety switch.

3.2 Operating panel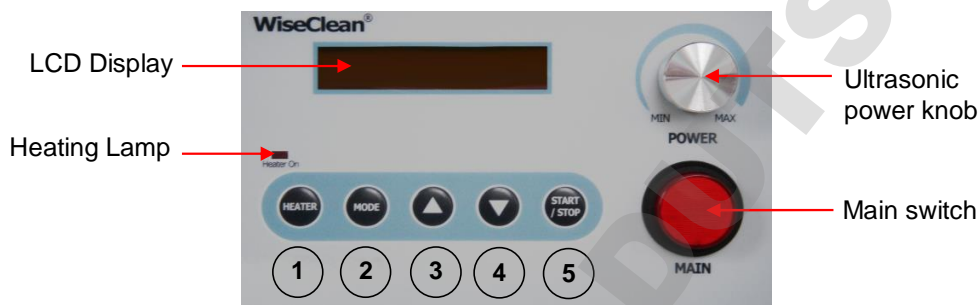

- 1 – HEATER button
- 2 – MODE button
- 3 – Arrow up button
- 4 – Arrow down button
- 5 – START/STOP button

Heating lamp

- Shines when heater is on.

Ultrasonic power knob (not matching for D03H)

- Power adjustment of ultrasonic cleaner.

Main switch

- To switch the device on or of.

HEATER button

- To switch the heater on or of.

MODE button

- To change menu item.

Arrow up button / Arrow down button

- To change timer and temperature values.

START/STOP button

- To switch the ultrasonic cleaning on or of.

4 Packing content

Article Amount	D03H	D06H	D10H	D22H
1	Ultrasonic Cleaner			
1	Stainless steel lid			
1	Power cord			
1	Operating manual			


5 Positioning the device and commissioning

1. Follow especially the safety instructions marked with following symbol:
2. Remove the package and perhaps soiling carefully.
3. If there are any unexpected transport damages, please inform your forwarder and your customer service.
4. Before you connect the device with the electrical outlet, please let some time laps until the device is acclimatised.
5. Make sure that you follow all safety instructions.
6. Fill enough cleaning liquid in. Minimum fill the half of the bath. Maximum filling level is 3 cm under bath edge.
7. Close the drain valve.
8. Connect the device with the power outlet.
9. Fill enough operating liquid in.
10. Switch the device on and start the process (look at job description).

6 Job description

6.1 Menu structure diagram

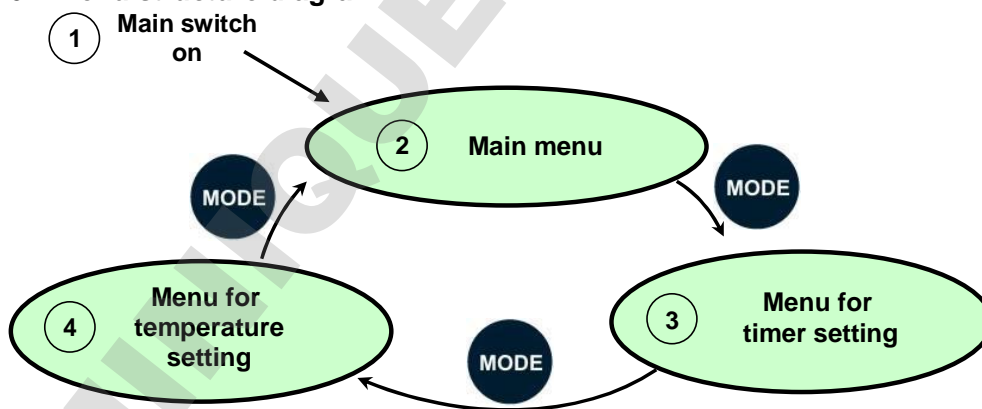6.2 Main menu

Main menu and the last used values are showed after switching the device on.
f.e.:

WAIT	10 m	13° Off
Status ultrasonic cleaner	Status Timer	Present temperature Heater is off

Oder z.B.:


WAIT	30 s	13° / 36 °
Status ultrasonic cleaner	Status Timer	Present temperature Heater is on Set value 36 °C

Switch heating on or off by HEATER button  in main menu.

Use START/STOP button  to switch ultrasonic cleaner on and off.

6.3 Menu for timer setting

Push MODE button in main menu to reach timer setting menu: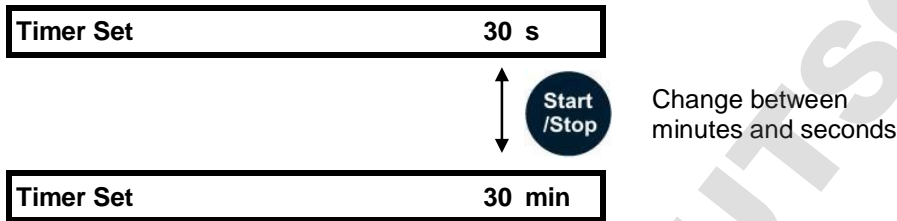


Use arrow buttons   to adjust time.

Ultrasonic cleaner will be automatically switched off when time is elapsed.

6.4 Menu for temperature setting

Push MODE button in menu for time setting (6.2) to open temperature setting menu item: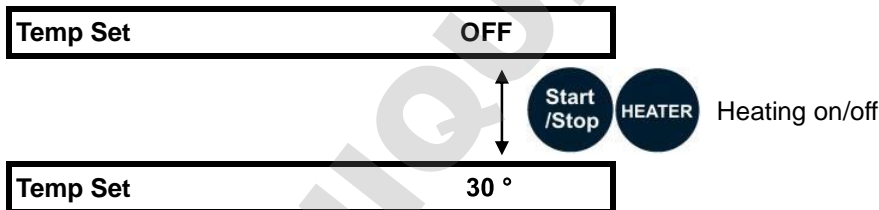


Switch heating off and use arrow buttons   to adjust temperature. Use Mode button to open main menu item.

6.5 Ultrasonic setting (not matching for D03H)

Use ultrasonic power knob to adjust the power of ultrasonic cleaning.

Use START/STOP button to switch ultrasonic on/off.

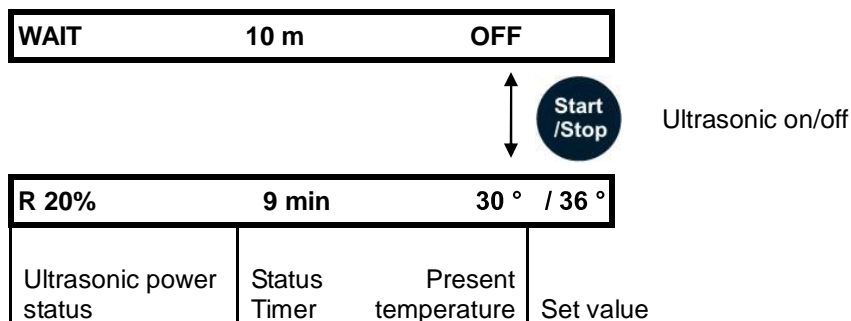7 Maintenance and cleaning

- Please read the whole safety instructions before you start to maintain or clean the device.
- Pull the power plug out of the electrical outlet.
- Only use care cleaning agents and soft cleaning rags.
- Organic lotions, strong chemicals and rough cleaning rags can damage the device.
- If you not want to use the device for a long time, pull of the power plug and keep it dry in the package.
- Only use original spare parts.
- There is no need to maintenance the device.

8 Technical data

	WUC-D03H	WUC-D06H	WUC-D10H	WUC-D22H
Capacity	3,3 l	6 l	10 l	22 l
Temperature	30-80 °C	30-80 °C	30-80 °C	30-80 °C
Capsule size	266x163x252	366x178x303	366x266x308	570x327x312
Bath size	240x140x100	290x150x150	290x240x150	500x300x150
Drain valve	no	no	yes	yes
HF-Power (output)	100 W	253 W	407 W	528 W
Heating power	172 W	172 W	258 W	344 W
Overall power	290 W	425 W	665 W	872 W
HF-Frequency	40 kHz			
Weight	3,7 kg	5,9 kg	8,1 kg	12,3 kg

9 Certificate

Certificate

for

The quality and all features were checked by the manufacturer before the shipment.

We grant from date of purchase

two years guaranty.

This certificate excludes damages by natural disasters or incorrect usages by the costumer.

Please look on your account and complete following table:

Article	WUC-D Ultrasonic Cleaner
Typ	
Serialno.	
Date	

witeg Labortechnik GmbH
Am Bildacker 16
D-97877 Wertheim
TEL: +49-9342-9301-0
FAX: +49-9342-9301-77
Email: info@witeg.de
www.witeg.de

DOMINIQUE DUTSCHER SAS

DOMINIQUE DUTSCHER SAS