
Finnpipette®

Stepper

Instructions for Use
 Bedienungsanleitung
 Guide d'utilisation
 Instrucciones de uso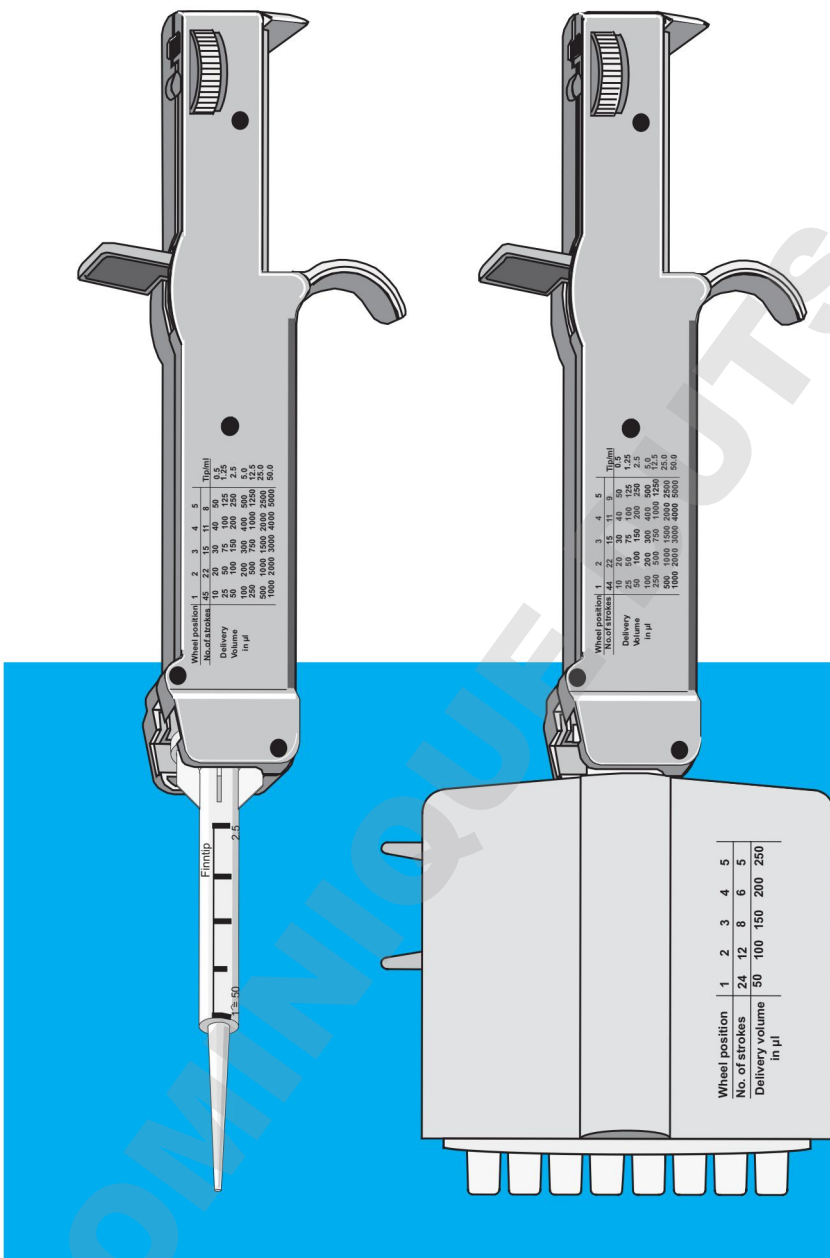


Thermo
 SCIENTIFIC


Thermo Fisher Scientific Oy
 Ratastie 2, P.O.Box 100,
 FI-01621 Vantaa, Finland
 info.pipettes@thermofisher.com
 www.thermo.com/finnpipette

Tip no. Tip Nr. Cône n° Punta n°	Pick up Volume ml Aufnahme-Volumen ml Volume prélevé ml Volumen de aspiracion ml	Delivery volume µl Dispensier-Volumen µl Volume distribué µl Volumen de dispensacion µl					
		1	0.5	10	20	30	40
2	1.25	25	50	75	100	125	
3	2.5	50	100	150	200	250	
4	5.0	100	200	300	400	500	
5	12.5	250	500	750	1000	1250	
6	25	500	1000	1500	2000	2500	
7	50	1000	2000	3000	4000	5000	
8-ch	1250	50	100	150	200	250	
Number of Strokes Anzahl Dispensierschritte Nombre de distributions N° de dispensaciones		1-ch	44	22	15	11	9
		8-ch	24	12	8	6	5
Wheel Position Volumeneinstellung Nr. Position de la molette de sélection Posicion N°		1	2	3	4	5	



TABLE 1, TABELLE 1, TABLEAU 1, TABLA 1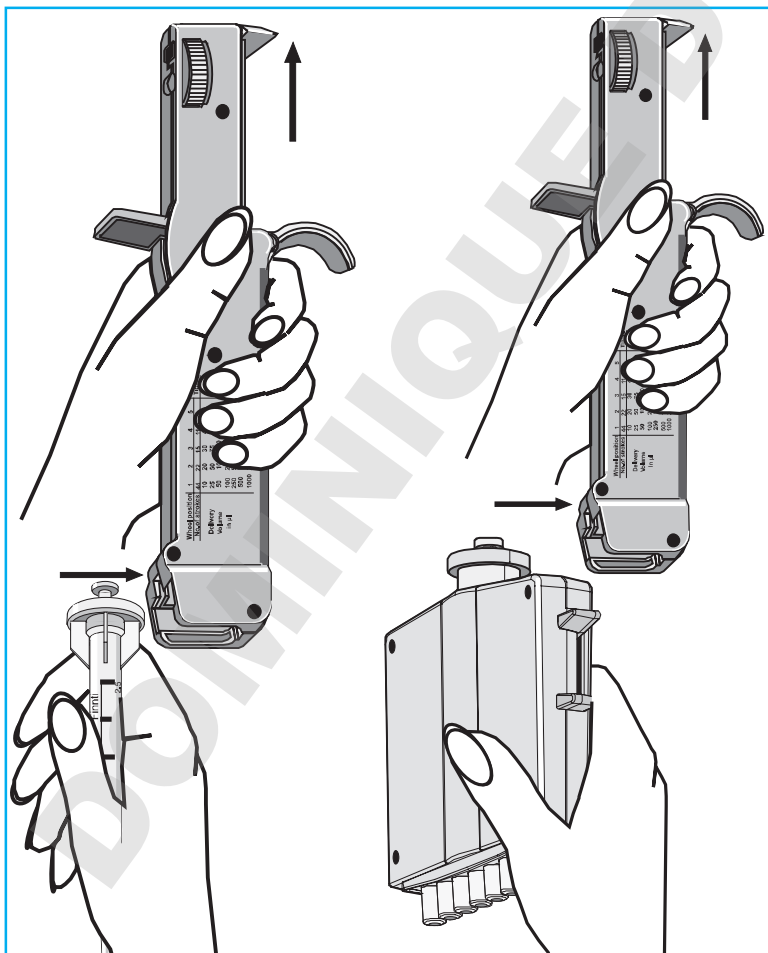

1

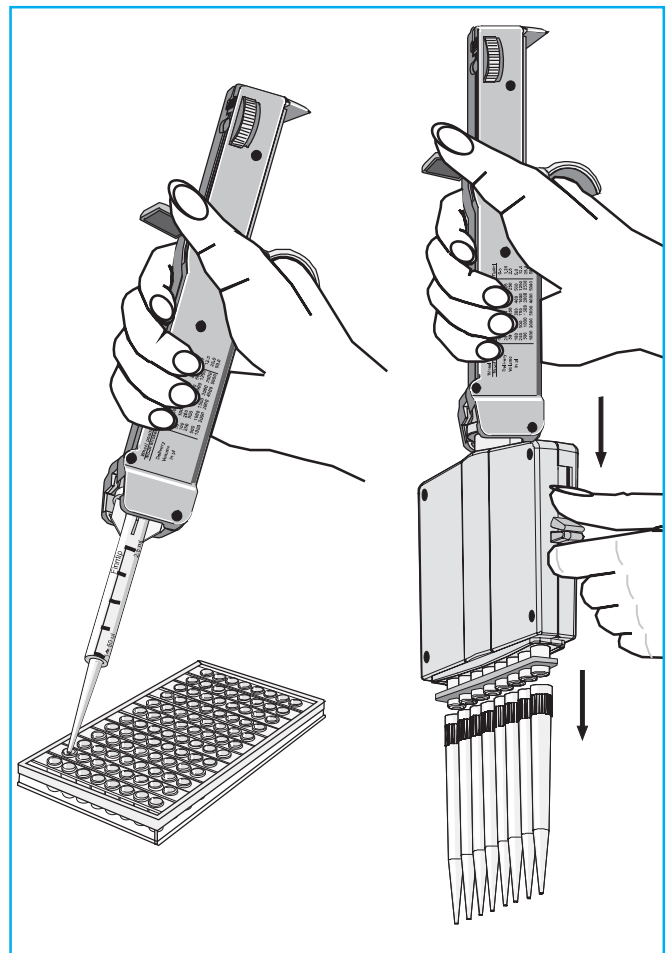2

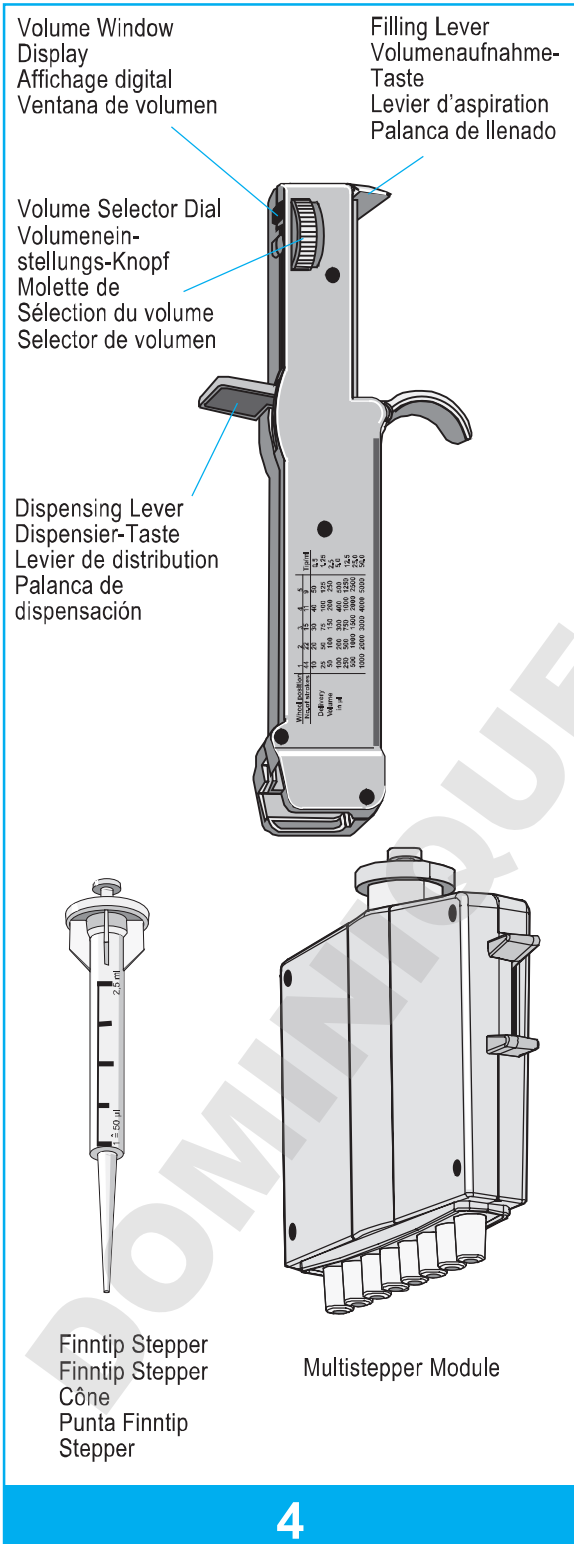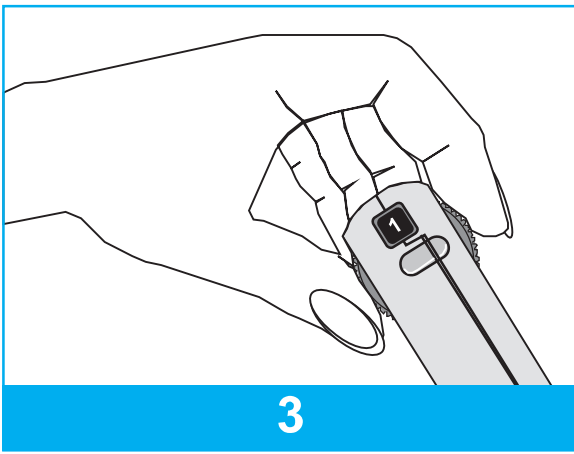CONTENTS

ENGLISH

DESCRIPTION	4
ADJUSTING THE DELIVERY VOLUME	4
MULTISTEPPEL MODULE	5
STORAGE	6
SERVICE	7
SPECIFICATIONS	7
PACKAGE	8
MULTISTEPPEL MODULE SPARE PARTS	23

INHALT

DEUTSCH

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DESCRIPTION

The Finnpiquette Stepper is an adjustable dispenser for rapid, one-handed repetitive dispensing.

This convenient repeater pipette dispenses up to 44 times in succession. Five delivery volumes are possible with each syringe size, for a volume range covering 10 μl to 5000 μl . The Finnpiquette Stepper consists of a handle and seven different tips.

The handle has three user-operated features:

- Dispensing lever
- Filling lever
- Volume selector dial.

It also has a volume selection window and an opening for the tips.

FINNTIP STEPPER

Table 1 lists the loading capacity and delivery volumes for each of the seven Finntip Stepper models. Delivery volumes are given for each position of the selector dial. The complete tip, with a piston inside a cylinder unit, is ready to attach to the handle. A special adapter must be used to attach tips 6 and 7. Insert the tips into the adapter, and turn to lock in place.

For your convenience, presterilized Finntip Steppers are also available.

ATTACHING THE TIP

1. Remove the appropriate tip from the box.
2. Make sure the filling lever is in the up position.
3. Insert the tip into the handle as shown in (Fig. 1)

Make sure that the cylinder flange and piston rod flange both slide into the appropriate grooves in the handle. Press the tip into the handle as far as it will go.

To detach the tip:

Make sure that the filling lever is in the up position. Pull the tip out of the handle.

ADJUSTING THE DELIVERY VOLUME

To obtain the desired delivery volume, select the appropriate Steptip and adjust the length of the piston stroke by turning the volume selector dial.

Example: For a delivery volume of 250 μl : **According to Table 1**, this volume can be delivered with either tip 3 (9 strokes) or tip 5 (44 strokes).

In this case, select tip 5 and insert it into the handle. Turn the volume selector dial to position 1 (250 μl) as shown in Fig. 2. Dispense with 44 strokes, as indicated in **Table 1**.

FILLING AND PRIMING THE TIP

1. The filling lever should be in the up position.

2. Dip the tip into the liquid and depress the filling lever steadily to the down position. Raise the filling lever completely and depress it again, repeating this procedure until no more air bubbles come out of the tip. Keep the tip below the liquid surface at all times.
3. The tip should now be filled. A small air bubble will remain under the plunger. This does not affect accuracy when the plunger approaches the bottom of the cylinder because dispensing stops.
4. Once the tip has been properly filled and primed, there is no need to repeat this procedure when refilling.

DISPENSING

1. For maximum hand comfort, hold the FinnpiPETTE Stepper lightly in your hand.
2. Fill the tip according to the instructions.
After filling the tip, the liquid delivered with the first stroke must be discarded.
3. Start dispensing by depressing the dispensing lever to the down position. Depress once for each dispensing stroke. Do not let the lever snap back.

CARE OF THE FINNPIPETTE STEPPER

Make sure your FinnpiPETTE Stepper is properly stored when not in use. The outside surfaces should be checked for dust and dirt every day. Do not allow liquids to enter the FinnpiPETTE Stepper's dispensing mechanism.

MULTISTEPPER MODULE

INTRODUCTION

The Multistepper module is an 8-channel cylinder unit for use with the FinnpiPETTE Stepper. This module has eight individual channels, each calibrated to simultaneously deliver the same amount of liquid. The 9 mm channel spacing is suitable for 96 well microplates.

With a single loading of the FinnpiPETTE Stepper, the Multistepper module can deliver $24 \times 50 \mu\text{l} = 192$ wells, $12 \times 100 \mu\text{l} = 96$ wells, $8 \times 150 \mu\text{l} = 64$ wells, $6 \times 200 \mu\text{l} = 48$ wells or $5 \times 250 \mu\text{l} = 40$ wells.

REAGENT

Pour the reagent into a clean reagent reservoir for easy loading.

ADJUSTING THE DELIVERY VOLUME

The volume selection table on the Multistepper module housing (Fig.1) describes how to obtain the desired delivery volume.

The volume selector dial provides a choice of five volume settings. Read vertically 1, 2, 3, 4 and 5.

Note that the maximum number of strokes for the Multistep module is not the same as for the single channel tips since the total stroke length is about 40% shorter. Therefore, the number of strokes is as follows : 24, 12, 8, 6 and 5.

Please note that the first stroke is not included in the total number of strokes. Select the volume by turning the volume selector dial (Fig. 1).

ATTACHING THE MODULE

The piston of the module should be in the lowest position before inserting the module to the handle. The filling lever should be in the upmost position.

NOTE:

If Multistep module is not in continuous use the dispensing or filling lever must be pressed harder at first, because the grease in piston may be sticky. If Multistep module is stored for long periods, the pistons may jam. When that happens and pistons are about 1 mm out of the tip cones, press the tip cones firmly against an even surface, in other case press the filling lever until piston are released.

TIPS

Attach eight tips firmly to the tip cone assembly. Insert the tips into the reagent.

Gently press the loading lever down. This fills the tips with reagent. All the tips should contain the same amount of reagent. This can be checked visually.

Withdraw the tips from the reagent, touching the edge of the reagent reservoir to remove excess liquid.

WARNING:

If the loading lever is pressed down too quickly, the liquid entering the tips may foam and enter the pipette. If this does happen, it will cause a loss of accuracy and possibly corrosion, contamination and malfunction.

DISPENSING

Dispense the first stroke back into the reagent reservoir in order to prime the tips properly. Align the tips with the first column of wells.

Press the dispensing lever down. This operation dispenses the pre-selected volume of reagent (either 50, 100, 150, 200 or 250 μ l).

Let the dispensing lever return to its original position.

Raise and gently tap the tips against the walls of the wells to remove any remaining drops from the tips.

Proceed to the next column of wells and repeat the dispensing procedure.

To avoid carry-over, tips should be changed before loading and dispensing a new reagent. Use the tip ejector (Fig. 3) to remove the tips.

QUICK UNLOADING OF THE TIPS

To quickly empty tips filled with reagent, raise the loading lever.

PARTIAL TIP FILLING

When necessary, the tips can be partially filled by depressing the loading lever part of the way down.

STORAGE

The Multistepper module should be stored with the filling lever in the upmost position. This is essential to ensure proper function of the pipette.

However, if MULTISTEPPER MODULE is stored for long periods, the pistons may jam. When that happens and pistons are about 1 mm out of the tip cones, press the tip cones firmly against an even surface, in other case press the filling lever until piston are released. (Fig. 4).

SERVICE

The Multistepper module requires a limited amount of maintenance and service. If used daily, clean the cylinders and O-ring base with a napless cloth once or twice a year. Similarly the O-rings and pistons should be lightly greased with the lubricant grease provided with the pipette.

There is no need to adjust the module.

In the case of malfunction, please return the Multistepper module to the nearest Thermo Fisher Scientific subsidiary or authorized agent.

If harmful substances have been handled with the Multistepper, please clean the module thoroughly before shipping it for service.

SPECIFICATIONS

Weight	162 g
Spacing	9 mm
Channels	8
Volume Uptake	1250 μ l/channel (maximum)
Volume Delivery	24x50 μ l/channel 12x100 μ l/channel 8 x150 μ l/channel 6x200 μ l/channel 5x250 μ l/channel

Accuracy	50 μ l: ± 10 %
	100 μ l: $\pm 6,25$ %
	150 μ l: $\pm 4,5$ %
	200 μ l: $\pm 3,25$ %
	250 μ l: $\pm 2,5$ %
Precision	50 μ l: ± 10 %
	100 μ l: $\pm 6,25$ %
	150 μ l: $\pm 4,5$ %
	200 μ l: $\pm 3,25$ %
	250 μ l: $\pm 2,5$ %

Specifications are subject to change without notice as part of Thermo Fisher Scientific continuous program of product improvement. The pipette is calibrated at the factory and cannot be recalibrated.

MATERIALS

The Finnpiquette Stepper is made of mechanically durable and chemically resistant materials.

PACKAGING

The Multistepper is shipped in a specially designed carton package. The package contains the following items:

- Finnpiquette Stepper
- Tube of grease
- Sample tips
- Instruction manual
- Multistepper module

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Product specifications are subject to change without prior notice. Finnipette® and Finntip® are registered trademarks of Thermo Scientific.
© 2007 Thermo Fisher Scientific Inc. All rights reserved.