


# 90 mm GLASS FILTER HOLDER KIT

## with stainless steel screen

### Description

The 90 mm Glass Filter Holder Kit allows the user to vacuum filter liquids for analysis of particulate or biological contamination retained on the filter surface, or to produce medium quantities of ultraclean filtrate. The user must aseptically install a sterile filter in the separately sterilized filter holder. Required accessories include: stainless smooth-tip filter forceps, a sidearm vacuum filtering flask; and a vacuum source.

### Kit Components

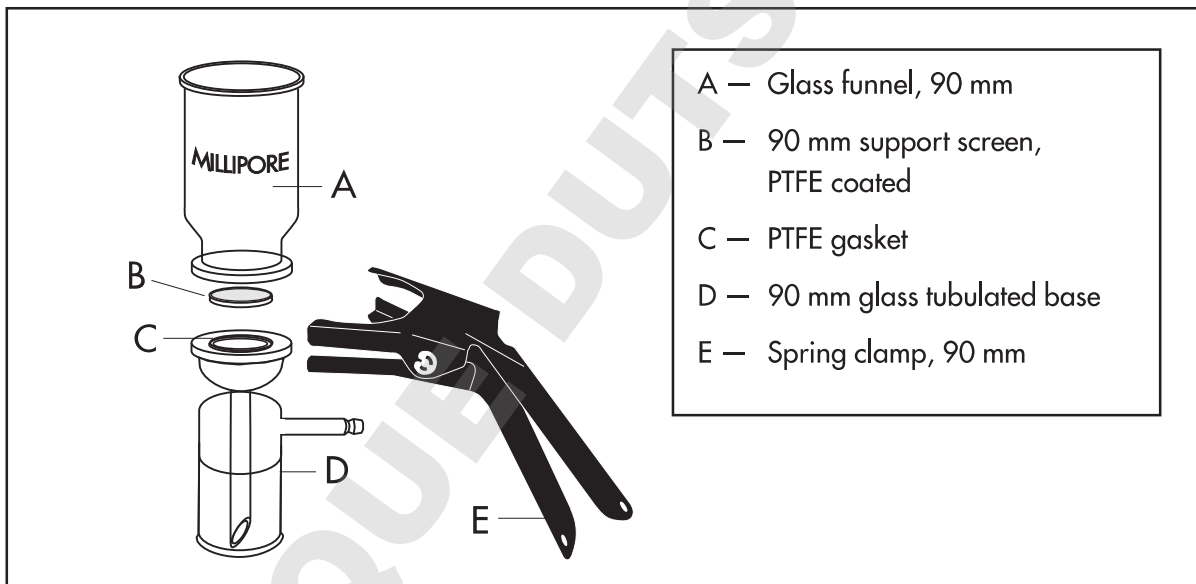

### Technical Assistance

For more information, contact the Millipore office nearest you. In the U.S., call **1-800-MILLIPORE** (1-800-645-5476). Outside the U.S., see your Millipore laboratory catalogue for the phone number of the office nearest you. You can reach us by e-mail at [tech\\_service@millipore.com](mailto:tech_service@millipore.com) or visit our web site ([www.millipore.com](http://www.millipore.com)).

# MILLIPORE

P36357, Rev. —, 02/01

© 2001 Millipore Corporation  
Printed in the U.S.A. All rights reserved.

Millipore is a registered trademark of the Millipore Corporation or an affiliated company.

## Product Ordering Information

This section lists the catalogue numbers for the 90 mm Glass Filter Holder Kit. See the Technical Assistance section for information about contacting Millipore. You can also buy Millipore products on-line at [www.millipore.com/purecommerce](http://www.millipore.com/purecommerce).

Description	Catalogue No.	Qty.	Diagram Key
90 mm Glass Filter Holder Kit consists of spring clamp, funnel, base, support screen and gaskets	XX10 090 20	1	
<b>Replacement Parts</b>			<b>Diagram Key</b>
Funnel, 90 mm, 1 ltr glass	XX10 090 00	1	A
Support screen, 90 mm, PTFE coated	XX10 090 02	1	B
Spring clamp, 90 mm, anodized	XX10 090 03	1	E
Glass base, 90 mm	XX10 090 04	1	D
Gasket, PTFE	XX10 090 10	10	C
<b>Accessories</b>			
Vacuum filtering flask, 1 liter	XX15 047 05	1	
Vacuum filtering flask, 2 liter	XX16 047 05	1	
Vacuum filtering flask, 5 liter	XX16 047 06	1	
Filter forceps	XX62 000 06	1	
Solvent filtering dispenser	XX66 025 00	1	
Vacuum Pressure Pump, 115V	XX55 000 00	1	
Vacuum Pressure Pump, 220V	XX55 220 50	1	
3/16" silicone tubing for vacuum use	XX71 000 04	1	

## Standard Warranty

**Millipore Corporation** ("Millipore") warrants its products will meet their applicable published specifications when used in accordance with their applicable instructions for a period of one year from shipment of the products. **MILLIPORE MAKES NO OTHER WARRANTY, EXPRESSED OR IMPLIED. THERE IS NO WARRANTY OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE.** The warranty provided herein and the data, specifications and descriptions of Millipore products appearing in Millipore's published catalogues and product literature may not be altered except by express written agreement signed by an officer of Millipore. Representations, oral or written, which are inconsistent with this warranty or such publications are not authorized and if given, should not be relied upon.

In the event of a breach of the foregoing warranty, Millipore's sole obligation shall be to repair or replace, at its option, the applicable product or part thereof, provided the customer notifies Millipore promptly of any such breach. If after exercising reasonable efforts, Millipore is unable to repair or replace the product or part, then Millipore shall refund to the customer all monies paid for such applicable product or part. **MILLIPORE SHALL NOT BE LIABLE FOR CONSEQUENTIAL, INCIDENTAL, SPECIAL OR ANY OTHER INDIRECT DAMAGES RESULTING FROM ECONOMIC LOSS OR PROPERTY DAMAGE SUSTAINED BY ANY CUSTOMER FROM THE USE OF ITS PRODUCTS.**

# MILLIPORE