


PureProteome™ Magnetic Stand, 8-well


Cat. No. LSKMAGS08

⚠ See Cautions on reverse side before using magnetic stand!

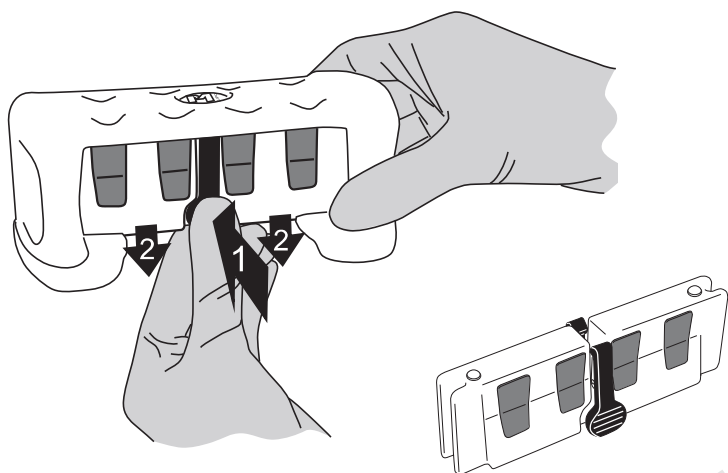


Disengage magnet

Hold magnetic stand securely and pinch both clips while pulling down on magnet bar. Place magnet bar upright on bench away from metal objects/ magnetic media.

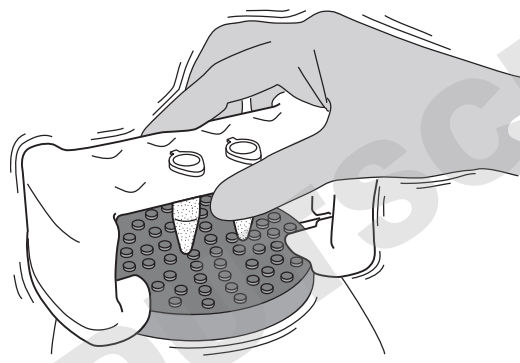

Resuspend beads

Grip magnetic stand firmly and align base of stand with 3" vortex platform. **Hold firmly while vortexing!**

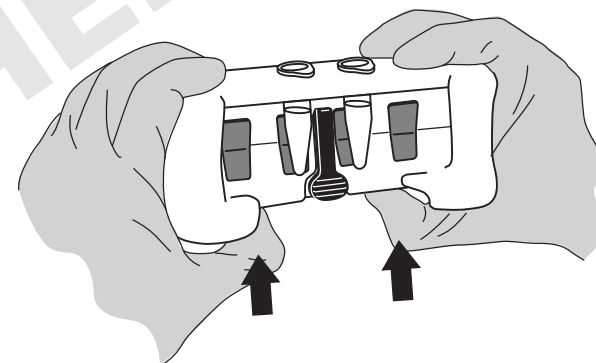


Re-engage magnet

Slide magnet bar up into stand until it clicks into place.



CAUTION – CONTAINS MAGNETS

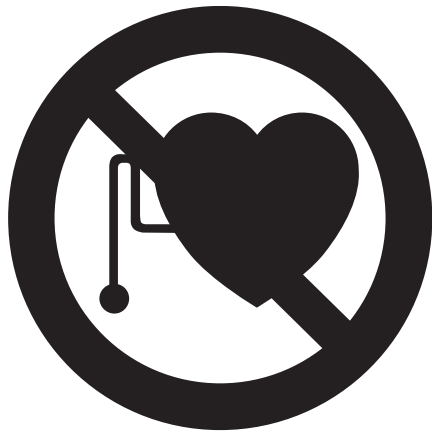

Strong magnets can cause serious injury to people with Pacemakers or Automatic Internal Cardiac Defibrillators.

Strong magnets can attract tools and metal objects, causing pinch, cut, or entrapment hazards.

Keep tools, metal objects, and magnetic media away from magnets.

