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B- CLOCK SET UP

- Press CLOCK key about 3 seconds until screen gleams.
- Press HR/MIN/SEC keys to set time. Leave press to accelerate the set up.
- Press CLOCK key to confirm

C- ALARM SET UP

- Press ALARM key about 3 seconds until screen gleams.
- Press HR/MIN/SEC keys to set time. Leave press to accelerate the set up.
- Press ALARM key to confirm

D- COUNT UP 1 AND 2 SET UP

- Press TIMER 1 or TIMER 2 key to set up count up 1 or 2
- Press CLEAR key to restore.
- Press START/STOP key to start the count up. Press again START/STOP to stop count down

Note : 422CA can count up on TIMER 1 and TIMER 2 simultaneously and independently

E- COUNT DOWN 1 AND 2 SET UP

- Press TIMER 1 or TIMER 2 key to set up count down 1 or 2
- Press CLEAR key to restore.
- Press HR/MIN/SEC keys to set time to count down. Leave press to accelerate the set up.
- Presser START/STOP key to start the count down.
- At count down's end, alarm rings during 1min. Time's up start and display until 23:59 99, then return to 0:00 00
- Press START/STOP key to stop the count down.
- Anytime, press CLOCK or ALARM key to return to CLOCK or ALARM display.

Note : 422CA can count down on TIMER 1 and TIMER 2 simultaneously and independently.

422CA can count up on one TIMER and count down on the other TIMER.

F- MEMORY FUNCTION

- Press MEMORY key to memorize the TIMER's set up.
- Press TIMER 1 or TIMER 2 key, then press MEMORY key to display the memory time.

G- SNOOZER FUNCTION

- Alarm rings during 1mn if it is not stopped by pressing any keys, then rings again after 5mn : SNOOZER function, « Zz » display on screen
- Press ALM ON/OFF key to definitely stop alarm.

Warning : deposit material and worn battery in an ecologically suitable place

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