

# When downstream quality matters, make sure your upstream tools are the best.

## Amicon® Ultra 2 mL Centrifugal Filters for Protein Purification and Concentration

The Amicon® Ultra 2 mL filter provides a fast and easy way to concentrate, desalt and exchange the buffer for both proteins and nucleic acids. Now you can concentrate 2 mL down to 15 – 70  $\mu$ L in an Amicon® Ultra centrifugal device.

By incorporating familiar Amicon® Ultra features such as the vertical membrane design and true dead stop, the Amicon® Ultra 2 mL filter delivers high recovery, short spin times and quality results even for the most dilute samples.

### Amicon® Ultra 2 mL Advantages

- True engineered dead stop provides consistent recovery without spinning to dryness
- Reverse spin aids in sample recovery, especially with small, dilute samples
- Vertical membrane design aligns with the filtrate for less clogging and faster filtering
- Ultracel® low-binding membrane combines with the innovative vertical housing to provide fast sample processing without sacrificing recovery
- Fast processing, with 10- to 30-minute spin times



## Easy to Choose!

Select the Amicon® Ultra Filter that best meets your sample needs. Choose your starting volume, identify the molecular weight of your molecule of interest, and identify final volume expected.


### Amicon® Ultra Centrifugal Filters

| Product                                   | Amicon® Ultra-0.5 | Amicon® Ultra-2 | Amicon® Ultra-4 | Amicon® Ultra-15 |
|-------------------------------------------|-------------------|-----------------|-----------------|------------------|
| Maximum initial sample volume (mL)        | 0.5               | 2               | 4               | 15               |
| Final concentrate (retentate) volume (µL) | 15 – 20           | 15 – 70         | 30 – 70         | 150 – 300        |

| MWCO         | Qty/Pk |           |           |           |           |
|--------------|--------|-----------|-----------|-----------|-----------|
| 3,000 MWCO   | 8      | UFC500308 |           | UFC800308 | UFC900308 |
|              | 24     | UFC500324 | UFC200324 | UFC800324 | UFC900324 |
|              | 96     | UFC500396 |           | UFC800396 | UFC900396 |
|              | 500    | UFC5003BK |           |           |           |
| 10,000 MWCO  | 8      | UFC501008 |           | UFC801008 | UFC901008 |
|              | 24     | UFC501024 | UFC201024 | UFC801024 | UFC901024 |
|              | 96     | UFC501096 |           | UFC801096 | UFC901096 |
|              | 500    | UFC5010BK |           |           |           |
| 30,000 MWCO  | 8      | UFC503008 |           | UFC803008 | UFC903008 |
|              | 24     | UFC503024 | UFC203024 | UFC803024 | UFC903024 |
|              | 96     | UFC503096 |           | UFC803096 | UFC903096 |
|              | 500    | UFC5030BK |           |           |           |
| 50,000 MWCO  | 8      | UFC505008 |           | UFC805008 | UFC905008 |
|              | 24     | UFC505024 | UFC205024 | UFC805024 | UFC905024 |
|              | 96     | UFC505096 |           | UFC805096 | UFC905096 |
|              | 500    | UFC5050BK |           |           |           |
| 100,000 MWCO | 8      | UFC510008 |           | UFC810008 | UFC910008 |
|              | 24     | UFC510024 | UFC210024 | UFC810024 | UFC910024 |
|              | 96     | UFC510096 |           | UFC810096 | UFC910096 |
|              | 500    | UFC5100BK |           |           |           |

Search with our online Amicon® Ultra Selector Tool and find the right filter to concentrate your sample. Visit: [www.millipore.com/FastEasy](http://www.millipore.com/FastEasy)

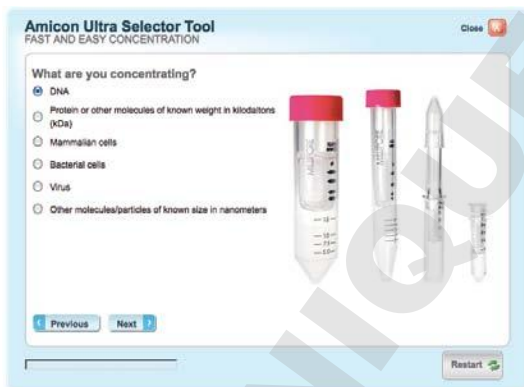

## To Place an Order or Receive Technical Assistance

In Europe, please call Customer Service:

France: 0825 045 645

Germany: 01805 045 645

Italy: 848 845 645

Spain: 901 516 645 Option 1

Switzerland: 0848 645 645

United Kingdom: 0870 900 4645

For other countries across Europe, please call: +44 (0) 115 943 0840

Or visit: [www.merckmillipore.com/offices](http://www.merckmillipore.com/offices)

For Technical Service visit:

[www.merckmillipore.com/techservice](http://www.merckmillipore.com/techservice)

## Get Connected!

Join Merck Millipore Bioscience on your favorite social media outlet for the latest updates, news, products, innovations, and contests!

 [facebook.com/MerckMilliporeBioscience](https://facebook.com/MerckMilliporeBioscience)

 [twitter.com/Merck4Bio](https://twitter.com/Merck4Bio)



[www.merckmillipore.com](http://www.merckmillipore.com)

Merck Millipore and the M logo are trademarks of Merck KGaA, Darmstadt, Germany. Amicon and Ultracel are registered trademarks of Merck KGaA, Darmstadt, Germany. Lit. No. DS3336ENEU 3/2012 LS-SBU-12-06120 Printed in the USA. © 2012 EMD Millipore Corporation, Billerica, MA, USA. All rights reserved.