

# When downstream quality matters, make sure your upstream tools are the best.

## Amicon® Ultra 2 mL Centrifugal Filters for Protein Purification and Concentration

The Amicon® Ultra 2 mL filter provides a fast and easy way to concentrate, desalt and exchange the buffer for both proteins and nucleic acids. Now you can concentrate 2 mL down to 15 – 70  $\mu$ L in an Amicon® Ultra centrifugal device.

By incorporating familiar Amicon® Ultra features such as the vertical membrane design and true dead stop, the Amicon® Ultra 2 mL filter delivers high recovery, short spin times and quality results even for the most dilute samples.

### Amicon® Ultra 2 mL Advantages

- True engineered dead stop provides consistent recovery without spinning to dryness
- Reverse spin aids in sample recovery, especially with small, dilute samples
- Vertical membrane design aligns with the filtrate for less clogging and faster filtering
- Ultracel® low-binding membrane combines with the innovative vertical housing to provide fast sample processing without sacrificing recovery
- Fast processing, with 10- to 30-minute spin times



## Easy to Choose!

Select the Amicon® Ultra Filter that best meets your sample needs. Choose your starting volume, identify the molecular weight of your molecule of interest, and identify final volume expected.


### Amicon® Ultra Centrifugal Filters

Product	Amicon® Ultra-0.5	Amicon® Ultra-2	Amicon® Ultra-4	Amicon® Ultra-15
Maximum initial sample volume (mL)	0.5	2	4	15
Final concentrate (retentate) volume (µL)	15 – 20	15 – 70	30 – 70	150 – 300

MWCO	Qty/Pk				
3,000 MWCO	8	UFC500308		UFC800308	UFC900308
	24	UFC500324	UFC200324	UFC800324	UFC900324
	96	UFC500396		UFC800396	UFC900396
	500	UFC5003BK			
10,000 MWCO	8	UFC501008		UFC801008	UFC901008
	24	UFC501024	UFC201024	UFC801024	UFC901024
	96	UFC501096		UFC801096	UFC901096
	500	UFC5010BK			
30,000 MWCO	8	UFC503008		UFC803008	UFC903008
	24	UFC503024	UFC203024	UFC803024	UFC903024
	96	UFC503096		UFC803096	UFC903096
	500	UFC5030BK			
50,000 MWCO	8	UFC505008		UFC805008	UFC905008
	24	UFC505024	UFC205024	UFC805024	UFC905024
	96	UFC505096		UFC805096	UFC905096
	500	UFC5050BK			
100,000 MWCO	8	UFC510008		UFC810008	UFC910008
	24	UFC510024	UFC210024	UFC810024	UFC910024
	96	UFC510096		UFC810096	UFC910096
	500	UFC5100BK			

Search with our online Amicon® Ultra Selector Tool and find the right filter to concentrate your sample. Visit: [www.millipore.com/FastEasy](http://www.millipore.com/FastEasy)

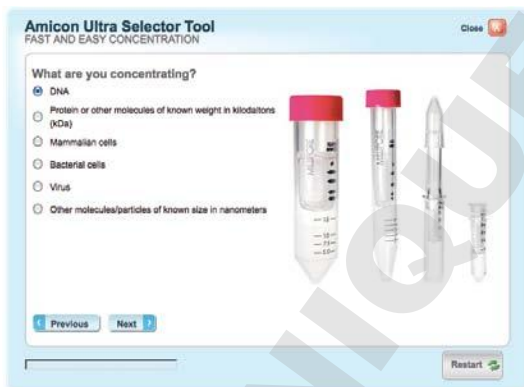

## To Place an Order or Receive Technical Assistance

In Europe, please call Customer Service:

France: 0825 045 645

Germany: 01805 045 645

Italy: 848 845 645

Spain: 901 516 645 Option 1

Switzerland: 0848 645 645

United Kingdom: 0870 900 4645

For other countries across Europe, please call: +44 (0) 115 943 0840

Or visit: [www.merckmillipore.com/offices](http://www.merckmillipore.com/offices)

For Technical Service visit:

[www.merckmillipore.com/techservice](http://www.merckmillipore.com/techservice)

## Get Connected!

Join Merck Millipore Bioscience on your favorite social media outlet for the latest updates, news, products, innovations, and contests!

 [facebook.com/MerckMilliporeBioscience](https://facebook.com/MerckMilliporeBioscience)

 [twitter.com/Merck4Bio](https://twitter.com/Merck4Bio)



[www.merckmillipore.com](http://www.merckmillipore.com)

Merck Millipore and the M logo are trademarks of Merck KGaA, Darmstadt, Germany. Amicon and Ultracel are registered trademarks of Merck KGaA, Darmstadt, Germany. Lit. No. DS3336ENEU 3/2012 LS-SBU-12-06120 Printed in the USA. © 2012 EMD Millipore Corporation, Billerica, MA, USA. All rights reserved.