

## CHEMICAL RESISTANCE DATA



# SHIELDskin XTREME™

## SHIELDskin Xtreme White Nitrile 300 DI



- Cleanroom glove - PPE Complex Design (Category III)
- 100% Nitrile glove - Powder Free
- Unique Skin Nitrile Technology
- Ambidextrous -Textured finger tips
- AQL 1.5 (EN374-2:2003 Level 2)
- Length 300 mm
- Reduced allergy risks (Type IV) - 100% Latex free
- Typical particle levels: less than 3000 per cm<sup>2</sup> more or equal 0,5µm

1239-45-8	Ethidium bromide 5%	Breakthrough time : 480 minutes	Class 6
1310-73-2	Sodium Hydroxide, 40%	Breakthrough time : 480 minutes	Class 6
50-00-0	Formaldehyde, 37%	Breakthrough time : 480 minutes	Class 6
75-59-2	Tetramethylammonium hydroxide 2.5%	Breakthrough time : 480 minutes	Class 6
7664-93-9	Sulphuric acid, 50%	Breakthrough time : 480 minutes	Class 6
79-06-1	Acrylamide 40%	Breakthrough time : 480 minutes	Class 6
7647-01-0	Hydrochloric Acid, 37%	Breakthrough time : 129 minutes	Class 4
67-63-0	Isopropanol 70%	Breakthrough time : 65 minutes	Class 3
67-63-0	Isopropanol 100%	Breakthrough time : 49 minutes	Class 2
71-36-3	Butanol 100%	Breakthrough time : 43 minutes	Class 2
7722-84-1	Hydrogen Peroxide, 30%	Breakthrough time : 39 minutes	Class 2
64-17-5	Ethanol, 70%	Breakthrough time : 33 minutes	Class 2
1336-21-6	Ammonium Hydroxide, 25%	Breakthrough time : 29 minutes	Class 1
64-17-5	Ethanol 100%	Breakthrough time : 17 minutes	Class 1
7664-39-3	Hydrofluoric Acid, 40%	Breakthrough time : 11 minutes	Class 1
64-19-7	Acetic Acid 100%	Breakthrough time : 7 minutes	Class 0
108-94-1	Cyclohexanone 99%	Breakthrough time : 6 minutes	Class 0