

CHEMICAL RESISTANCE DATA



SHIELDskin XTREME™

SHIELDskin Xtreme White Nitrile 300 DI++



- Cleanroom glove - PPE Complex Design (Category III)
- 100% Nitrile glove - Powder Free
- Unique Skin Nitrile Technology
- Ambidextrous -Textured finger tips
- AQL 1.5 (EN374-2:2003 Level 2)
- Length 300 mm
- Reduced allergy risk (Type IV) - Latex Free glove
- Typical particle levels: Less than 650 per cm² more or equal 0,5µm

1239-45-8	Ethidium bromide 5%	Breakthrough time : 480 minutes	Class 6
1310-58-3	Potassium Hydroxyde 40%	Breakthrough time : 480 minutes	Class 6
1310-73-2	Sodium Hydroxide, 40%	Breakthrough time : 480 minutes	Class 6
50-00-0	Formaldehyde, 37%	Breakthrough time : 480 minutes	Class 6
75-59-2	Tetramethylammonium hydroxide 2.5%	Breakthrough time : 480 minutes	Class 6
7664-93-9	Sulphuric acid, 50%	Breakthrough time : 480 minutes	Class 6
79-06-1	Acrylamide 40%	Breakthrough time : 480 minutes	Class 6
7647-01-0	Hydrochloric Acid, 37%	Breakthrough time : 134 minutes	Class 4
67-63-0	Isopropanol 70%	Breakthrough time : 76 minutes	Class 3
67-63-0	Isopropanol 100%	Breakthrough time : 62 minutes	Class 3
7722-84-1	Hydrogen Peroxide, 30%	Breakthrough time : 52 minutes	Class 2
64-17-5	Ethanol, 70%	Breakthrough time : 35 minutes	Class 2
1336-21-6	Ammonium Hydroxide, 25%	Breakthrough time : 32 minutes	Class 2
64-17-5	Ethanol 100%	Breakthrough time : 20 minutes	Class 1
7664-39-3	Hydrofluoric Acid, 40%	Breakthrough time : 13 minutes	Class 1
67-56-1	Methanol 99,9%	Breakthrough time : 7 minutes	Class 0
1330-20-7	Xylene 98,5%	Breakthrough time : 1 minutes	Class 0