

## CHEMICAL RESISTANCE DATA



# SHIELDskin XTREME™

## SHIELDskin Xtreme White Nitrile 400 DI+



- Cleanroom glove - PPE Complex Design (Category III)
- 100% Nitrile glove - Powder Free
- Unique Skin Nitrile Technology
- Ambidextrous - Textured finger tips
- AQL 1.5 (EN374-2:2003 Level 2)
- Length 400 mm
- Reduced allergy risks (Type IV) - Latex free glove
- Virus & Microorganism resistant (ASTM F1671-97b and EN374-2:2003)
- Typical Particle level: less than 950 per cm<sup>2</sup> more or equal 0.5µm

1239-45-8	Ethidium bromide 5%	Breakthrough time : 480 minutes	Class 6
1310-58-3	Potassium Hydroxyde 40%	Breakthrough time : 480 minutes	Class 6
1310-73-2	Sodium Hydroxide, 40%	Breakthrough time : 480 minutes	Class 6
50-00-0	Formaldehyde, 37%	Breakthrough time : 480 minutes	Class 6
7664-93-9	Sulphuric acid, 50%	Breakthrough time : 480 minutes	Class 6
79-06-1	Acrylamide 40%	Breakthrough time : 480 minutes	Class 6
7647-01-0	Hydrochloric Acid, 37%	Breakthrough time : 157 minutes	Class 4
67-63-0	Isopropanol 70%	Breakthrough time : 95 minutes	Class 3
67-63-0	Isopropanol 100%	Breakthrough time : 72 minutes	Class 3
7722-84-1	Hydrogen Peroxide, 30%	Breakthrough time : 59 minutes	Class 2
64-17-5	Ethanol, 70%	Breakthrough time : 46 minutes	Class 2
1336-21-6	Ammonium Hydroxide, 25%	Breakthrough time : 32 minutes	Class 2
64-17-5	Ethanol 100%	Breakthrough time : 22 minutes	Class 1
7664-39-3	Hydrofluoric Acid, 40%	Breakthrough time : 15 minutes	Class 1