

## CHEMICAL RESISTANCE DATA



# SHIELDskin XTREME™

## SHIELDskin Xtreme Sterile Nitrile 330 DI+



- Cleanroom glove - PPE Complex Design (Category III)
- 100% Nitrile glove - Powder Free - Sterile glove
- Hand Specific - Fully Textured
- AQL 0.65 (EN374-2:2003 Level 3)
- Length 330 mm
- Reduced allergy risk (Type IV) - Latex Free glove
- Virus and Microorganism resistant (ASTM F1671-97b and EN374-2:2003)
- Typical particle levels: less than 950 per cm<sup>2</sup> more or equal 0.5µm

## CHEMICAL RESISTANCE DATA



# SHIELDskin XTREME™

### SHIELDskin Xtreme Sterile Nitrile 330 DI+

111-30-8	Glutaraldehyde 25%	Breakthrough time : 480 minutes	Class 6
1239-45-8	Ethidium bromide 5%	Breakthrough time : 480 minutes	Class 6
1310-73-2	Sodium Hydroxide, 40%	Breakthrough time : 480 minutes	Class 6
50-00-0	Formaldehyde, 37%	Breakthrough time : 480 minutes	Class 6
7664-93-9	Sulphuric acid, 50%	Breakthrough time : 480 minutes	Class 6
7681-52-9	Sodium Hypochlorite 13%	Breakthrough time : 480 minutes	Class 6
79-06-1	Acrylamide 40%	Breakthrough time : 480 minutes	Class 6
67-63-0	Isopropanol 70%	Breakthrough time : 57 minutes	Class 2
67-63-0	Isopropanol 100%	Breakthrough time : 44 minutes	Class 2
7722-84-1	Hydrogen Peroxide, 30%	Breakthrough time : 41 minutes	Class 2
64-17-5	Ethanol, 70%	Breakthrough time : 37 minutes	Class 2
1336-21-6	Ammonium Hydroxide, 25%	Breakthrough time : 27 minutes	Class 1
142-82-5	Heptane	Breakthrough time : 25 minutes	Class 1
64-17-5	Ethanol 100%	Breakthrough time : 16 minutes	Class 1
7664-93-9	Sulphuric acid, 95-98%	Breakthrough time : 6 minutes	Class 0
1330-20-7	Xylene 98,5%	Breakthrough time : 3 minutes	Class 0
67-64-1	Acetone 100%	Breakthrough time : 1 minutes	Class 0
75-05-8	Acetonitrile 99,9%	Breakthrough time : 1 minutes	Class 0
80-62-6	Methyl Methacrylate 99%	Breakthrough time : 1 minutes	Class 0
109-89-7	Diethylamine 99,5%	Breakthrough time : 0 minutes	Class 0
75-09-2	Dichloromethane 99%	Breakthrough time : 0 minutes	Class 0