

Level 0
0 - <10 min

Level 1
10 - 30 min

Level 2
30 - 60 min

Level 3
60 - 120 min

Level 4
120 - 240 min


Level 5
240 - 480 min

Level 6
< 480 min

SHIELDskin XTREME* Eco Nitrile 300 DI+



0120

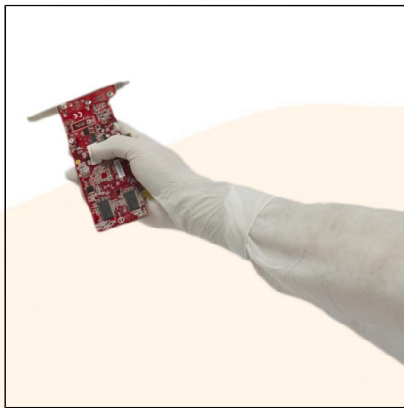

- Category III PPE glove (PPE Directive 89/686/EEC)
- Complex Design - For mortal & irreversible risks
- Powder-free white nitrile glove
- Ambidextrous
- 300 mm / 0.10 mm (EN 420:2003 + A1:2009)
- AQL 1.5 (EN 374-2:2014 Level 2)
- Viral penetration test (ISO 16604:2004 Procedure B & ASTM F1671-97b)
- Waterproof and for low chemical protection
- Tested for chemical permeation (EN 16523-1:2015 supersedes EN 374-3:2003)
- Typical particle levels: less than 900 per cm² more or equal 0.5µm
- Tested according to EN 1149-1-2-3 & 5
- Type I hypersensitivity eliminated - Type IV hypersensitivity reduced

7664-93-9
Sulphuric acid 50%

LEVEL 6
480 min

108-95-2
Phenol 0.1% solution

LEVEL 6
480 min

75-59-2
Tetramethylammonium hydroxide 2.5%

LEVEL 6
480 min

1310-58-3
Potassium Hydroxide 40%

LEVEL 6
480 min

1310-73-2
Sodium Hydroxide, 50%

LEVEL 6
480 min

7722-84-1
Hydrogen Peroxide 30%

LEVEL 6
480 min

50-00-0
Formaldehyde 37%

LEVEL 6
480 min

1239-45-8 Ethidium bromide 5%	LEVEL 6 480 min
7647-01-0 Hydrochloric Acid conc.	LEVEL 3 79 min
7647-01-0 Hydrochloric Acid 37%	LEVEL 3 74 min
79-06-1 Acrylamide 40%	LEVEL 6 480 min
67-63-0 Isopropanol 70%	LEVEL 2 45 min
142-82-5 n-Heptane 99%	LEVEL 2 41 min
67-63-0 Isopropanol 100%	LEVEL 2 35 min
64-17-5 Ethanol 70%	LEVEL 1 22 min
64-17-5 Ethanol 99.8%	LEVEL 1 11 min
67-68-5 Dimethyl Sulfoxide 99% (DMSO)	LEVEL 0 9 min
7664-39-3 Hydrofluoric Acid 40%	LEVEL 0 8 min
7697-37-2 Nitric acid 65%	LEVEL 0 5 min
1330-20-7 Xylene 98,5%	LEVEL 0 2 min
1336-21-6 Ammonium Hydroxide 25%	LEVEL 1 16 min

68-12-2 Dimethyl Formamide 99,8%	LEVEL 0 1 min
67-64-1 Acetone 99,8%	LEVEL 0 0 min
67-56-1 Methanol 99,9%	LEVEL 0 0 min
141-78-6 Ethyl Acetate 99.8%	LEVEL 0 0 min
109-89-7 Diethylamine 99,5%	LEVEL 0 0 min
67-66-3 Chloroform 99,8%	LEVEL 0 0 min
75-05-8 Acetonitrile 99,9%	LEVEL 0 0 min

DOMINIQUE DUTSCHER SAS