


# pipetman<sup>®</sup>

DIAMOND TIPS

## Refill for TOWERPACK boxes

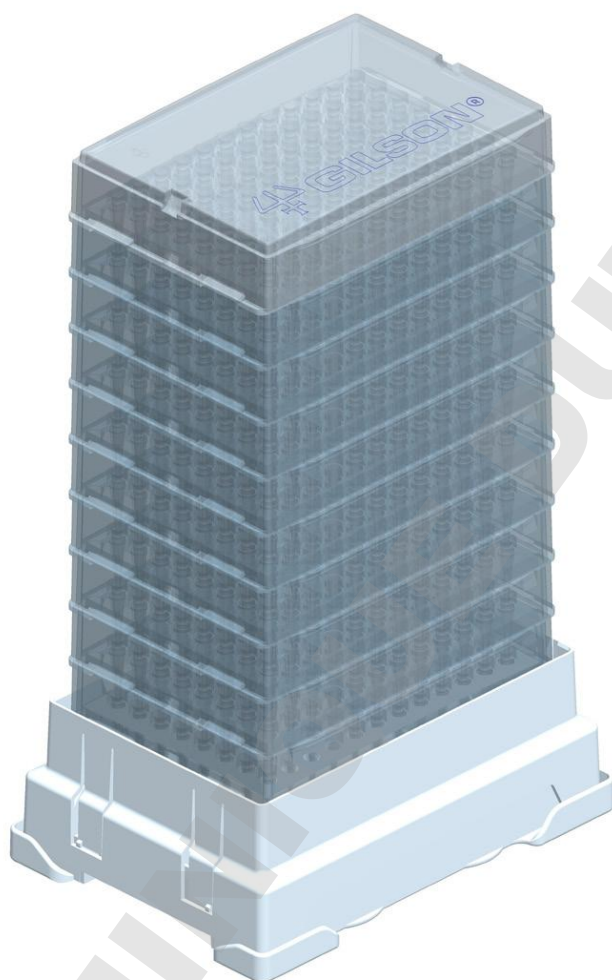

- Uses 60 % less space and 70 % less plastic.
- One reload box for the entire TOWERPACK range.
- Easy-to-use one-handed system.
- Protective base to prevent contamination from environmental pollution and dust.
- Available in standard and sterilized versions.

Loading instructions on the reverse side.



Made in France

Download certificates at [www.gilson.com](http://www.gilson.com)

**Gilson S.A.S.**


19, avenue des Entrepreneurs BP 145 • F-95400  
Villiers-le-Bel, France  
T +33 (0) 1 34 29 50 00 • F +33 (0) 1 34 29 50 20

[www.gilson.com/contactus](http://www.gilson.com/contactus)

 **GILSON<sup>®</sup>**

## Loading a TOWER PACK Box

1. Remove the TOWER PACK dispenser from the protective support.
2. Place the TOWER PACK on the opened reload box.
3. Press down firmly.
4. The rack snaps into place.
5. When a rack is empty, remove it from the reload box, and load a new rack.

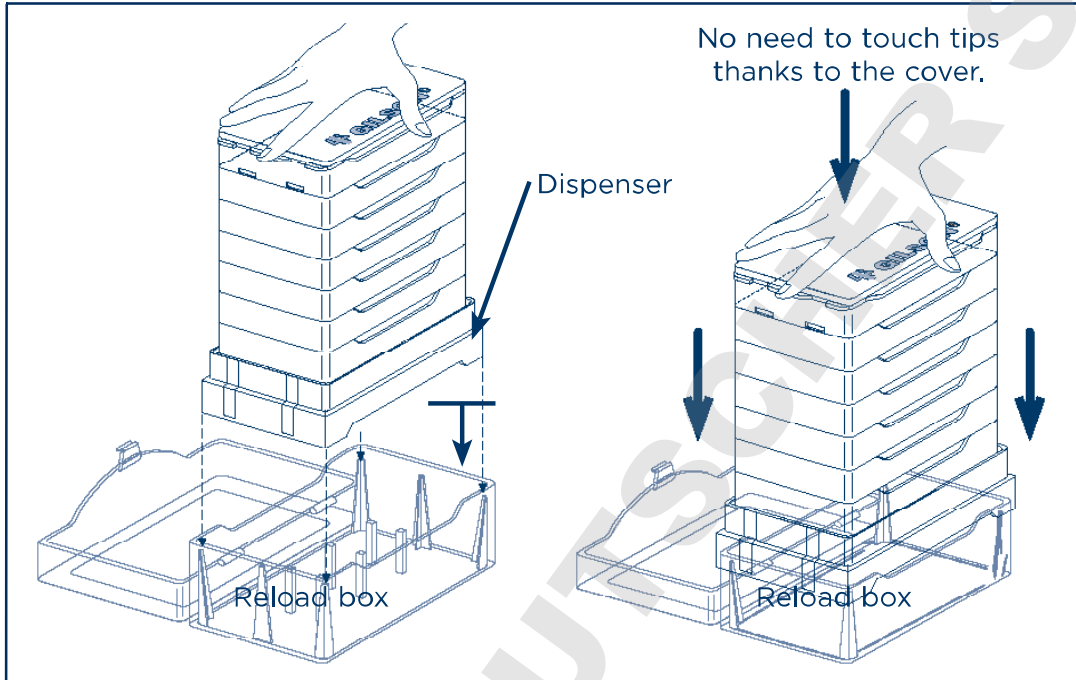

## Autoclaving

The entire TOWER PACK can be autoclaved or just the reload box filled with one rack of tips.

Autoclaving conditions: moist heat, 121°C, 0.1 MPa for 20 minutes).

### NOTE

When autoclaving PIPETMAN DIAMOND D10 tips, take care to keep the box upright to prevent these tips from falling out of the rack.

## Loading a Multichannel PIPETMAN

The patented Rocky Rack™ makes it extremely easy to fit PIPETMAN DIAMOND tips on a multichannel PIPETMAN.

Exert a light vertical force to obtain a perfect seal on all eight or twelve channels.

