

The JIMMY™

Your thumb's best friend!

FR téléchargeable sur www.gilson.com

DU von www.gilson.com downloaden

IT scaricabili su www.gilson.com

SP se puede telecargar desde www.gilson.com



PIPETMAN® ergonomic and hands-free microtube opener:

- ▮ Opens both snap-cap and screw-cap tubes.
- ▮ Increases pipetting efficiency.
- ▮ Lessens the fatigue associated with manual tube opening.

Product ref. F144983

 **GILSON** 

QUICK GUIDE

This simple yet highly effective device eliminates the need to manually handle the caps of tubes when opening.

Clipping The JIMMY™

- 1 Rest the pipette underneath the two tabs on the upper surface of the prongs.



- 2 While holding the pipette stationary, rotate The JIMMY™ down and toward the pipette until it «snaps» into place.



- 3 To remove, pull straight off.

Cleaning and Sterilization

Made of high quality polyacetal, The JIMMY™ can be cleaned with alcohol or sterilized by autoclaving.

Opening Snap-cap Tubes



1 Insert the tab into the slot.



2 Move the pipette upwards to open the tube.

Opening Screw-cap Tubes



1 Insert the cap into the socket.



2 Once the cap is secure, rotate the tube to remove the cap.

3 To close, return the tube to the secured cap, and rotate the tube until it is tight and remove.

Key features and benefits

2 TOOLS IN 1:

- You can open both snap-cap tubes and screw-cap tubes.
- You can leave the screw-cap in the opener and still open snap-cap tubes with The JIMMY's™ 2 separate tools in 1.

PIPETTE ADAPTABILITY:

- You can fit all PIPETMAN and PIPETMAN Neo models: P2 to P10 mL, P2N to P1000N.

CONTAMINATION RISK FREE:

- You can keep your fingers and pipette free of microtube contents, due to The JIMMY's™ specific design and position on the pipette.

PERSONALIZATION FACILITIES:

- You can personalize The JIMMY™ with a permanent ink, due to its high quality polyacetal plastic.
- You can identify your applications with The JIMMY's™ set of 3 colors: orange, yellow and white.



www.gilson.com | sales@gilson.com | service@gilson.com | training@gilson.com

Gilson, Inc. World Headquarters

3000 Parmenter Street | P.O. Box 620027 | Middleton, WI 53562-0027 USA
Tel: (1) 800-445-7661 or (1) 608-836-1551 | Fax: (1) 608-831-4451

Gilson S.A.S. | 19, avenue des Entrepreneurs | BP 145, F-95400 Villiers-le-Bel, France

Tel: (33-1) 34 29 50 00 | Fax: (33-1) 34 29 50 20