



CORIO™ CD



english

Immersion / Open Bath / Refrigerated Circulators

Operating Instructions

Important: keep original operating manual for future use.

1.951.0800-V1

Table of contents

1 CORIO® CD product overview	5
2 Intended use	5
3 Description	6
4 Explanation of safety information	7
4.1 Explanation of other information	7
5 Safety instructions	8
6 Operator's responsibility - safety instructions	10
6.1 Requirements for the operating personnel.....	10
6.2 Operating and ambient conditions for using the unit.....	10
6.3 Operating the unit.....	11
7 Control and function elements	13
7.1 Installation of the circulator	16
8 Preparations for operating the device	16
8.1 Securing the immersion circulator.....	16
8.2 Open bath tanks.....	19
9 Closed stainless steel bath tanks	22
9.1 Bath fluids.....	23
Water bath fluid.....	23
9.2 Temperature control for external connected systems.....	25
9.3 Tubing.....	26
10 Commissioning	27
10.1 Excess temperature and low level safety devices	28
10.2 Filling	28
10.3 Pump settings	29
10.4 Switching on / Start - Stop	29
10.5 Adjusting of temperature setpoint	30
10.6 ATC - Absolute Temperature Calibration (1-point calibration)	31
10.7 Switch Autostart on and off	32
11 Error messages / Possible causes of faults	33

12 Operation of the CORIO CD via the USB interface	36
13 Emptying the bath tank.....	38
14 Technical data	39
14.1 Technical data for circulator	39
14.2 Technical data for refrigeration circulation circulator.....	40
15 Materials of parts in contact with the bath fluid.....	42
15.1 Circulator	42
16 Accessories	43
16.1 For external connection.....	43
16.2 For open baths	43
17 Maintenance, cleaning, storage.....	43
17.1 Maintain the refrigeration capacity.....	43
17.2 Cleaning.....	44
18 Storage	45
19 Repair service.....	45
19.1 Warranty	45
20 Waste disposal	46
20.1 Packaging.....	46
20.2 Unit	46
20.3 Refrigerant	46
21 EC conformity.....	47

Congratulations.

You have made an excellent choice.

JULABO thanks you for the trust you have placed in us.

This operating manual is designed to familiarize you with the operation of our units and their possible applications. Please read the operating manual carefully.

Please call us if you have any questions about the operation of the unit or about the operating manual.

Contact:

JULABO GmbH
Eisenbahnstrasse 45
77960 Seelbach / Germany
Tel. +49 (0) 7823 / 51-0
Fax +49 (0) 07823 / 2491
info.de@julabo.com
www.julabo.com

The JULABO quality management system

The standards for the development, production and distribution of temperature control devices for laboratory and industry use satisfy the requirements of ISO 9001 and ISO 14001.
Registration certificate No. 01 100044846.



Unpacking and inspection

If the packaging is damaged or if you discover any concealed transport damage when you have unpacked the devices and the accessories, please notify the supplier in the form of a statement of damage.

NOTICE

The operating manual

- should be kept for future use.
- must be available to operating personnel at all times.

DOMINIQUE DUTSCHER SAS