

Economy Counter

HS5360A / HS5360B

Operating Instructions:

- 1) Place counter on even surface.
- 2) Turn knob to zero all counters.
- 3) Press key firmly for each count. Each key has independent counter. Bell will ring once the Total window reaches each hundred total.
- 4) Turn knob to re-zero counter for next project.

FAQ

- 1) The counter won't reset!
Try turning the knob slowly.
- 2) The values on the counter are not turning!
Press the buttons firmly all the way down.
- 3) The casing arrived damaged!
We have extras in the sample room! Do not order a whole counter!