STERILBIO

3000788

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

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CONTENTS

GENERAL INFORMATION
 CONTENTS
 PACKING LISTS
 TECHNICAL FEATURES
 EQUIPMENT DESCRIPTION
 INSTALLATION
 OPERATION
 GUARANTEE
 "EC" CONFORMITY DECLARATION

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PACKING LIST

The standard equipment consist of the following components:

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TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Potencia eléctrica consumida / Consumption	Peso / Weight	Medidas exteriores cm / Overall dimensions cm			Medidas interiores tubo cuarzo cm / Internal dimensions quartz tube cm	
		Alto / Height	Ancho / Length	Fondo / Width	∅	Largo / Length
110 W	1.5 Kg	20	13	18	1.4	15

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EQUIPMENT DESCRIPTION

Infra-red incinerators STERILBIO, are made of ceramic parts with internal tube of high quality quartz. No oxygen required and can be used in anaerobics environments. Eliminates problems of aerosols. Avoids bacterial contamination resulting from flame methods of sterilization. Side stand with four housing include for Kolle handles, needles, lances, dissecting forceps, etc...

Effective incineration of culture handles, needles, culture tubes and small instruments and of all organic material in 5/8 seconds at a temperature of 900°C.

INSTALLATION

It is not necessary any special installation, simply plug it to the mains.

Important:

During the first starting up, the apparatus can gives off some fume because of the initial warming of the materials. That circumstance is completely normal and will disappear in a period of time.

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Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the apparatus.

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Do not use the apparatus if it is NOT earthed.

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If you have to change the plug bear in mind the following:

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For the starting up, you only have to press the main switch.

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After 7 or 8 minutes a temperature of 900°C will be reached inside the tube and the apparatus is ready for use.

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To sterilize an instrument, introduce it by the central tube, and after 5/8 seconds, the end of the instrument will be sterilized because of the high temperature.

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The sterilizer can be on without stopping for about 3 or 4 hours, because of the temperature stops raising when reaches 950°C. But if it is not necessary, we recommend disconnect it after using and foresee the connection 10 minutes before the use. Doing that, the live of the sterilizer will be longer.

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Caution:

The ceramic tube envelope can reach temperatures of about 80°C

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GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of J.P. SELECTA, S.A..

Any manipulation of the apparatus by unauthorized personnel by J.P. SELECTA, S.A. cancels the guarantee automatically.

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DOMINIQUE DUTSCHER SAS