PLAC-CENTER

44 000 501

PLAQUE REFRIGERANTE

COOLING PLATE

INFORMATION GENERALE

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

SOMMAIRE

INFORMATION GENERALE
SOMMAIRE
LISTE D'EMBALLAGE
SPECIFICATIONS TECHNIQUES
DESCRIPTION DE L'EQUIPEMENT
INSTALLATION
MODE OPERATOIRE
ENTRETIEN
GARANTIE
DECLARATION DE CONFORMITE "CE"

CONTENTS

GENERAL INFORMATION
CONTENTS
PACKING LISTS
TECHNICAL FEATURES
EQUIPMENT DESCRIPTION
INSTALLATION
OPERATION
MAINTENANCE
GUARANTEE
"EC" CONFORMITY DECLARATION

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PACKING LIST

The standard equipment consist of the following components:

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TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Code / Code		44000501
Dimensions utiles plaque cm / Plate dimensions cm.	Ancho / Length	30
	Fondo / Width	15
Dimensions extérieures cm. / Overall dimensions cm	Haut. / Height	27
	Larg. / Length	33
	Prof. / Width	50
Puissance compresseur H/P Compressor power HP		1 / 3
Consommation W. Consumption W		160
Poids Kg. / Weight Kg.		15

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EQUIPMENT DESCRIPTION

The cooling plate PLAC-CENTER is specially suitable for maintaining substances in solid state.

For adjustable temperatures on plate from 5°C up to -10°C, with electronic regulation and digital reading. AISI 304 stainless steel outer casing and cooling plate in black anodized aluminium. Hermetically-sealed compressor installed on an anti-vibration mounting with ventilated evaporator, regulating valve and condenser.

Control panel equipped with main switch with indicator lamp, analogue electronic temperature regulator and digital reading (2 digits) with PTC probe and resolution ±1 digit.

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INSTALLATION

Place the cooling plate on a flat horizontal, level surface, leaving a space at the back and sides of the apparatus of about 10cm.

**ATTENTION ! IMPORTANT POUR VOTRE SECURITE
CAUTION!!! IMPORTANT FOR YOUR SAFETY!**

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Do NOT use the plate for thermic treatment of products which made form explosive gases, not even in special protection aerea.

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Be sure that the voltage supply is the same as the one indicated on the characteristics plate.

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Do not use the apparatus if it is NOT earthed.

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If you have to change the plug bear in mind the following:

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*Blue cable: Neutral.
Brown cable: Phase.
Yellow/green cable: Earth.*

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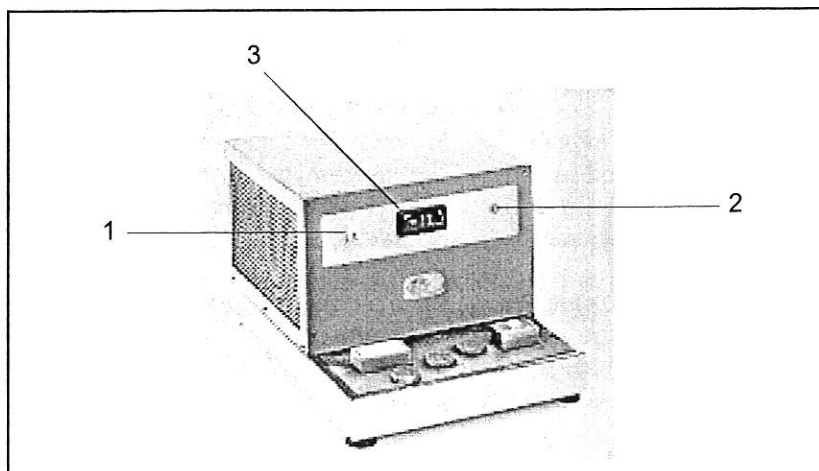
1. Main switch with indicator lamp.
2. Compressor operation indicator lamp.
3. Temperature controller.

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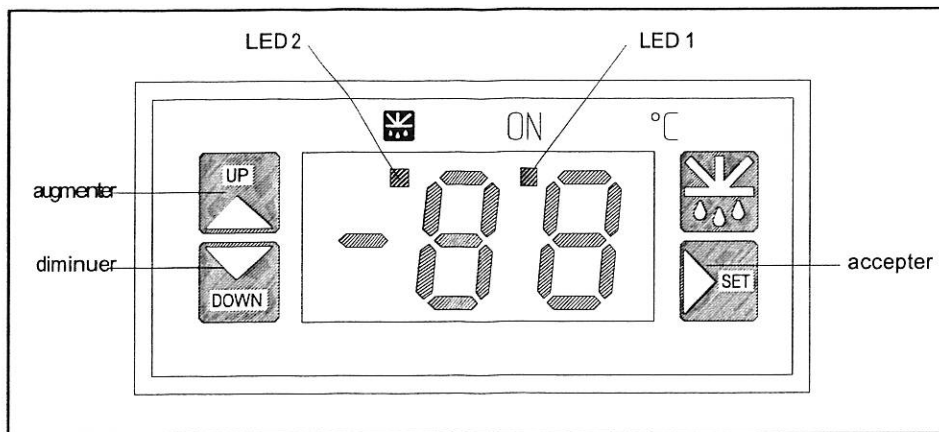
STARTING UP:

Before the introduction of the first charge, it is advisable to leave the chest working for about two hours at the preset temperature.

OPERATION:

1. Connect the plate to the mains.
2. Press the main switch (1).
3. Set point adjust:

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3.1 Push  during 3 seconds until LED 2 starts flashing.

3.2 Select set point by means of keys  

3.3 Accept set point pushing key .

Display will show 3 dashes (---) indicating that the controller is storing new set point.

Les indicateurs / Indicator Leds:

LED 1: Indicateur de relai activé / Indicator of actived relay.

LED 2: Indicateur de désactivation activé / Indicator of actived defrosting.

LED 2 : Intermittent / Intermittent: Phase de programmation / Programming phase.

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3. The stabilization time at the desired temperature will vary according to the ambient temperature.

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MAINTENANCE

Before removing the side covers disconnect the apparatus from the mains.

The guarantee will not have any effect if the apparatus is manipulated by somebody who has not been expressly authorized by J.P.SELECTA.

CLEANING:

Remove the metallic casing and clean the condenser wings with the help of a brusher or compressed air. It is recommended to do it periodically.

GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of J.P. SELECTA,S.A..

Any manipulation of the apparatus by unauthorized personnel by J.P. SELECTA,S.A. cancels the aguarantee automatically.

DECLARATION DE CONFORMITE "CE" "EC" CONFORMITY DECLARATION

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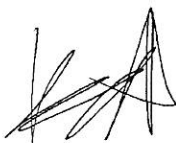
PROGRAMME DE FABRICATION MANUFACTURING PROGRAMME

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Agent en France de J.P. SELECTA S.A