



sartorius

# Tacta®

## Mechanical Pipettes

### Product Information

Tacta® is a premium mechanical pipette featuring superb comfort and reliability. Designed with comfort in mind, Tacta® makes pipetting effortless and safe, while producing accurate and reliable results time after time. Tacta® is made from carefully selected materials, with each component designed to meet the highest standards.



#### Description

Tacta® is a premium mechanical pipette available in volume range from 0.1 to 10 000 µl in single channel models, and from 0.5 µl to 300 µl in multichannel models.

#### Features

- Low pipetting cycle forces reduce the risk of WRULD (Work Related Upper Limb Disorder)
- Ergonomic design for comfortable pipetting
- Optiload feature in both single and multichannel models for easy and light tip loading with perfect tip sealing
- Optijet soft tip ejection feature detaches the tip in a smooth, controlled manner
- Optilock's dual-function volume locking prevents accidental volume changes
- Easy-to-read, four-digit volume display helps to set exact volumes

- Colour-coding of volumes to ease the selection of corresponding pipette tips
- Safe-Cone Filters available for models above 10 µl with convenient built-in filter ejection mechanism
- Fully autoclavable without disassembling
- Simple to clean and maintain with only three parts to disassemble
- Easy calibration adjustment with adjustment scale ensures accurate results, also in cases where factory calibration does not apply, for example when pipetting viscous liquids
- Materials have high chemical and UV-resistance to ensure a long life span for the pipette

#### Intended Use

Manual liquid-handling of volumes ranging from 0.1 µl to 10 ml.



## Technical Data

### Technical Specifications

Weight	79 g (1-ch, 1000 µl) 138 g (8-ch, 300 µl)	Stand options	Carousel Stand for 6 pipettes Linear Stand Pipette Holder for one pipette (included in the pipette package)
Length	225 mm (1-ch, 1000 µl) 240 mm (8-ch, 300 µl)	Safe-Cone Filter ejector	Built-in ejector
Autoclavability	Fully autoclavable (121°C, 20 min, 1 bar/100 kPa)	Limitations	The accuracy and precision values are valid under the controlled conditions described in ISO 8655. Test temperature: +15°C to +30°C +/-0.5°C Relative humidity: >50%
UV-resistance	UV resistance: Tacta <sup>®</sup> pipettes are made of UV-resistant materials. Sartorius pipettes tolerate temporary exposure to UV radiation. Take note that prolonged or frequent exposure to UV radiation may cause yellowing and brittling of the pipette.	Instrument use outside +20°C to +25°C	When using instruments outside this temperature range, the correction factor for water density (Z-value µl/mg) must be used.
Pipetting principle	Air displacement	Transportation conditions	Temperature: -50°C to +50°C Humidity: 10-90% Shocks: <10 G
Pipetting modes	Pipetting (P) Reverse pipetting (rP) Post-delivery mixing	Long-term storage conditions	Temperature: -20°C to +40°C Humidity: 10-60%
		Warranty	2 years

### Volume Adjustment

The calibration of Tacta<sup>®</sup> is based on ISO 8655-6: gravimetric test method for volumetric instruments. It has been factory-checked using the forward-pipetting technique and certified according to ISO 3696 at 22°C using grade-3 distilled water.

The pipette calibration should be adjusted when necessary, for example when reverse pipetting, when pipetting liquids other than water, when using the pipette at high altitudes, or when the geometry of the tip you use clearly differs from standard tip geometry.

For accurate adjustment results, please refer to the table below.

### Volume Adjustment

Model [µl]	Channels	Increment [µl]	-45	-10	-5	-1	0	1	5	10	45
0.1-3	1	0.002	-0.09	-0.02	-0.01	-0.002	0	0.002	0.01	0.02	0.09
0.5-10	1	0.01	-0.45	-0.1	-0.05	-0.01	0	0.01	0.05	0.1	0.45
2-20	1	0.02	-0.9	-0.2	-0.1	-0.02	0	0.02	0.1	0.2	0.9
10-100	1	0.1	-4.5	-1	-0.5	-0.1	0	0.1	0.5	1	4.5
20-200	1	0.2	-9	-2	-1	-0.2	0	0.2	1	2	9
100-1000	1	1	-45	-10	-5	-1	0	1	5	10	45
500-5000	1	5	-225	-50	-25	-5	0	5	25	50	225
1000-10000	1	10	-450	-100	-50	-10	0	10	50	100	450
0.5-10	8 & 12	0.01	-0.45	-0.1	-0.05	-0.01	0	0.01	0.05	0.1	0.45
5-100	8 & 12	0.1	-4.5	-1	-0.5	-0.1	0	0.1	0.5	1	4.5
30-300	8 & 12	0.2	-9	-2	-1	-0.2	0	0.2	1	2	9

## Ordering Information

Tacta® Order Code	Channels	Volume Range (µl)	Increment (µl)	Test Volume (µl)	Systematic Error <sup>N</sup> Limit ±		Random Error <sup>N</sup> Limit	
					(%)	(µl)	(%)	(µl)
LH-729010	1	0.1–3	0.002	3	1.4	0.042	0.8	0.024
				1.5	2.6	0.039	1.6	0.024
				0.3	10.0	0.030	6.0	0.018
LH-729020	1	0.5–10	0.01	10	1.0	0.100	0.6	0.060
				5	1.5	0.075	1.0	0.050
				1	3.0	0.030	2.0	0.020
LH-729030	1	2–20	0.02	20	1.0	0.200	0.5	0.100
				10	1.4	0.140	0.9	0.090
				2	4.0	0.080	3.0	0.060
LH-729050	1	10–100	0.10	100	0.8	0.80	0.2	0.20
				50	1.0	0.50	0.4	0.20
				10	3.0	0.30	1.0	0.10
LH-729060	1	20–200	0.20	200	0.6	1.20	0.2	0.40
				100	0.8	0.80	0.3	0.30
				20	2.3	0.46	0.9	0.18
LH-729070	1	100–1,000	1.00	1,000	0.7	7.0	0.2	2.0
				500	0.8	4.0	0.2	1.0
				100	2.5	2.5	0.6	0.6
LH-729080	1	500–5,000	10.0	5,000	0.6	30	0.2	10
				2,500	0.7	17.5	0.3	7.5
				500	2.4	12	0.6	3
LH-729090	1	1,000–10,000	20.0	1,000	0.6	60	0.2	20
				5,000	1.2	60	0.3	15
				1,000	3.0	30	0.6	6
LH-729120	8	0.5–10	0.01	10	1.5	0.150	1.0	0.100
LH-729220	12			5	2.5	0.125	2.0	0.100
				1	5.5	0.055	4.0	0.040
LH-729130	8	5–100	0.10	100	0.9	0.90	0.4	0.40
LH-729230	12			50	1.2	0.60	0.7	0.35
				10	4.0	0.40	2.0	0.20
LH-729140	8	30–300	0.20	300	0.6	1.80	0.25	0.75
LH-729240	12			150	1.0	1.50	0.5	0.75
				30	2.5	0.75	1.0	0.30

<sup>N</sup> Note: The listed systematic and random error values can be achieved only under strictly controlled conditions during type tests per ISO 8655. Due to the continuous product development by Sartorius, the systematic and random error values are subject to change without prior notice.

# Sales and Service Contacts

For further contacts, visit [www.sartorius.com](http://www.sartorius.com)

## Europe

**Germany**  
Sartorius Lab Instruments  
GmbH & Co. KG  
Otto-Brenner-Strasse 20  
37079 Goettingen  
Phone +49.551.308.0  
Fax +49.551.308.3289

**France & Suisse Romande**  
Sartorius France  
2, rue Antoine Laurent de Lavoisier  
ZA de la Gaudriée  
91410 Dourdan  
Phone +33.1.70.62.50.00  
Fax +33.1.64.59.76.39

**Austria**  
Sartorius Austria GmbH  
Modecenterstrasse 22  
1030 Vienna  
Phone +43.1.7965760.0  
Fax +43.1.7965760.24

**Belgium**  
Sartorius Belgium N.V.  
Rue Colonel Bourg 105  
1030 Bruxelles  
Phone +32.2.756.06.90  
Fax +32.2.481.84.11

**Finland & Baltics**  
Sartorius Biohit Liquid Handling Oy  
Laippatie 1  
00880 Helsinki  
Phone +358.9.755.951  
Fax +358.9.755.95.200

**Hungary**  
Sartorius Hungária Kft.  
Kagyló u. 5.  
2092 Budakeszi  
Phone +3623.457.227  
Fax +3623.457.147

**Ireland**  
Sartorius Ireland Ltd.  
Unit 41, The Business Centre  
Stadium Business Park  
Ballycoolin Road  
Dublin 11  
Phone +353.1.8089050  
Fax +353.1.8089388

**Italy**  
Sartorius Italy S.r.l.  
Via Torino 3/5  
20814 Varedo (MB)  
Phone +39.0362.5557.11  
Fax +39.0362.5557.99

**Netherlands**  
Sartorius Netherlands B.V.  
Phone +31.30.60.53.001  
Fax +31.30.60.52.917  
[info.netherlands@sartorius.com](mailto:info.netherlands@sartorius.com)

**Poland**  
Sartorius Poland sp.z o.o.  
ul. Wrzesinska 70  
62-025 Kostrzyn  
Phone +48.61.6473830  
Fax +48.61.6473839

**Russian Federation**  
LLC "Sartorius RUS"  
Vasilyevsky Island  
5<sup>th</sup> line 70, Lit. A  
199178 St. Petersburg  
Phone +7.812.327.53.27  
Fax +7.812.327.53.23

**Spain & Portugal**  
Sartorius Spain, S.A.  
Avda. de la Industria, 32  
Edificio PAYMA  
28108 Alcobendas (Madrid)  
Phone Spain +34.913.586.095  
Phone Portugal +351.800.855.800  
Fax Spain +34.913.589.623  
Fax Portugal +351.800.855.799

**Switzerland**  
Sartorius Mechatronics Switzerland AG  
Ringstrasse 24a  
8317 Tagelswangen (ZH)  
Phone +41.44.746.50.00  
Fax +41.44.746.50.50

**U.K.**  
Sartorius UK Ltd.  
Longmead Business Centre  
Blenheim Road, Epsom  
Surrey KT19 9QQ  
Phone +44.1372.737159  
Fax +44.1372.726171

**Ukraine**  
LLS "Sartorius RUS"  
Post Box 440 "B"  
01001 Kiev, Ukraine  
Phone +380.44.411.4918  
Fax +380.50.623.3162

## Americas

**USA**  
Sartorius Corporation  
5 Orville Drive, Suite 200  
Bohemia, NY 11716  
Phone +1.631.254.4249  
Toll-free +1.800.635.2906  
Fax +1.631.254.4253

**Argentina**  
Sartorius Argentina S.A.  
Int. A. Ávalos 4251  
B1605ECS Munro  
Buenos Aires  
Phone +54.11.4721.0505  
Fax +54.11.4762.2333

**Brazil**  
Sartorius do Brasil Ltda  
Avenida Senador Vergueiro 2962  
São Bernardo do Campo  
CEP 09600-000 - SP- Brasil  
Phone +55.11.4362.8900  
Fax +55.11.4362.8901

**Canada**  
Sartorius Canada Inc  
1173 North Service Road West, D4  
Oakville, ON L6M 2V9  
Phone +1.905.569.7977  
Toll-Free +1.800.668.4234  
Fax +1.905.569.7021

**Mexico**  
Sartorius de México, S.A. de C.V.  
Libramiento Norte de Tepotzotlán s/n,  
Colonia Barrio Ilacateco,  
Municipio de Tepotzotlán,  
Estado de México,  
C.P. 54605  
Phone +52.55.5562.1102  
Fax +52.55.5562.2942  
[leadsmex@sartorius.com](mailto:leadsmex@sartorius.com)

**Peru**  
Sartorius Peru S.A.C.  
Avenue Alberto del Campo 411  
Floor 12 - The Office  
15076 - San Isidro, Lima  
Phone +51.1.441 0158  
Fax +51.1.422 6100

## Asia | Pacific

**Australia**  
Sartorius Australia Pty. Ltd.  
Unit 5, 7-11 Rodeo Drive  
Dandenong South Vic 3175  
Phone +61.3.8762.1800  
Fax +61.3.8762.1828

**China**  
Sartorius (Shanghai) Trading Co., Ltd.  
3rd Floor, North Wing, Tower 1  
No. 4560 Jinke Road  
Zhangjiang Hi-Tech Park  
Pudong District  
Shanghai 201210, P.R. China  
Phone +86.21.6878.2300  
Fax +86.21.6878.2882

**Hong Kong**  
Sartorius Hong Kong Ltd.  
Unit 1012, Lu Plaza  
2 Wing Yip Street  
Kwun Tong  
Kowloon, Hong Kong  
Phone +852.2774.2678  
Fax +852.2766.3526

**India**  
Sartorius Weighing India Pvt. Ltd.  
#69/2-69/3, NH 48, Jakkasandra,  
Nelamangala Tq  
562 123 Bangalore, India  
Phone +91.80.4350.5250  
Fax +91.80.4350.5253

**Japan**  
Sartorius Japan K.K.  
4th Fl, Daiwa Shinagawa North Bldg.  
8-11, Kita-Shinagawa 1-chome  
Shinagawa-ku, Tokyo, 140-0001 Japan  
Phone +81.3.3740.5408  
Fax +81.3.3740.5406

**Malaysia**  
Sartorius Malaysia Sdn. Bhd  
Lot L3-E-3B, Enterprise 4  
Technology Park Malaysia  
Bukit Jalil  
57000 Kuala Lumpur, Malaysia  
Phone +60.3.8996.0622  
Fax +60.3.8996.0755

**Singapore**  
Sartorius Singapore Pte. Ltd  
10 Science Park Rd  
The Alpha #02-13/14  
Singapore Science Park II  
Singapore 117684  
Phone +65.6872.3966  
Fax +65.6778.2494

**South Korea**  
Sartorius Korea Ltd.  
8th Floor, Solid Space B/D,  
PanGyoYeok-Ro 220, Bundang-Gu  
SeongNam-Si, GyeongGi-Do, 463-400  
Phone +82.31.622.5700  
Fax +82.31.622.5799

**Thailand**  
Sartorius (Thailand) Co. Ltd.  
129 Rama 9 Road,  
Huaykwang  
Bangkok 10310  
Phone +66.2643.8361-6  
Fax +66.2643.8367



◀ [www.sartorius.com](http://www.sartorius.com)