





USE OF THE BIOHAZARD SPILL KIT

FOR BLOOD AND BLOOD-STAINED BODY FLUID SPILLS

BEFORE YOU START

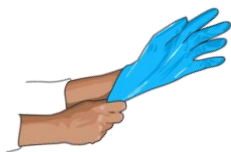

Always wear gloves and an apron. Use eye protection if required.

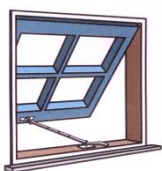

Wherever possible ensure good ventilation when using all chlorine products.




Be aware of the safety notes on the product labels and always check the expiry date.




Sprinkle Haz-Tab Granules over spill until all moisture is completely absorbed.

Leave for no more than two minutes.

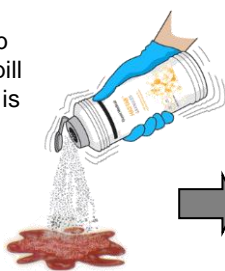

Peel backing away from adhesive strip on yellow disposal bag and attach bag to wall or table and allow to hang open.

Retain backing to act as a tie later.

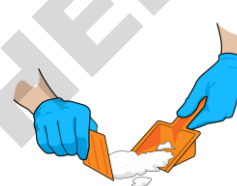

Collect spilt material and granules mixture using scoop and scraper and discard all into the yellow disposal bag.

When you need to make up a chlorine solution for splashes or drips on vertical surfaces, walls, etc.



Fill the 1 litre diluter to the line with cold water then add 4 Haz-Tab Tablets.

Set aside for 2 minutes to allow the tablets to dissolve.



When tablets have dissolved screw down the diluter cap and mix gently by inversion.

DO NOT SHAKE THE DILUTER



Use the solution on disposable towels to wipe splashes on walls and remove any smears.




Discard all remaining solutions by flushing down the sink or sluice

REPLACE THE DILUTER IN THE YELLOW BOX



Place all used materials, gloves last, into the disposal bag. Secure with the tie.

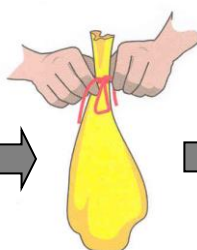


The bag and its contents should be destroyed unopened by incineration.

DON'T FORGET TO WASH YOUR HANDS!



Important Points:

- **Never** use on chemical spills or mix with any other chemicals, including cleaning products.
- **Never** use hot or warm water to dissolve the tablets.
- **Never** use the solution in a spray bottle.
- Chlorine products may bleach soft furnishings and corrode metal (not stainless steel).
- Refer to the Infection Control Policy for full details on the use of these products.
- If in doubt contact your Infection Control Nurse.

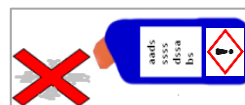

Issue Date: March 2021

* **HAZ-TAB** products are manufactured by **Guest Medical Limited** of Aylesford, Kent. **01622 791895**