



661 Route 23 South
Wayne, New Jersey 07470
Tel: 973-694-0500 * Fax: 973-694-7199
Internet Address: <http://www.bel-art.com>

Velveteen Squares Cleaning

Although Velveteen Squares can be washed / cleaned, it is recommended to use new sterile squares from the sterile package. Remove one square at a time with either a gloved hand or forceps to prevent contamination, close the package tightly immediately after removing the Velveteen Square.

If you choose to clean the squares for reuse, hand wash the Velveteen Squares with a mild detergent, rinse thoroughly with warm clean water. Squares can then be stacked on Kraft paper or aluminum foil and autoclaved. The Velveteen Squares may shrink during drying cycle of the autoclave. Not all the squares may become dry during autoclave drying cycle and should be allowed to dry prior to use.





661 Route 23 South
Wayne, New Jersey 07470
Tel: 973-694-0500 * Fax: 973-694-7199
Internet Address: <http://www.bel-art.com>

Dominique Dutscher SAS