Instrucciones

INSTALACIÓN

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Instructions

Four Channel Timer with White Board

Installation

1. Open the battery compartment by pushing tab back
2. Install 2 AAA batteries following the correct polarity
3. Close the battery compartment by snapping the battery cover on

Operation

Count Down Function

- Press the [EVENT] button to move from timer 1 (T1) to timer 2 (T2) to timer 3 (T3) to timer 4 (T4)
- Press the [HR] button to set the hours to count down
- Press the [MIN] button to set minutes to count down
- Press the [START/STOP] to start the timer counting down
- The alarm will go off when the display counts down to 00:00
- The alarm will continue to sound for five minutes as the time elapsed is counted up
- Press the [START/STOP] to stop counting - the remaining time will be displayed
- Press the [START/STOP] to continuing the counting
- Press the [HR] and [MIN] buttons simultaneously to reset the display value to "00:00"
- Press the [START/STOP] button to start counting up
- When display counts up to maximum range the timer will stop

Count Up Function

- Press the [EVENT] button to move from timer 1 (T1) to timer 2 (T2) to timer 3 (T3) to timer 4 (T4)
- Press the [START/STOP] to start the timer counting up
- Press the [START/STOP] to stop counting - the elapsed time will be displayed
- Press the [START/STOP] to continuing the counting
- When display counts up to maximum range the timer will stop
- Press the [HR] and [MIN] buttons simultaneously to reset the display value to "00:00"

Battery Handling/Packaging

- Remove batteries from equipment immediately after they fail to operate the equipment.
- If equipment uses two or more batteries, always replace batteries in complete sets.
- Do not attempt to recharge alkaline batteries, they are not rechargeable.
- Do not handle hot or warm batteries.
- Do not heat, incinerate, crush, puncture or mutilate batteries.
- Wear personal protective equipment if batteries show signs of leakage, bulging, swelling or deformity.

Warning: If potassium hydroxide electrolyte from an alkaline battery comes in contact with the skin, do not try to neutralize the electrolyte with vinegar or any other acidic solutions. Neutralization may trap electrolyte on the skin. Flush the affected skin area with copious amounts of water. If the battery electrolyte gets into your eyes, it can cause severe damage and/or blindness.

To register this product for a one year limited warranty, visit www.belart.com/warranty

H-B Instrument – SP Scienceware
Triple Accredited/Registered
ISO 9001:2015 Registered Company
ISO 14001:2004 Registered Company
ISO/IEC 17025:2017 Accredited Laboratory

For more information call (973) 694-0500 • Fax (973) 694-7199
info@belart.com • www.belart.com/hb

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