



PREPSAFE™ MINI FLOATING MCT RACK

Patent Pending

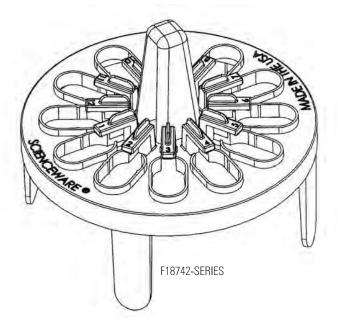

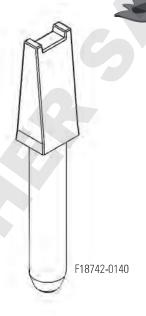

INSERTING MICROCENTRIFUGE TUBES:

- 1. Hold rack securely in one hand.
- 2. Insert a microcentrifuge tube into the desired location.
- 3. Hold the tube underneath rack and slide it completely under the immobilizing tab until the tube snaps in place securing it in the rack.

REMOVING MICROCENTRIFUGE TUBES:

- 1. Hold rack securely in one hand.
- 2. Place your finger underneath the rack behind the tube you wish to remove and slide it out from under the immobilizing tab.

INSERTING VORTEXING ATTACHMENT:

- 1. Holding the rack in one hand, invert rack so that the legs are pointing upward.
- 2. Insert the larger end of the vortexing attachment into the hollow space within the handle of the rack until fully engaged.

REMOVING VORTEXING ATTACHMENT:

- 1. Holding the rack in one hand, invert rack so that the legs are pointing upward.
- 2. Pull vortexing attachment away from rack.

| CAIALUG NO. | DESCRIPTION |
|-------------|--|
| F18742-0120 | Clear PrepSafe™ Mini MCT Rack without Vortexing Attachment |
| F18742-0121 | Turquoise PrepSafe™ Mini MCT Rack without Vortexing Attachment |
| F18742-0122 | Fuchsia PrepSafe™ Mini MCT Rack without Vortexing Attachment |
| F18742-0123 | Neon Yellow PrepSafe™ Mini MCT Rack without Vortexing Attachment |
| F18742-0124 | Clear PrepSafe™ Mini MCT Rack with Vortexing Attachment |
| F18742-0125 | Turquoise PrepSafe™ Mini MCT Rack with Vortexing Attachment |
| F18742-0126 | Fuchsia PrepSafe™ Mini MCT Rack with Vortexing Attachment |
| F18742-0127 | Neon Yellow PrepSafe™ Mini MCT Rack with Vortexing Attachment |
| F18742-0140 | Replacement Vortexing Attachment |

Ecienceware Bel-Art Products



TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

918742028

