

NEW 10 ML DURAN® LABORATORY GLASS BOTTLE

The original DURAN® laboratory glass bottles are ideal for a huge range of uses in the laboratory thanks to their incredible properties, which include a very high chemical resistance, nearly inert behaviour and a high usage temperature.

The trend for **miniaturisation** allows for a decrease in the quantities of material or fluid required for testing in precise analytical methods. If less material is used, procurement costs are reduced and the amount of waste is lowered.



Areas of application

- Storing expensive/sensitive media, reference samples etc.
- Packaging bottle for reference samples, stability tests, round robin tests and deuteration solutions
- Shipping bottle for small sample volumes
- Storing proteins, vitamins, DNA

Target groups

- Pharmaceutical industry
- Bioengineering industry
- Food and beverage industry



Best cleaning method

The best way to clean and rinse DURAN® bottles is in a cleaning machine with injector nozzles. Insert the bottles with the opening facing downwards. To prevent damage during cleaning, ensure that the bottles are not placed too close together. The injector nozzle should not come into contact with the base of the bottle.

The cap is rinsed and cleaned using a rotating spray arm. Place and cover the caps in a layer with the opening facing downwards on the tray with perforated plate or mesh.

Product features

- Glass type I/Neutral glass in accordance with USP⁽¹⁾/EP⁽²⁾ and JP⁽³⁾
- Tried-and-tested DURAN® properties
- The threaded Polypropylene cap with plug seal offers a reliable sealing system
- Popular thread type GL 25: existing caps can be used
- Single-handed opening allows easy pipetting
- Reduced ullage - minimises air available to react with media
- Autoclavable/sterilisable
- Also available in amber or with a plastic coating

¹United States Pharmacopoeia / ²European Pharmacopoeia / ³Japanese Pharmacopoeia

Order information

Order no.	Description	Capacity (ml)	PU/quantity
21 801 08	DURAN® GL 25 laboratory glass bottle, clear, without cap	10	10
21 801 08 5	DURAN® GL 25 laboratory glass bottle, clear, with cap	10	10
21 806 08	DURAN® GL 25 laboratory glass bottle, amber, without cap	10	10
21 806 08 5	DURAN® GL 25 laboratory glass bottle, amber, with cap	10	10
21 805 08	DURAN® protect GL 25 laboratory glass bottle, without cap	10	10

Accessories	Description	Thread	PU/quantity
11 296 00	Premium cap (TpCh260)	25	5
89 299 13 04	Premium cap STERICLIN®-packed	25	20
29 240 13	Screw cap (PBT)	25	10
29 239 13	Screw cap (PP)	25	10
11 832 50	Membrane screw cap (PP)	25	5

These products are available from your specialist laboratory dealer.



Premium cap



PBT screw cap



PP screw cap



Membrane screw cap