

Notice d'emploi – Réf 0706CA – Compteur / Décompteur 100 mn / Façade métallique

- Capacité : 100 mn à la seconde
- Précision : +/- 1,7 sec /24heures
- Alarme : 1 mn / Mémoire
- Écran : 47x20 mm / Ch : 17 mm
- Façade alu brossé
- Béquille / Aimant / Fixation murale
- D : 72x65x19 mm / 70 g
- Pile : AAA (isolée)

FONCTION COMPTEUR

En mode Comptage (affichage 00 :00), appuyer sur START pour commencer le comptage.
Appuyer sur STOP pour arrêter le comptage.
Appuyer simultanément sur MIN + SEC pour remettre à zéro le compteur.

REGLAGE DU DECOMPTEUR

Lorsque le compteur affiche 00 :00, rester appuyé sur les touches MIN ou SEC pour incrémenter et programmer le temps souhaité.
Appuyer sur START pour débiter le décompte.
Appuyer sur STOP pour arrêter le décompte.
A échéance, un signal sonore de 1 min retenti. Le temps programmé reste en mémoire.

CHANGEMENT DE LA PILE

Lorsque l'affichage ou l'alarme devient faible, il faut changer la pile.
Ouvrir la trappe arrière et insérer une pile neuve LR03/AA en veillant à respecter la polarité.

Note : Veiller à déposer le matériel et les piles usagés dans un endroit écologiquement approprié

Operating instructions - Ref 0706CA - 100 mn count up/down / Metal front panel

- Capacity : 100 min with 1 second accuracy
- Precision : +/- 1.7 sec /24 hours
- Alarm : 1 mn / Memory
- Display : 47x20 mm / Dgt : 17 mm
- Brushed aluminium front panel
- Lever handle / Magnet / Wall mounting
- D : 72x65x19 mm / 70 g
- Battery : AAA (insulated)

COUNT UP FUNCTION

In count up mode (display 00 :00), press START to start counting.
Press STOP to stop counting.
Press together MIN + SEC to reset the counter.

SETTING COUNT DOWN TIMER

When the counter displays 00 :00, press and maintain MIN or SEC key to increase and set the desired time.
Press START to start counting down.
Press STOP to stop the countdown.
At the end of the countdown, a 1 min signal sounds. The programmed time remains in memory.

CHANGING THE BATTERY

When the display or alarm becomes weak, change the battery.
Open the rear cover and insert a new LR03/AA battery, making sure that the polarity is correct.

Note: Take care to dispose of used equipment and batteries in an environmentally appropriate place.