
Ref 098494

Notice d'emploi / Instruction / Anweisung / Instrucción / Instrução / Instructie

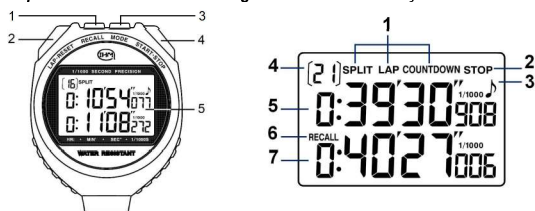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1 - DESCRIPTION

Features:

- 2 Display lines
- Ability 10h to 1 / 1000th sec
- 99 SPLIT time (elapsed from the start) or LAP (between intervals)
- 30 SPLIT or LAP memories
- Repetitive Countdown
- Buttons Colors: yellow, red, black
- Bip On / Off
- Automatic power off
- Batterie : LR44

Description Stopwatch :

Screen display :

- | | |
|----------------------|--|
| 1. RECALL Button | 1. Icon Split / Lap / Countdown |
| 2. Button LAP/RESET | 2. Stop Icon |
| 3. MODE Button | 3. Beep Icon |
| 4. START/STOP button | 4. No. Split / Lap / repeat Number |
| 5. LCD Display | 5. Time Split / LAP / programmed Countdown |
| | 6. Recall Icon |
| | 7. Current Time |

2 - OPERATING

Select the Stopwatch Mode SPLIT or LAP countdown by pressing the MODE button.

a) Stopwatch SPLIT or LAP

- Press START/STOP to begin timing.
- Press LAP/RESET to store a time. The number of SPLIT/LAP increments on the top left of the screen.
- Press START/STOP to stop timing.
- Press LAP/RESET to reset the stopwatch.

b) Time recall SPLIT/LAP

- In Timing mode, current time or stopped, press RECALL to enter recall memories.
- Press RECALL again to cycle the fastest time (FS LAP), slowest (SL LAP) and average time (AV LAP).
- Press LAP/RESET or START/STOP to display the top / bottom, different saved times.
- Press MODE to return to clock mode.

Note: The stopwatch can measure up to 99 times SPLIT/LAP but can not remember 30.

c) Countdown

- Press And hold RECALL to begin setting. Hours flash.
- Press START/STOP to change value of fSPLIT / LAP flashing digits.
- Press LAP/RESET to set the other values in this order: Hours → Minutes → Seconds
- Press MODE to confirm.

- Press START/STOP to start or stop the countdown. The number of repeat increments
- Press And hold RECALL to put the time to zero count.
- Press LAP/RESET to set the countdown and the number of repetitions to zero

Note: If programmed time exceeds 5 seconds, a beep sounds and the last 5 seconds due.

Note: The countdown is repeated 99 times. In addition, the number of repeat again from 0.

3 - SETTING THE TONE / BEEP

- Press And hold MODE and START/STOP for 2 seconds to enable / disable the tone. The symbol ♪ appears on the screen.

4 - POWER SAVE / BATTERY REPLACEMENT

- In Order to ensure long battery life, the LCD turns off automatically after 30 minutes without use, when stationary. Press any button to turn the display.
- This Timer is equipped with 2 batteries in parallel. 1 single cell is in operation, at a time. If the screen or the sound becomes weak, change the / stack (s).
- Unscrew housing base, insert a new LR44 battery and screw the bottom case making sure to replace the gasket.

Note: ensuring waste batteries and deposit the materials in an environmentally appropriate place.

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