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DOMINIQUE DUTSCHER SAS

Contents

| | |
|---|----|
| 1. Preface | 4 |
| 2. Introduction | 5 |
| 3. Safety precautions | 6 |
| 4. EEC - Declaration of Conformity | 7 |
| 5. Unpacking and checking the equipment supplied | 8 |
| 6. Description of the MD8 airscan Remote-Control | 9 |
| 7. Specifications of the MD8 airscan Remote-Control | 10 |
| 8. Switching the Remote-Control on | 11 |
| 9. How the Remote-Control functions | 12 |
| 10. Terminal assignment | 13 |
| 11. Warranty | 14 |
| 12. Information and instruction on disposal | 15 |
| Non-hazardous Declaration | 17 |

Version 1.0 - as of: June 2009

Please read these operating instructions carefully before switching the instrument on!

1. Preface

Dear new user

Congratulations!

With your decision to purchase the MD8 airscan Remote Control 1ZE--0003, a newly developed accessory for our practice-proven microbiological air sampler, you have selected a product which is manufactured according to the highest quality German electrical and precision engineering standards.

We are sure that it will provide you with satisfying and trouble-free service for many years.

Please do not hesitate to contact us, should you need further information or assistance on installation or on your specific applications.

Goettingen, 2009

2. Introduction

The MD8 airscan Remote Control 1ZE---0003 is an interface which allows the user to carry out automatic air sampling with the MD8 airscan air sampler at several collection locations in succession in a process line via a PLC.

These operating instructions are mainly intended to provide a technical description on how the Remote Control is to be connected to and operated with an available process logic controller (PLC).

Therefore, a complete function test of the Remote Control is only possible in conjunction with a PLC and a connected MD8 airscan. These Operating Instructions have been primarily written for technicians.

As these air monitoring systems are customer-specific solutions, further details concerning installation of this equipment must be discussed on site with the user and the equipment needs to be installed together with the user.

DOMINIQUE DUTSCHER SAS

3. Safety precautions

1. Switch off the instrument before wiring the MD8 airscan Remote-Control 1ZE---0003.
2. Disconnect the instrument from the electrical supply (unplug it) before changing the fine fuse.
3. Only switch on the MD8 airscan Remote Control when it is connected to an MD8 airscan air sampling instrument and a PLC unit.
4. Any use other than for the intended purpose, or improper operation, voids any and all manufacturer warranties.
5. Before starting the system, make sure that the available electrical supply matches that which is printed on the label on the bottom of the instrument.
6. Do not open or make changes to the instrument. Damage resulting from such actions or changes are not covered by the manufacturer's warranty.

4. EEC – Declaration of Conformity

We hereby declare that this version of the MD8 airscan Remote Control (order number 1ZE---0003) conforms, in its design and function, to the requirements of the following standards:

EN 50081-1
EN 55022 Cl. B
EN 50082-2
EN 61000-4-2, -3, -4, -5, -6,-11
EN 50204

As of: Oct. 2000

This declaration becomes invalid if unauthorized changes have been made to the instrument.

DOMINIQUE DUTSCHER SAS

5. Unpacking and checking the equipment supplied

Sartorius Stedim Biotech products are packaged very carefully. However, despite this, damage may still occur during transport, so please check the equipment supplied for any damage while unpacking it, and at the same time check each piece against the following list, to ensure that everything has been correctly delivered and that nothing is left in the packaging:

- MD 8 airscan Remote Control (check the voltage)
- a) Line cord (country specific) (see Fig.1)
- b) Connecting cable to the MD8 airscan
- c) Connecting cable to the PLC
- Operating instructions

Fig. 1

c

b

a

6. Description of the MD8 airscan Remote Control

The MD8 airscan Remote Control has a metal housing which can be wiped down and disinfected using common cleaning agents and disinfectants.

The operating elements (on | off switch (a) and status indicators (b)) for the various instrument functions, which are described in the section on how the Remote Control functions, are located on the front of the instrument (Fig. 2).

The power supply connection (a), fuse (0.5 A T, (b)) and connections for the PLC (c) and for an MD8 airscan air sampler (d) are on the back of the instrument (Fig.3).

Fig. 2

Fig 3

7. MD8 airscan Remote Control specifications

| | |
|---------------------------|-------------------------------|
| Order number | 1ZE---0003 |
| Supply voltage | 90 – 260 V ~ 50 60 Hz |
| Power consumption | 5 W |
| Safety fuse | 500 mA slow |
| Connecting voltage output | 24 V = 0.3 Amp |
| Weight | Approx. 2.2 kg |
| Dimensions | 258 × 159 × 94 mm (L × W × H) |
| Total width | 215 mm with connecting plug |

8. Switching on the Remote Control

Use the on | off switch on the front of the instrument to switch the MD8 airscan Remote Control on (Fig. 2 (a) + Fig. 4 (a)). The switch lights up red to show that the instrument is on, and at the same time, the instrument carries out a self test in which it checks its functions. The status indicator LED flashes during this test.

When the test is successfully completed, and provided the instrument is connected to an MD8 airscan air sampler, the green LED stays on ("on") (Fig. 4, b). This shows that the instrument is ready for use and the measurement (air sampling) can be started.

Should the test not be successfully completed, the green LED ("on") is not lit. Possible cause of the fault: Remote Control and MD8 airscan are not connected.

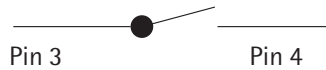
9. How the Remote Control functions

A measurement can be started when the green LED ("on") of the MD8 airscan Remote Control is on.

To start the measurement procedure, the PLC control must trigger a relay with a closing contact for about 100 ms.

The closing contact of the relay must be connected at pins 3 and 4, potential free.

As soon as the MD8 airscan air sampler and the MD8 airscan Remote Control are supplied with a 220 V operating voltage and switched on, the measurement procedure can be started. At the end of a measurement, provided no fault has occurred, a signal is output at pin 5 (signal: Ready = yellow LED (Fig. 4, c)).

After the measurement procedure has been initiated with the start pulse, a 24 V return signal is output at pin 2 (during the complete measuring phase). After the measuring phase, the instrument must be reset to start again by an acknowledgement signal at pin 1 (signal: Running = yellow LED, (Fig. 4, d))

If the measurement was not successfully carried out, a fault is reported by a flashing signal which alternates between pin 2 and pin 5, and by a 24 V return signal at pin 6. In this case, the instrument can be reset to start again by a double acknowledgement at pin 1 (signal: Error = red LED, (Fig. 4,e)).

Pin 1 (Enter | Acknowledge) must be externally switched to the MINUS of the supply voltage.

Pin 7 is connected to the external supply voltage for the galvanic separated outputs + 24 V =.

The PLC connecting plug

The signal plug is a midget round plug connector made by Binder.

Order number: 09 - 0327 - 00 - 07

Flange plug, solderable, can be screwed from the back

Number of pins: 7

10. Terminal assignment


Socket pin assignment | Colors of the connecting wires

| Pin no. | Color | Function | Type |
|---------|--------|--|--------|
| 1 | Blue | Enter Acknowledge | Input |
| 2 | Yellow | Message Measurement running High Signal = 24 V | Output |
| 3 | White | Start Stop Signal Measurement Potential free | Input |
| 4 | Green | Start Stop Signal Measurement Potential free | Input |
| 5 | Brown | Message Measurement finished High Signal = 24 V | Output |
| 6 | Gray | Message Error High Signal = 24 V | Output |
| 7 | Pink | 24 V Supply voltage external | |

Shielding: brown, thick

7 Pin socket External

| Connection | Cable | Note |
|------------|--------|----------------------|
| 1 | Blue | Enter Acknowledge |
| 2 | Yellow | Measurement running |
| 3 | White | Start Stop |
| 4 | Green | Start Stop |
| 5 | Brown | Measurement finished |
| 6 | Gray | Error |
| 7 | Pink | + 24V DC external |

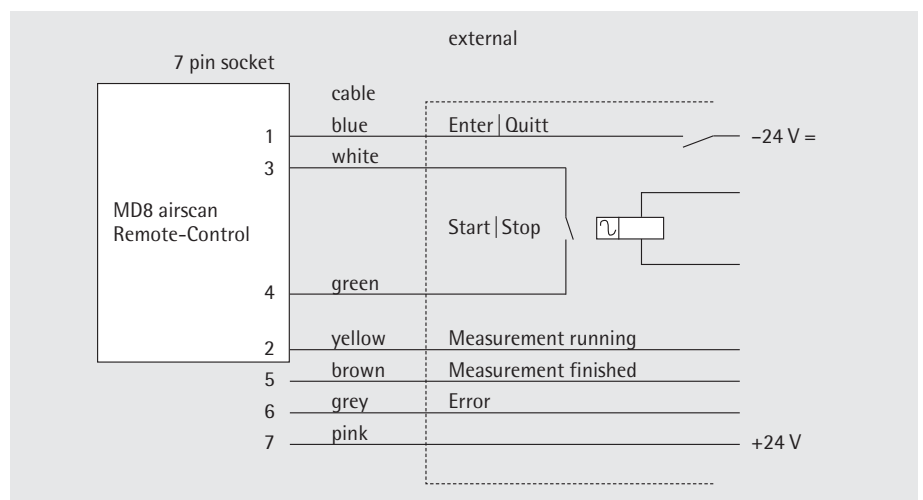Specifications subject to change without notice.

11. Warranty

The warranty conditions can differ somewhat in various countries. Should no warranty card be included with the instrument, please contact the Sartorius Stedim Biotech subsidiary in your country at the address given on the back of this publication.

In Germany, the Remote Control has a factory warranty of 2 years from the date of purchase.

DOMINIQUE DUTSCHER SAS

12. Information and instructions on disposal

Attention must be paid to the applicable laws, rules and regulations, when disposing of the contaminated components.

The packaging is to be taken to a local waste disposal site if no longer required. The packaging is made of environmentally friendly materials that can be used as secondary raw materials.

The equipment, including accessories and batteries, does not belong in your regular household waste. EU legislation requires its Member States to collect electrical and electronic equipment and disposed of it separately from other unsorted municipal waste with the aim of recycling it.

In Germany and several other countries, Sartorius Stedim Biotech GmbH itself assumes responsibility for the return and conformant disposal of its electronic and electrical products. These products may not be placed with household waste or brought to collection centers run by local public disposal operations – not even by small commercial operators.

For disposal in Germany and in the other member nations of the European Economic Area (EEA), please contact our local service technicians or our Service Center in Goettingen, Germany:

Sartorius Stedim Biotech GmbH
Servicezentrum Biotechnologie
August-Spindler-Strasse 11
37079 Goettingen, Germany

WEEE-Reg.-No. DE 89907997

In countries that are not members of the European Economic Area (EEA) or where no Sartorius Stedim Biotech subsidiaries or dealerships are located, please contact your local authorities or a commercial disposal operator.

Prior to disposal and |or scrapping of the equipment, any batteries should be removed and disposed of in local collection boxes.

Sartorius Stedim Biotech will not take back equipment contaminated with hazardous materials (ABC contamination) – either for repair or disposal. Please refer to the package leaflet | the included CD ROM or to our web site (www.sartorius-stedim.com) for more detailed information and addresses for repair service or disposal of your device.

Sales and Service Contacts

For further contacts, visit www.sartorius-stedim.com

Europe

Germany
Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen
Phone +49.551.308.0
Fax +49.551.308.3289
www.sartorius-stedim.com

Sartorius Stedim Systems GmbH
Schwarzenberger Weg 73-79
34212 Melsungen
Phone +49.5661.71.3400
Fax +49.5661.71.3702
www.sartorius-stedim.com

France
Sartorius Stedim Biotech S.A.
ZI Les Paluds
Avenue de Jouques – BP 1051
13781 Aubagne Cedex
Phone +33.442.845600
Fax +33.442.845619

Sartorius Stedim France SAS
ZI Les Paluds
Avenue de Jouques – CS 71058
13781 Aubagne Cedex
Phone +33.442.845600
Fax +33.442.846545

Austria
Sartorius Stedim Austria GmbH
Franzosengraben 12
A-1030 Vienna
Phone +43.1.7965763.18
Fax +43.1.796576344

Belgium
Sartorius Stedim Belgium N.V.
Leuvensesteenweg, 248/B
1800 Vilvoorde
Phone +32.2.756.06.80
Fax +32.2.756.06.81

Denmark
Sartorius Stedim Nordic A/S
Hoerskaetten 6D, 1.
DK-2630 Taastrup
Phone +45.7023.4400
Fax +45.4630.4030

Italy
Sartorius Stedim Italy S.p.A.
Via dell'Antella, 76/A
50012 Antella-Bagno a Ripoli (FI)
Phone +39.055.63.40.41
Fax +39.055.63.40.526

Netherlands
Sartorius Stedim Netherlands B.V.
Edisonbaan 24
3439 MN Nieuwegein
Phone +31.30.6025080
Fax +31.30.6025099

Spain
Sartorius Stedim Spain SA
C/Isabel Colbrand 10,
Oficina 70
Polígono Industrial de Fuencarral
28050 Madrid
Phone +34.90.2110935
Fax +34.91.3589623

Switzerland
Sartorius Stedim Switzerland GmbH
Lerzenstrasse 21
8953 Dietikon
Phone +41.44.741.05.00
Fax +41.44.741.05.09

U.K.
Sartorius Stedim UK Limited
Longmead Business Park
Blenheim Road, Epsom
Surrey KT19 9 QQ
Phone +44.1372.737159
Fax +44.1372.726171

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Asia | Pacific

China
Sartorius Stedim Beijing
Representative Office
No. 33, Yu'an Road,
Airport Industrial Zone B, Shunyi District
Beijing 101300
Phone +86.10.80426516
Fax +86.10.80426580

Sartorius Stedim Shanghai
Representative Office
Room 618, Tower 1, German Centre,
Shanghai, PRC., 201203
Phone +86.21.28986393
Fax +86.21.28986392.11

Sartorius Stedim Guangzhou Office
Room 704, Broadway Plaza,
No. 233-234 Dong Feng West Road
Guangzhou 510180
Phone +86.20.8351.7921
Fax +86.20.8351.7931


India
Sartorius Stedim India Pvt. Ltd.
10, 6th Main, 3rd Phase Peenya
KIADB Industrial Area
Bangalore – 560 058
Phone +91.80.2839.1963|0461
Fax +91.80.2839.8262

Japan
Sartorius Stedim Japan K.K.
KY Building, 8-11
Kita Shinagawa 1-chome
Shinagawa-ku
Tokyo 140-0001
Phone +81.3.3740.5407
Fax +81.3.3740.5406

Malaysia
Sartorius Stedim Malaysia Sdn. Bhd.
Lot L3-E-3B, Enterprise 4
Technology Park Malaysia
Bukit Jalil
57000 Kuala Lumpur
Phone +60.3.8996.0622
Fax +60.3.8996.0755

Singapore
Sartorius Stedim Singapore Pte. Ltd.
10, Science Park Road, The Alpha
#02-25, Singapore Science Park 2
Singapore 117684
Phone +65.6872.3966
Fax +65.6778.2494

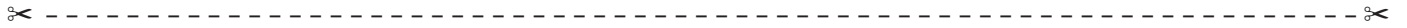Australia
Sartorius Stedim Australia Pty. Ltd.
Unit 5, 7-11 Rodeo Drive
Dandenong South Vic 3175
Phone +61.3.8762.1800
Fax +61.3.8762.1828

Non-hazardous Declaration

Declaration of Safety and Acceptability as Non-Hazardous to Health

The safety and health of our employees, the legal regulations governing the handling of materials that pose hazards to health and the environment and the regulations governing occupational safety mandate and require that this Non-hazardous Declaration be filled out completely for all equipment returned to the respective Sartorius Stedim Biotech affiliate, subsidiary, dealer or distributor listed for your country. No repair or return is possible unless a completed Non-hazardous Declaration form has been submitted.

Fax or mail a copy of a completed Non-hazardous Declaration form in advance to the respective Sartorius Stedim Biotech affiliate, subsidiary, dealer or distributor listed for your country (for the fax number or address, please refer to our list of countries). We require this information before the equipment | component arrives. An additional copy must accompany the equipment | component. If necessary, the shipping company must be notified.

Incomplete entries or non-compliance with this procedure will automatically lead to considerable delays in processing.

In the event of return, we do not differentiate between

- still utilizable and | or repairable used electrical and electronic equipment and no longer utilizable and | or repairable used electrical and electronic equipment or
- between waste that is subject to monitoring by the supervisory authorities and waste that is not subject to any special monitoring by supervisory authorities (used electrical and electronic equipment).

| Model | Serial No. |
|--|------------|
| <p>I we guarantee that the following measures have been taken:</p> <ul style="list-style-type: none"><input type="checkbox"/> The equipment component has been freed of hazardous materials to ensure that persons involved with the handling repair are not exposed to any hazard or danger whatsoever.<input type="checkbox"/> The equipment has been packaged for safety and the complete labeling has been affixed to it.<input type="checkbox"/> The shipping agent has been notified (as prescribed) about the hazardous nature of the shipment. | |
| <p>The person sending in the equipment component is hereby notified that said person shall be liable to Sartorius Stedim Biotech, its affiliates, subsidiaries, dealers and distributors as well as to any third parties – in particular to any of employees of Sartorius, its affiliates, its subsidiaries, dealers or distributors involved with the handling repair of the equipment component – for any damage, caused by entries that are incomplete or incorrect on account of negligence, gross negligence or willful intent.</p> | |
| <p>Name _____</p> <p>Position _____</p> <p>Date _____</p> <p>Signature _____</p> <p>Company stamp _____</p> | |

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Sartorius Stedim Nordic A/S
Hoerskaetten 6D, 1.
DK-2630 Taastrup

Telefon +45.7023.4400
Fax +45.4630.4030

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Sartorius Stedim UK Limited
Longmead Business Park
Blenheim Road, Epsom
Surrey KT19 9 QQ

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Asien | Pazifik

China

Sartorius Stedim Beijing
Representative Office
No. 33, Yu'an Road,
Airport Industrial Zone B, Shunyi District
Beijing 101300

Telefon +86.10.80426516
Fax +86.10.80426580

Sartorius Stedim Shanghai
Representative Office
Room 618, Tower 1, German Centre,
Shanghai, PRC., 201203

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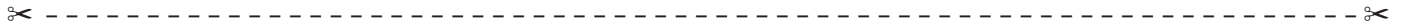
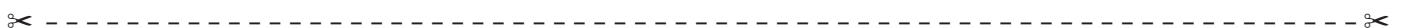
Sartorius Stedim Biotech GmbH
August-Spindler-Str. 11
37079 Goettingen, Germany

Phone +49.551.308.0
Fax +49.551.308.3289
www.sartorius-stedim.com

Copyright by
Sartorius Stedim Biotech GmbH,
Goettingen, Germany.
All rights reserved. No part of
this publication may be reprinted
or translated in any form or by any
means without the prior
written permission of Sartorius
Stedim Biotech GmbH.
The status of the information,
specifications and illustrations
in this manual is indicated by
the date given below.
Sartorius Stedim Biotech GmbH
reserves the right to make changes
to the technology, features,
specifications and design of the
equipment without notice.

Status:
November 2009,
Sartorius Stedim Biotech GmbH,
Goettingen, Germany

DOMINIQUE DUTSCHER SAS