



sartorius

Directions for Use

Vivaflow 50|50R|200

DOMINIQUE DUTSCHER SAS



85030-519-25

## Contents

|                                          |         |
|------------------------------------------|---------|
| <b>Vivaflow 50</b>                       | 3       |
| Single Module                            | 3       |
| Linked Modules                           | 4       |
| Diafiltration   Desalting                | 5       |
| Concentration                            | 6       |
| Recovery                                 | 6       |
| System Components                        | 7       |
| Technical Specifications                 | 7       |
| Performance Characteristics              | 8       |
| Ordering Information                     | 9       |
| Chemical Compatibility                   | 10      |
| <br><b>Vivaflow 50R and Vivaflow 200</b> | <br>11  |
| Operation of Single Module               | 11      |
| Operation of Two Modules                 | 12 – 13 |
| Diafiltration   Desalting                | 14      |
| Concentration                            | 14      |
| Recovery                                 | 14      |
| Technical Specifications                 | 15      |
| Cleaning                                 | 16      |
| Storage                                  | 16      |
| Performance Characteristics              | 17 – 18 |
| Ordering Information                     | 19 – 20 |
| Chemical Compatibility                   | 21      |

# Vivaflow 50

Designed for **single use** only. Anything else constitutes improper use.

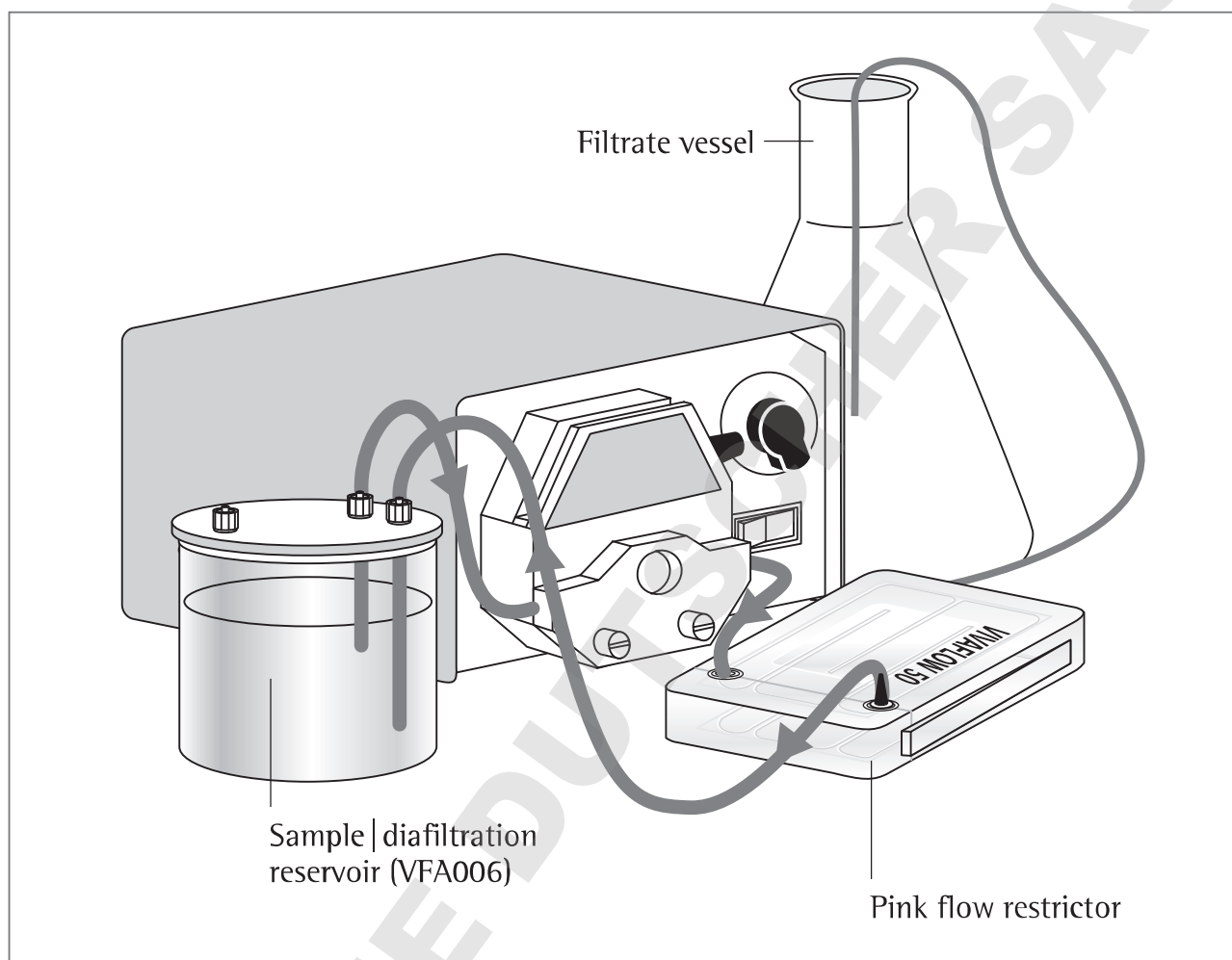


Diagram 1: Vivaflow 50 set up

## Single module

1. Set up the system as illustrated in Diagram 1. Note the positioning of the flow restrictor on the return line.
2. Vivaflow membranes contain trace amounts of glycerine and sodium azide. To remove these chemicals and to check the security of the tube connections, it is recommended to rinse the module, and to test the system at full pressure before introducing the sample.
3. Place 500 ml deionised water in a suitable reservoir. The Vivaflow sample | diafiltration reservoir (prod. no. VFA006) is recommended.
4. Pump liquid through the system to purge any air pockets. The recirculation rate should be in the range 200–400 ml/min, and suitable flow should exit the filtrate line. If used, the pressure indicator (prod. no. VFA020), should read approximately 2.5 bar.
5. Allow 400 ml to pass into the filtrate vessel. Check for any leakage at tubing connection points. Drain the system and empty or replace the filtrate vessel, (see recovery section). The system is now ready for use.

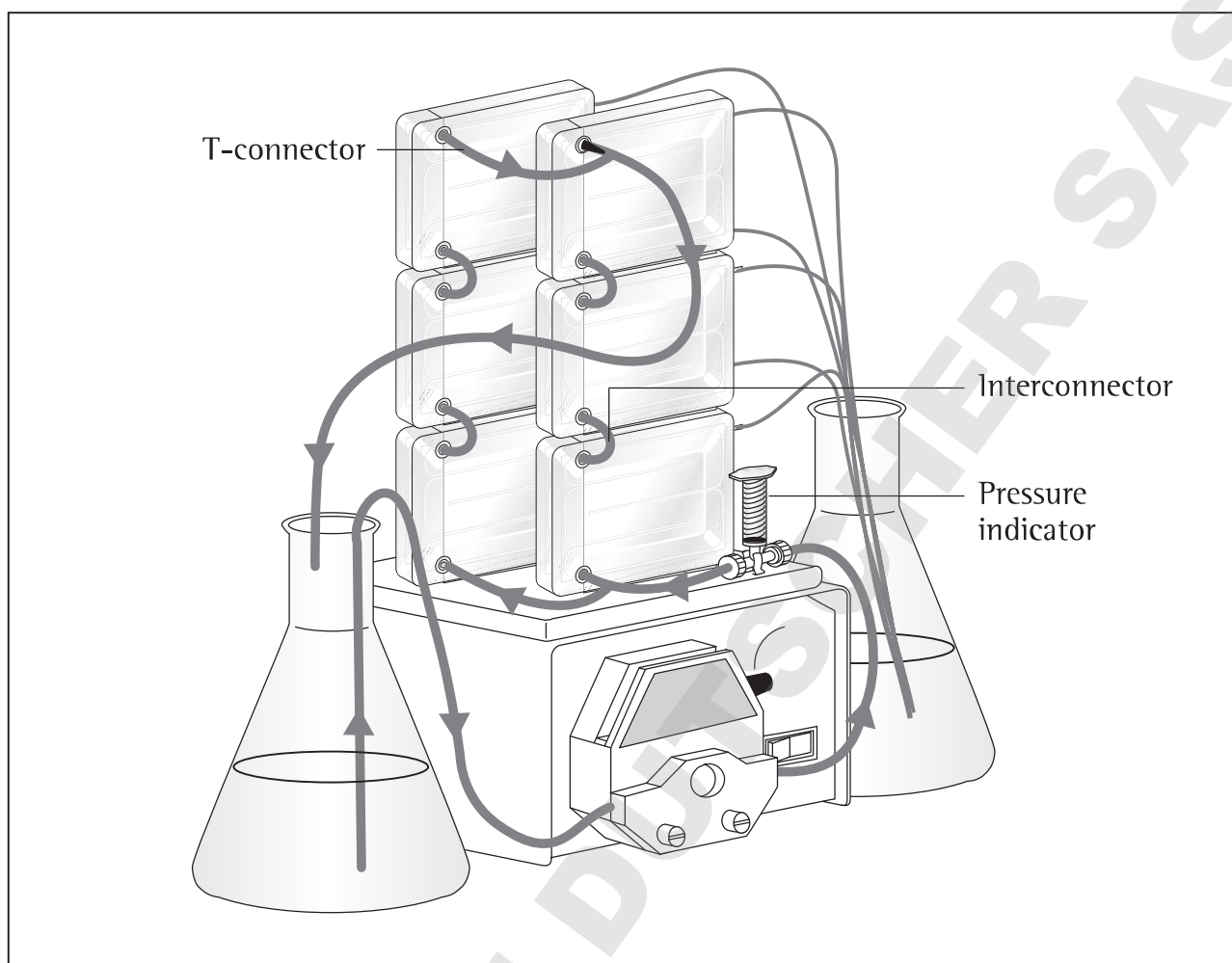


Diagram 2: Setting up multiple Vivaflow 50 cassettes

### Linked modules

1. Set up the system as illustrated in Diagram 2. Link the required number of modules by sliding the tongue and groove edges together. Note the positioning of the flow restrictor on the return line.
2. Use series interconnectors, (prod. no. VFA031), to connect a single row of modules. Detailed ordering information for system components to be found on page 7, Diagram 4.
3. In addition, use T-connectors, (prod. no. VFA030), to connect two rows of modules in parallel.
4. The flow path of the system can be configured to suit the membrane|sample combination. Most solutions are better suited by modules connected in parallel. Very low viscosity solutions favour modules connected only in series. Rinse the system as detailed under single module operation.

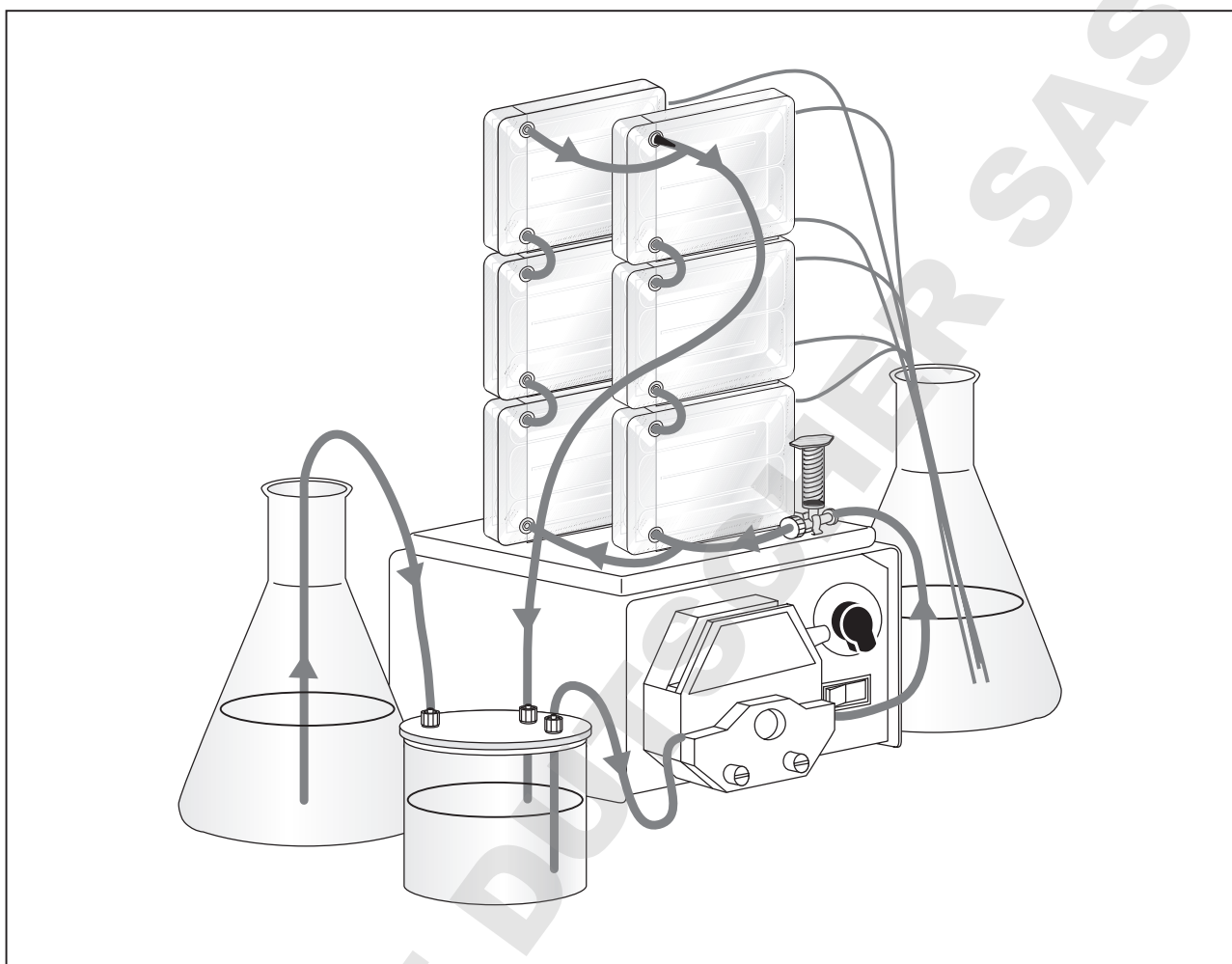


Diagram 3: Buffer exchange | diafiltration with Vivaflow 50

### **Diafiltration | Desalting**

1. Set up the system as illustrated in Diagram 3 for linked modules. If only one module is required, connect it in place of the linked modules.
2. Rinse the system as detailed previously.
3. Either place the solution to be purified in the 500 ml sample | diafiltration reservoir (prod. no. VFA006), or concentrate a larger volume to 500 ml or less so that it is fully contained in the 500 ml reservoir. Ensure the reservoir lid is firmly closed.
4. Fill a large feed reservoir with exchange solvent and pump sample through the system as for concentration.
5. As the volume in the 500 ml reservoir decreases, the vacuum created draws exchange solvent through the feed line from the larger reservoir. Over 99% solvent exchange can be accomplished with an exchange volume approximately 5 times the volume of the sample.

## Concentration

1. Fill the feed reservoir with sample solution. When initial volumes larger than 500 ml are required, place 500 ml in the sample | diafiltration reservoir, the remaining volume in another suitable container and connect the vessels as detailed in the diafiltration section. Alternatively, use a larger container for the entire sample volume and immerse the feed and return lines directly into the liquid.
2. Pump liquid through the system. The recirculation rate should be in the range 200–400 ml/min, and suitable flow should exit the filtrate line. If used, the pressure indicator, (prod. no. VFA020), should read approximately 2.5 bar.
3. Concentrate the sample to the desired volume.

⚠ **Warning:** Do not run the same section of tubing through the pump head for longer than six hours, over use of tubing will result in significant pressure drop and ultimately, failure.

4. When the desired volume has been reached, reduce the recirculation rate to 20–40 ml/min and recirculate the concentrated sample for 1–2 minutes to maximise recovery.

## Recovery

1. Disconnect the feed line from the lid of the 500 ml reservoir or when using a different container, remove the feed line from the sample.
2. Pump residual system volume back into the reservoir | container. (When parallel modules are used with viscous solutions, ensure that all modules are empty by pinching the tubing between each of the stacks of modules in turn).
3. For a more complete sample recovery rinse approximately 5–10 ml per module of water or sample buffer through the system, and recover as before.

## System Components

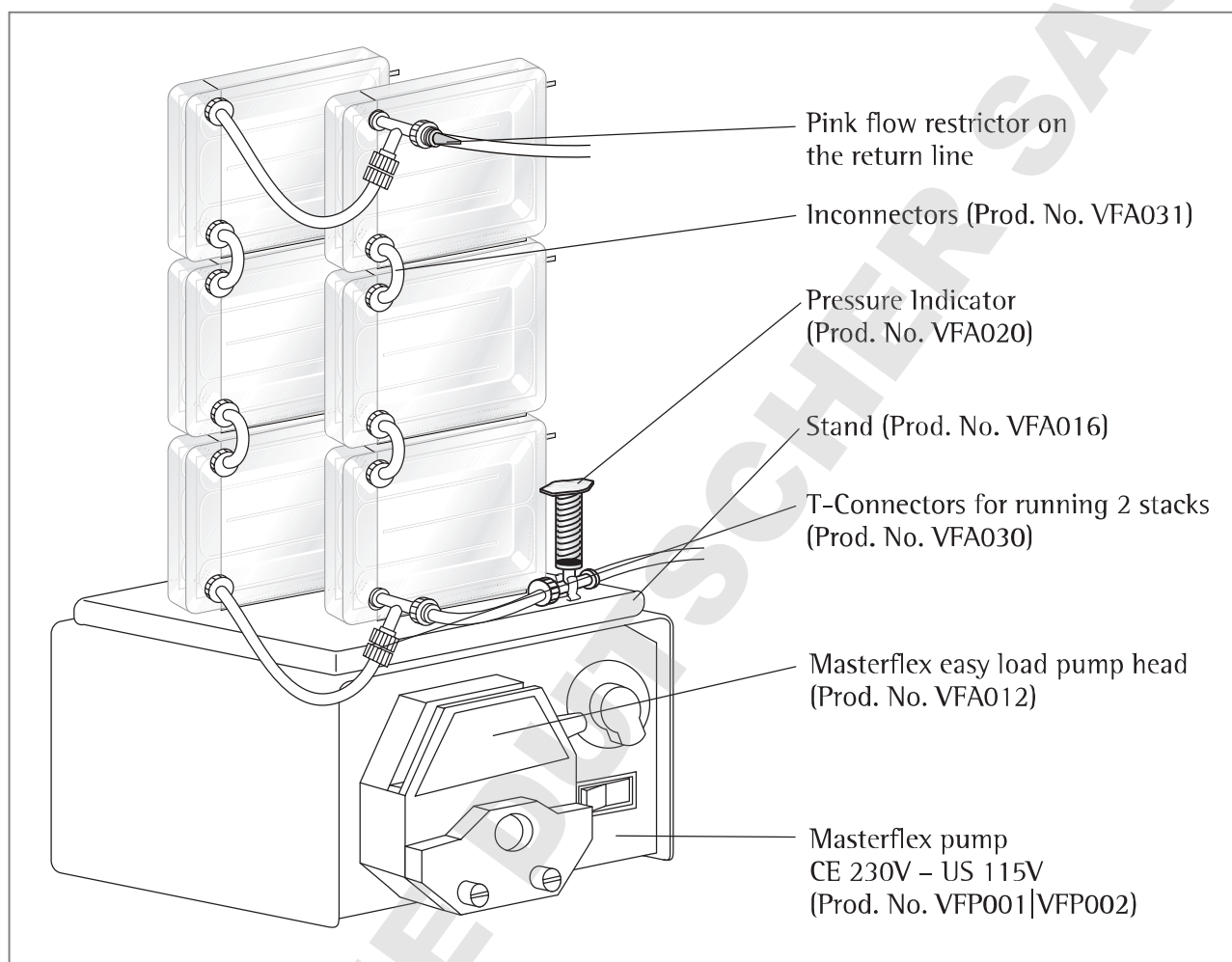


Diagram 4: System components for setting up multiple Vivaflow 50 units

**Table 1: Technical Specifications**

### Dimensions

|                          |                    |
|--------------------------|--------------------|
| Overall L   H   W        | 107   84   25 mm   |
| Channel W   H            | 15 mm   0.3 mm     |
| Active membrane area     | 50 cm <sup>2</sup> |
| Hold up volume (module)  | 1.5 ml             |
| Min recirculation volume | < 10 ml            |
| Non recoverable hold-up  | < 0.5 ml           |

### Operating Conditions

|                     |                |
|---------------------|----------------|
| Pump flow           | 200–400 ml/min |
| Maximum pressure    | 3 bar (45 psi) |
| Maximum temperature | 60 °C          |

## Materials of Construction

|                    |                           |
|--------------------|---------------------------|
| Main housing       | Polycarbonate             |
| Flow channel       | TPX (PMP)                 |
| Membrane support   | TPX (PMP)                 |
| Seals and O rings  | Silicone                  |
| Pressure indicator | Polypropylene, SS Spring, |
| Flow restrictor    | Polypropylene             |
| Fittings           | Nylon                     |
| Tubing             | PVC (medical grade)       |

**Table 2: Performance characteristics**

|                                       | Time to concentrate up to 20x (min.) at 3 bar inlet pressure, 20°C |                     |                   |             |
|---------------------------------------|--------------------------------------------------------------------|---------------------|-------------------|-------------|
|                                       | Single Device                                                      | Three Devices       | Solute Recovery % |             |
|                                       | 250 ml Start Vol. min.                                             | 1 L Start Vol. min. | Direct            | 10 ml rinse |
| <b>BSA 1.0 mg/ml (66,000 MW)</b>      |                                                                    |                     |                   |             |
| 5,000 MWCO PES                        | 34                                                                 | 49                  | 96%               | > 99%       |
| 10,000 MWCO PES                       | 22                                                                 | 32                  | 94%               | > 99%       |
| 30,000 MWCO PES                       | 22                                                                 | 32                  | 92%               | 99%         |
| 50,000 MWCO PES                       | 20                                                                 | 29                  | 92%               | 98%         |
| <b>γ Globulins 1.0 mg/ml</b>          |                                                                    |                     |                   |             |
| 100,000 MWCO PES                      | 43                                                                 | 62                  | 92%               | 98%         |
| 100,000 MWCO RC                       | 40                                                                 | 58                  | 92%               | 98%         |
| 100,000 MWCO Hydrosart                | 40                                                                 | 58                  | 92%               | 98%         |
| <b>Yeast 1.0 mg/ml (S.Cerevisiae)</b> |                                                                    |                     |                   |             |
| 0.2 μm PES                            | 33                                                                 | 47                  | 92%               | 98%         |

## Ordering Information

| <b>Vivaflow 50 include filtrate tube, size 16 peristaltic tubing, flow restrictor and fittings and 1 × Series Interconnector</b> | <b>Pack Size</b> | <b>Prod. No.</b> |
|----------------------------------------------------------------------------------------------------------------------------------|------------------|------------------|
| 3,000 MWCO (PES)                                                                                                                 | 2                | VF05P9           |
| 5,000 MWCO (PES)                                                                                                                 | 2                | VF05P1           |
| 10,000 MWCO (PES)                                                                                                                | 2                | VF05P0           |
| 30,000 MWCO (PES)                                                                                                                | 2                | VF05P2           |
| 50,000 MWCO (PES)                                                                                                                | 2                | VF05P3           |
| 100,000 MWCO (PES)                                                                                                               | 2                | VF05P4           |
| 0.2 µm (PES)                                                                                                                     | 2                | VF05P7           |
| 100,000 MWCO (RC)                                                                                                                | 2                | VF05C4           |

### Vivaflow 50 complete system comprises

|                                                                                                                                                                      |   |        |
|----------------------------------------------------------------------------------------------------------------------------------------------------------------------|---|--------|
| Pump (240 V), easy load pump head (size 16), tubing, 500 ml sample   diafiltration reservoir, module stand, pressure indicator, T connectors, series interconnectors | 1 | VFS502 |
| Pump (115 V), easy load pump head (size 16), tubing, 500 ml sample   diafiltration reservoir, module stand, pressure indicator, T connectors, series interconnectors | 1 | VFS504 |

### Vivaflow 50 PVC tubing and fittings

|                                                                                                                                                            |        |
|------------------------------------------------------------------------------------------------------------------------------------------------------------|--------|
| Size 16 PVC pump tubing (3 metres, 3.2 × 1.6 mm)                                                                                                           | VFA004 |
| Flow restrictor set (2 × 0.4, 0.6, 0.8 mm)                                                                                                                 | VFA009 |
| T connectors for running 2 stacks (2 pieces)                                                                                                               | VFA030 |
| Series interconnectors (6 pieces)                                                                                                                          | VFA031 |
| Female luer fittings (10 pieces)                                                                                                                           | VFA032 |
| VF50 tubing kit (2 × 1 m size 16 PVC tubing with inlets fittings, 2 × 50 cm size 16 PVC tubing with 0.6 mm flow restrictors, 1 × series inlet connections) | VFA034 |
| Flow restrictor 0.6 mm (6 pieces)                                                                                                                          | VFA035 |

### Vivaflow 50 accessories

|                                                                  |        |
|------------------------------------------------------------------|--------|
| Masterflex economy drive variable speed peristaltic pump (230 V) | VFP001 |
| Masterflex economy drive variable speed peristaltic pump (115 V) | VFP002 |
| 500 ml sample and   or diafiltration reservoir                   | VFA006 |
| Masterflex easy load pump head – size 15                         | VFA013 |
| Masterflex easy load pump head – size 16                         | VFA012 |
| Vivaflow 50 stand                                                | VFA016 |
| Pressure indicator (1–3 bar)                                     | VFA020 |

**Table 3: Chemical Compatibility  
for the Vivaflow 50**

| <b>Solution</b>           | <b>PES</b> | <b>RC</b> |
|---------------------------|------------|-----------|
| Acetic acid (25%)         | OK         | OK        |
| Acetone                   | NO         | NO        |
| Ammonium hydroxide (5%)   | OK         | OK        |
| Ammonium sulfate sat.     | OK         | OK        |
| DMEM                      | OK         | OK        |
| Ethanol (70%)             | OK         | OK        |
| Ethyl acetate             | NO         | NO        |
| Formaldehyde (30%)        | OK         | OK        |
| Formic acid (5%)          | OK         | OK        |
| Guanidine HCl (6 M)       | OK         | OK        |
| Hydrocarbons, Aromatic    | NO         | ?         |
| Hydrocarbons, Chlorinated | NO         | NO        |
| Hydrochloric acid (1 M)   | OK         | NO        |
| Isopropanol               | NO         | NO        |
| Lactic acid (5%)          | OK         | OK        |
| Mercaptoethanol (10 mM)   | OK         | OK        |
| Methanol (60%)            | OK         | OK        |
| n-Butanol (70%)           | OK         | OK        |

| <b>Solution</b>             | <b>PES</b> | <b>RC</b> |
|-----------------------------|------------|-----------|
| Peracetic acid (0.2%)       | OK         | OK        |
| Phenol (1%)                 | OK         | ?         |
| Phosphate Buffer (1 M)      | OK         | OK        |
| Pyridine                    | NO         | NO        |
| RPMI-1640                   | OK         | OK        |
| Sodium azide                | OK         | OK        |
| Sodium deoxycholate (5%)    | OK         | OK        |
| Sodium hydroxide (1 M)      | NO         | NO        |
| Sodium hydroxide (0.1 M)    | NO         | NO        |
| Sodium hypochlorite (0.02%) | ?          | ?         |
| Sodium nitrate              | NO         | NO        |
| Sulfamic acid (5%)          | OK         | NO        |
| Surfactants (0.1%)          | OK         | OK        |
| Toluene                     | NO         | NO        |
| Trichloroacetic acid (10%)  | NO         | ?         |
| Trifluoroacetic acid (10%)  | OK         | NO        |
| Urea (8 M)                  | OK         | OK        |

OK = Acceptable

? = Questionable

NO = Not recommended

## Vivaflow 50R and Vivaflow 200

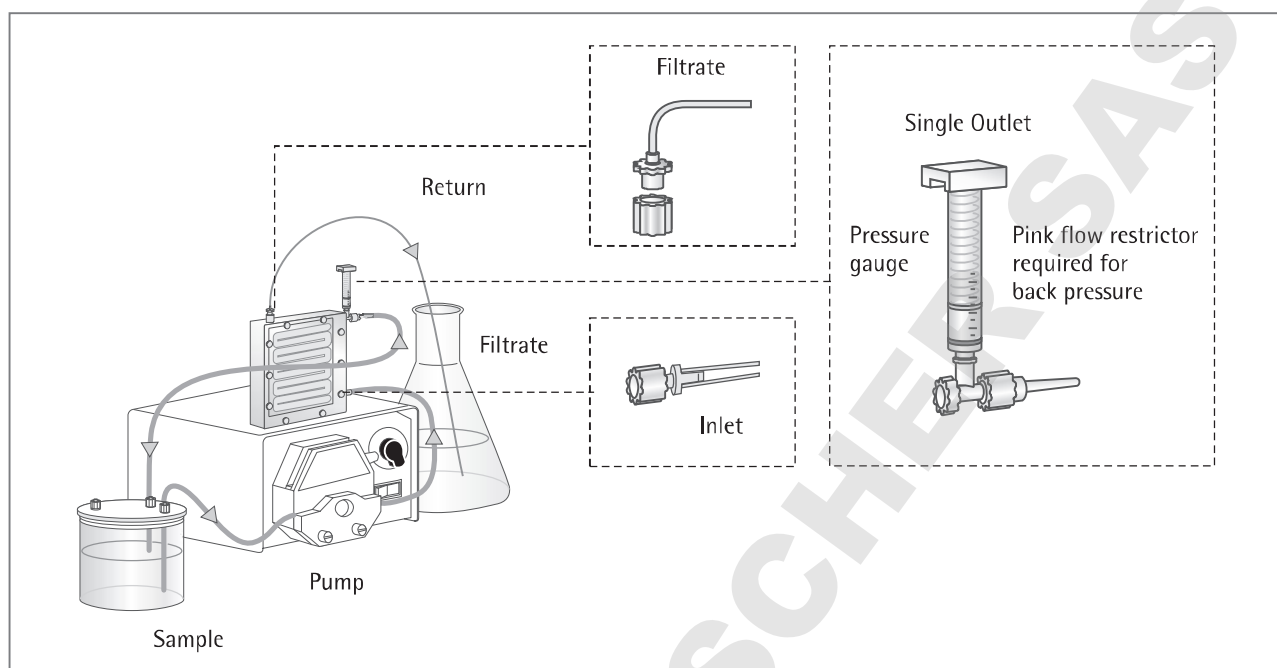


Diagram 5: Vivaflow 50R and Vivaflow 200 set up

### Operation Single Vivaflow 50R or Vivaflow 200 Modules

#### Assembly and rinsing single module

1. Set up the system as illustrated above in Diagram 5.. Note the positioning of the flow restrictor on the return line. A set of flow restrictors of different gauges (Product No. VFA009), is available for use with solutions of unusually high or low viscosity.
2. Vivaflow membranes contain trace amounts of glycerine and sodium azide. To remove these chemicals and to check the security of the tube connections, it is recommended to rinse the module, and to test the system at full pressure before introducing the sample.
3. Place 500 ml deionised water in a suitable reservoir.
4. Pump liquid through the system to purge any air pockets. The recirculation rate should be 200–400 ml/min, and suitable flow should exit the filtrate line. If used, the pressure indicator should read approximately 2.5 bar.
5. Allow 400 ml to pass into the filtrate vessel. Check for any leakage at tubing connection points.
6. Drain the system and empty or replace the filtrate vessel (see Recovery section). The system is now ready for use.

**⚠ Warning:** Ensure Luer connections are secure before operation.

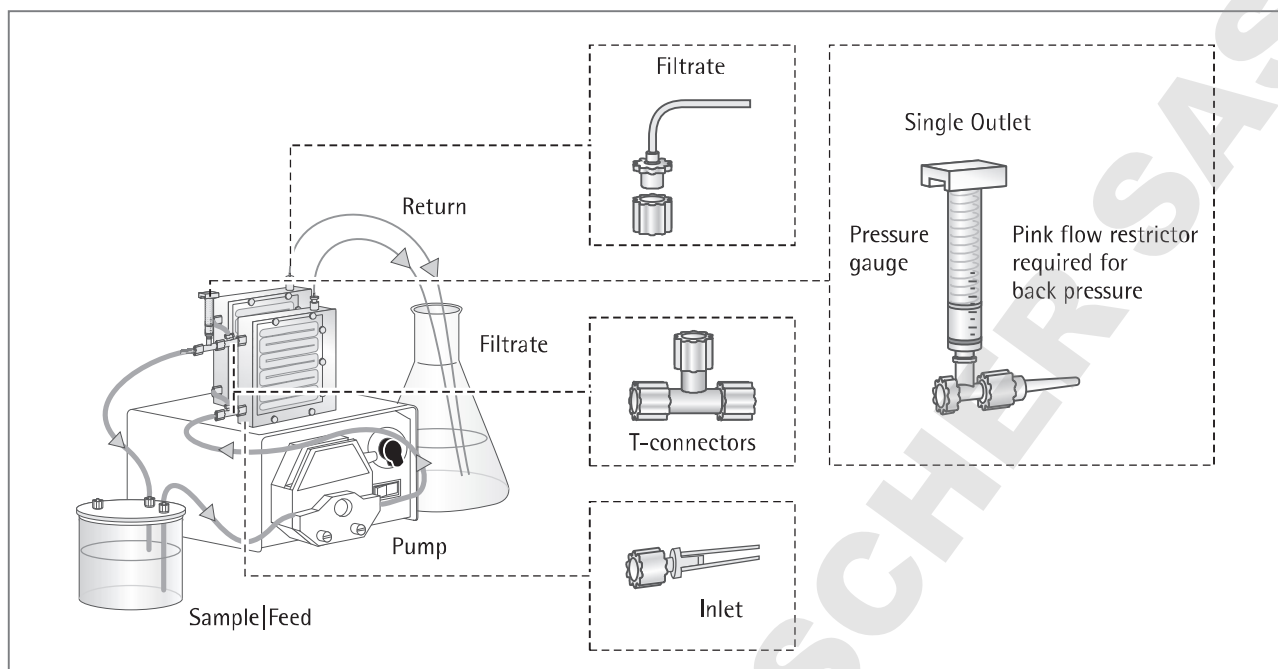


Diagram 6: Setting up two Vivaflow 50R units

## Operating two Vivaflow 50R or two Vivaflow 200 Modules

### Assembly and rinsing two modules

1. Set up the Vivaflow 50R system as illustrated in Diagram 6 and the Vivaflow 200 System as illustrated in Diagram 7 system as illustrated above.

**⚠ Warning:** Ensure Luer connections are secure before operation.

### 2. Vivaflow 50R:

Use two T-connectors (VFA030) to connect to the inlets of both Vivaflow 50R modules.

### Vivaflow 200:

Use the Y-connector, (VFA005) to connect to the inlets of both Vivaflow 200 modules.

The size 15 tubing connected to the Y-connector is the inlet tubing for the pump. Use a size 15 pump head.

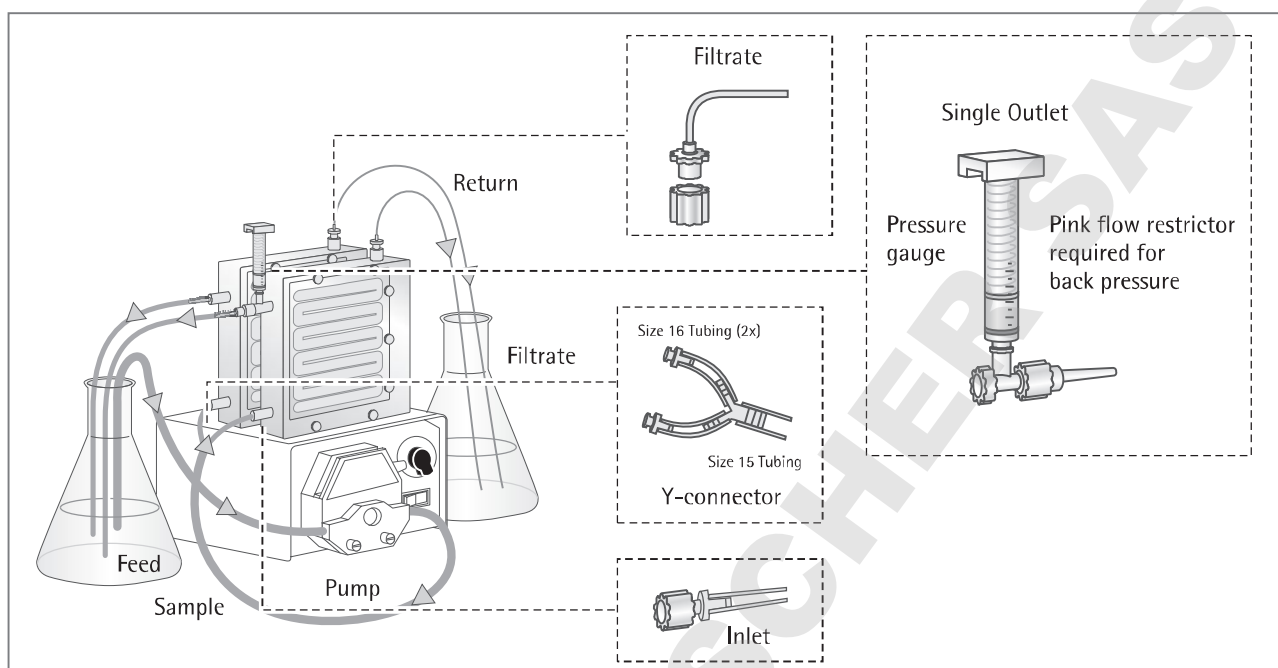


Diagram 7: Setting up two Vivaflow 200 units

3. Note the positioning of a flow restrictor on the return line from each Vivaflow 200 module. Vivaflow membranes contain trace amounts of glycerine and sodium azide. To remove these chemicals and to check the security of the tube connections, it is recommended to rinse the module, and to test the system at full pressure before introducing the sample.
4. Place 1 litre of deionised water in a suitable reservoir.
5. Pump liquid through the system to purge any air pockets. The recirculation rate should be 500–900 ml/min, and suitable flow should exit the filtrate line. If used, the pressure indicator should read approximately 2.5 bar.
6. Allow 800 ml to pass into the filtrate vessel. Check for any leakage at tubing connection points.
7. Drain the system and empty or replace the filtrate vessel (see Recovery section). The system is now ready for use.

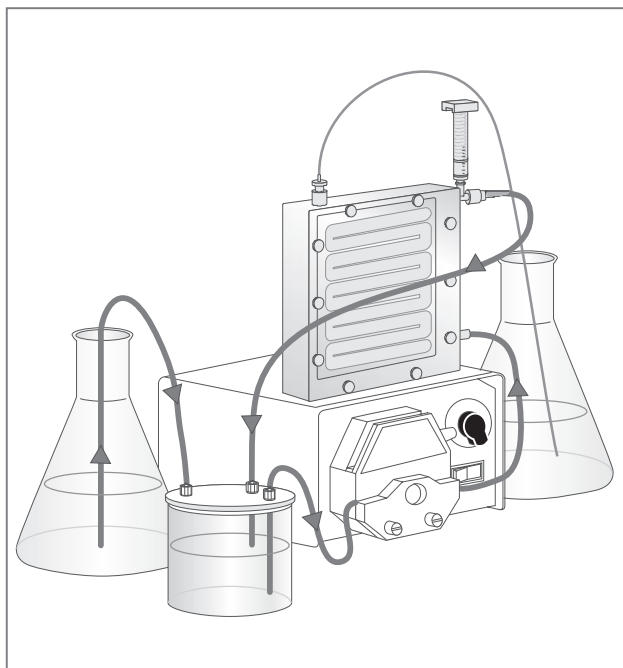


Diagramm 8: Buffer exchange | Desalting with Vivaflow 50R and Vivaflow 200

### Diafiltration | Desalting

1. Set up the system as illustrated in Diagram 8.
2. Rinse the system as detailed previously.
3. Place the solution to be purified in a suitable diafiltration reservoir. For volumes up to 500 ml the Vivaflow 500 ml sample | diafiltration reservoir (VFA006), is recommended. Ensure the lid is firmly closed.
4. Fill the large feed reservoir with exchange solvent and pump sample through the system as for concentration.
5. As liquid passes through the membrane, the vacuum created in the diafiltration reservoir draws in exchange solvent through the feed line from the large reservoir. Over 99% solvent exchange can be accomplished with an exchange volume of approximately 5 times the sample volume.

### Concentration

1. Fill the feed reservoir with sample solution.
2. Pump liquid through the system. The recirculation rate should be 200–400 ml/min, (500–900 ml/min for two modules), and suitable flow should exit the filtrate line. If used, the pressure indicator should read approximately 2.5 bar.

**⚠ Warning:** Do not run the same section of tubing through the pump-head for longer than 6 hours, overuse of tubing will result in significant pressure drop and ultimately, failure.

3. Concentrate the sample.
4. When nearing the desired volume, reduce the recirculation rate to 20–40 ml/min and recirculate the concentrated sample for 1–2 minutes to increase sample recovery.

### Recovery

1. Remove the feed line from the sample.
2. Pump residual system volume back into the reservoir | container.
3. For a more complete recovery, rinse 25–50 ml of water or sample buffer through the system, and recover as before.

**Table 4: Technical specifications Vivaflow 50R and Vivaflow 200**

| <b>Dimensions</b>                | <b>Vivaflow 50 R</b>                                                  | <b>Vivaflow 200</b> |
|----------------------------------|-----------------------------------------------------------------------|---------------------|
| Overall L   H   W                | 100   100   24 mm                                                     | 126   138   38 mm   |
| Channel W   H                    | 7.5   0.4 mm                                                          | 10 mm   0.4 mm      |
| Active membrane area             | 50 cm <sup>2</sup>                                                    | 200 cm <sup>2</sup> |
| Hold up volume (module)          | 1.7 ml                                                                | 5.3 ml              |
| Min. recirculation volume        | 10 ml                                                                 | < 20 ml             |
| Non recoverable hold-up          | < 0.5 ml                                                              | < 1 ml              |
| <b>Materials of construction</b> | <b>Vivaflow 50R and Vivaflow 200</b>                                  |                     |
| Main housing                     | Acrylic                                                               |                     |
| Flow channel                     | Acrylic                                                               |                     |
| Membrane support                 | Polypropylene                                                         |                     |
| Seals and O rings                | Silicone                                                              |                     |
| Pressure indicator               | Polypropylene, SS spring                                              |                     |
| Flow restrictor                  | Polypropylene                                                         |                     |
| Fittings                         | Nylon                                                                 |                     |
| Tubing                           | PVC (medical grade)                                                   |                     |
| <b>Operating conditions</b>      | <b>Vivaflow 50R and Vivaflow 200</b>                                  |                     |
| Pump flow                        | 200–400 ml/min                                                        |                     |
| Maximum pressure                 | 4 bar (60 psi)<br>Pressure drop across inlet   outlet 0.5 bar (7 psi) |                     |
| Maximum temperature              | 60 °C                                                                 |                     |

## Cleaning and storage information only for Vivaflow 50R and Vivaflow 200

### Cleaning

Vivaflow 50R and 200 modules may be used several times if cleaning and storage instructions are followed.

1. Flush the system with 200 ml of deionised water with the filtrate going to waste.
2. Place the feed, return and filtrate lines in a suitable container for the cleaning solution.
3. Prepare cleaning solutions suitable for the membrane.
  1. Polyethersulfone membranes:  
250 ml of 0.5 mM NaOCl in 0.5 M NaOH
  - II. Hydrosart membranes: 250 ml of 0.5 M NaOH
4. Recirculate at 50–100 ml/min for 30–40 minutes.
5. Drain the system and recirculate 250 ml of deionised water through the system for 5–10 minutes.
6. Drain and discard rinse solution and flush with a further 500 ml of deionised water with the filtrate going to waste. The system is now ready for further use.

### Storage

To store Vivaflow 50R and Vivaflow 200 after cleaning, fill module with deionised water and 10% ethanol. Seal inlet, outlet and filtrate ports and refrigerate at approximately 4°C.

**Table 5: Performance characteristics Vivaflow 50R**

|                                                                                           | Time to concentrate up to 20x (min)<br>at 3.0 bar inlet   2.5 bar outlet pressure, 20°C |              |            |             |
|-------------------------------------------------------------------------------------------|-----------------------------------------------------------------------------------------|--------------|------------|-------------|
|                                                                                           | start volume                                                                            | Average flux | Recovery % |             |
|                                                                                           | 250 ml                                                                                  | ml/min       | Direct     | 25 ml rinse |
| <b>Lysozyme 0.25 mg/ml (14,000 MW)</b>                                                    |                                                                                         |              |            |             |
| 5,000 MWCO Hydrosart                                                                      | 70                                                                                      | 3.4          | 96%        | 98%         |
| 10,000 MWCO Hydrosart                                                                     | 23                                                                                      | 10.3         | 94%        | 96%         |
| <b>BSA 1.0 mg/ml (66,000 MW)</b>                                                          |                                                                                         |              |            |             |
| 10,000 MWCO Hydrosart                                                                     | 24                                                                                      | 9.9          | 98%        | > 99%       |
| 30,000 MWCO Hydrosart                                                                     | 15                                                                                      | 15.8         | 97%        | > 99%       |
| <b>γ Globulins 1.0 mg/ml (150,000 MW)</b>                                                 |                                                                                         |              |            |             |
| 100,000 MWCO Hydrosart                                                                    | 46                                                                                      | 5.2          | 97%        | > 99%       |
| <b>Start volume 1 litre (one Vivaflow 50R at 3 bar) 10,000 MWCO Hydrosart</b>             |                                                                                         |              |            |             |
| BSA 1.0 mg/ml                                                                             | 95                                                                                      | 10.0         | 98%        | > 99%       |
| <b>Start volume 1 litre (two Vivaflow 50R in parallel at 3 bar) 10,000 MWCO Hydrosart</b> |                                                                                         |              |            |             |
| BSA 1.0 mg/ml                                                                             | 48                                                                                      | 19.8         | 98%        | > 99%       |

**Table 6: Performance characteristics Vivaflow 200**

|                                                                                      | Time to concentrate up to 20x (min.) at 3 bar inlet pressure, 20°C |              |            |             |
|--------------------------------------------------------------------------------------|--------------------------------------------------------------------|--------------|------------|-------------|
|                                                                                      | start volume                                                       | Average flux | Recovery % |             |
|                                                                                      | 1 litre                                                            | ml/min       | Direct     | 25 ml rinse |
| <b>Lysozyme 0.25 mg/ml (14,000 MW)</b>                                               |                                                                    |              |            |             |
| 2,000 MWCO Hydrosart                                                                 | 160                                                                | 6            | 97%        | > 99%       |
| 3,000 MWCO PES                                                                       | 180                                                                | 5            | 97%        | > 99%       |
| <b>BSA 1.0 mg/ml (66,000 MW)</b>                                                     |                                                                    |              |            |             |
| 5,000 MWCO PES                                                                       | 29                                                                 | 33           | 98%        | > 99%       |
| 5,000 MWCO Hydrosart                                                                 | 70                                                                 | 14           | 98%        | > 99%       |
| 10,000 MWCO PES                                                                      | 23                                                                 | 41           | 96%        | > 99%       |
| 10,000 MWCO Hydrosart                                                                | 35                                                                 | 27           | 98%        | > 99%       |
| 30,000 MWCO PES                                                                      | 25                                                                 | 38           | 96%        | 99%         |
| 30,000 MWCO Hydrosart                                                                | 20                                                                 | 48           | 96%        | > 99%       |
| 50,000 MWCO PES                                                                      | 22                                                                 | 43           | 96%        | 98%         |
| <b>γ Globulins 1.0 mg/ml (average 160,000 MW)</b>                                    |                                                                    |              |            |             |
| 100,000 MWCO PES                                                                     | 54                                                                 | 18           | 96%        | 99%         |
| <b>Yeast 1.0 mg/ml (S. Cerevisiae)</b>                                               |                                                                    |              |            |             |
| 0.2 µm PES                                                                           | 11                                                                 | 86           | 92%        | 98%         |
| <b>Dilute solute concentration, start volume 1 litre at 3 bar, 10,000 MWCO PES</b>   |                                                                    |              |            |             |
| BSA 0.001 mg/ml                                                                      | 18                                                                 | 52           | 90%        | 98%         |
| BSA 0.01 mg/ml                                                                       | 20                                                                 | 47           | 92%        | 98%         |
| BSA 0.1 mg/ml                                                                        | 21                                                                 | 45           | 94%        | 99%         |
| <b>Start volume 5 litres (two Vivaflow 200 in parallel at 3 bar) 10,000 MWCO PES</b> |                                                                    |              |            |             |
| BSA 1.0 mg/ml (66,000 MW)                                                            | 67                                                                 | 70           | 97%        | > 99%       |

## Ordering Information

| <b>Vivaflow 50R modules include pressure indicator, flow restrictor and size 16 pvc peristaltic tubing and fittings</b> | <b>Qty per box</b> | <b>Prod. no.</b> |
|-------------------------------------------------------------------------------------------------------------------------|--------------------|------------------|
| 5,000 MWCO Hydrosart                                                                                                    | 1                  | VF05H1           |
| 10,000 MWCO Hydrosart                                                                                                   | 1                  | VF05H0           |
| 30,000 MWCO Hydrosart                                                                                                   | 1                  | VF05H2           |
| 100,000 MWCO Hydrosart                                                                                                  | 1                  | VF05H4           |
| <b>Vivaflow 200 modules include pressure indicator, flow restrictor and size 16 pvc peristaltic tubing and fittings</b> | <b>Qty per box</b> | <b>Prod. no.</b> |
| 3,000 MWCO PES                                                                                                          | 1                  | VF20P9           |
| 5,000 MWCO PES                                                                                                          | 1                  | VF20P1           |
| 10,000 MWCO PES                                                                                                         | 1                  | VF20P0           |
| 30,000 MWCO PES                                                                                                         | 1                  | VF20P2           |
| 50,000 MWCO PES                                                                                                         | 1                  | VF20P3           |
| 100,000 MWCO PES                                                                                                        | 1                  | VF20P4           |
| 0.2 µm PES                                                                                                              | 1                  | VF20P7           |
| 2,000 MWCO Hydrosart                                                                                                    | 1                  | VF20H9           |
| 10,000 MWCO Hydrosart                                                                                                   | 1                  | VF20H0           |
| 30,000 MWCO Hydrosart                                                                                                   | 1                  | VF20H2           |
| 100,000 MWCO Hydrosart                                                                                                  | 1                  | VF20H4           |
| <b>Vivaflow 50R   200 complete system comprises</b>                                                                     |                    |                  |
| Pump (240 V), Easy load pump head (size 16), tubing, 500 ml sample   diafiltration reservoir                            | 1                  | VFS202           |
| Pump (115 V), Easy load pump head (size 16), tubing, 500 ml sample   diafiltration reservoir                            | 1                  | VFS204           |

**Vivaflow 50R | 200 accessories**

|                                                                                   |        |
|-----------------------------------------------------------------------------------|--------|
| Masterflex economy drive variable speed peristaltic pump (230V)                   | VFP001 |
| Masterflex economy drive variable speed peristaltic pump (115V)                   | VFP002 |
| 500 ml sample and/or diafiltration reservoir                                      | VFA006 |
| Masterflex easy load pump head – size 16                                          | VFA012 |
| Masterflex easy load pump head – size 15<br>(only for operating two Vivaflow 200) | VFA013 |
| Size 15 pvc pump tubing and Luer fittings (3 m, 4.8 × 2.6 mm)                     | VFA003 |
| Size 16 pvc pump tubing and Luer fittings (3 m, 3.2 × 1.6 mm)                     | VFA004 |
| Flow restrictor set (2 × 0.4, 0.6, 0.8 mm)                                        | VFA009 |
| Female luer fittings size 16 (10 pieces)                                          | VFA032 |
| Flow restrictors 0.6 mm (6 pieces)                                                | VFA035 |
| Female luer fittings size 15 (10 pieces)                                          | VFA036 |
| Flow restrictor set (2 + 0.4, 0.6, 0.8 mm)                                        | VFA009 |
| Female luer fittings size 16 (10 pieces)                                          | VFA032 |
| Flow restrictors 0.6 mm (6 pieces)                                                | VFA035 |
| Female luer fittings size 15 (10 pieces)                                          | VFA036 |

**Vivaflow 200 tubing and fittings for operation two modules**

|                                                     |        |
|-----------------------------------------------------|--------|
| Y connector (size 15 to 2 × size 16, Luer fittings) | VFA005 |
| Masterflex easy load pump head – size 15            | VFA013 |

**Vivaflow 50R tubing and fittings for operation of two modules**

|             |        |
|-------------|--------|
| T-Connector | VFA030 |
|-------------|--------|

**Table 7: Chemical Compatibility  
for Vivaflow 50R and Vivaflow 200**

| <b>Solution</b>           | <b>PES</b> | <b>HY</b> |
|---------------------------|------------|-----------|
| Acetic acid (25%)         | OK         | OK        |
| Acetone                   | NO         | NO        |
| Ammonium hydroxide (5%)   | OK         | OK        |
| Ammonium sulfate sat.     | OK         | ?         |
| DMEM                      | OK         | OK        |
| Ethanol (70%)             | OK         | OK        |
| Ethyl acetate             | NO         | NO        |
| Formaldehyde (30%)        | OK         | OK        |
| Formic acid (5%)          | OK         | OK        |
| Guanidine HCl (6 M)       | OK         | OK        |
| Hydrocarbons, Aromatic    | NO         | NO        |
| Hydrocarbons, Chlorinated | NO         | NO        |
| Hydrochloric acid (1 M)   | OK         | OK        |
| Isopropanol               | NO         | NO        |
| Lactic acid (5%)          | OK         | OK        |
| Mercaptoethanol (10 mM)   | OK         | OK        |
| Methanol (60%)            | OK         | OK        |
| n-Butanol (70%)           | OK         | OK        |

| <b>Solution</b>             | <b>PES</b> | <b>HY</b> |
|-----------------------------|------------|-----------|
| Peracetic acid (0.2%)       | OK         | NO        |
| Phenol (1%)                 | OK         | NO        |
| Phosphate Buffer (1 M)      | OK         | OK        |
| Pyridine                    | NO         | NO        |
| RPMI-1640                   | OK         | OK        |
| Sodium azide                | OK         | OK        |
| Sodium deoxycholate (5%)    | OK         | ?         |
| Sodium hydroxide (1 M)      | OK         | OK        |
| Sodium hydroxide (0.1 M)    | OK         | OK        |
| Sodium hypochlorite (0.02%) | OK         | NO        |
| Sodium nitrate              | NO         | OK        |
| Sulfamic acid (5%)          | OK         | ?         |
| Surfactants (0.1%)          | OK         | OK        |
| Toluene                     | NO         | NO        |
| Trichloroacetic acid (10%)  | NO         | OK        |
| Trifluoroacetic acid (10%)  | OK         | OK        |
| Urea (8 M)                  | OK         | OK        |

OK = Acceptable

? = Questionable

NO = Not recommended

**DOMINIQUE DUTSCHER SAS**

**DOMINIQUE DUTSCHER SAS**

Sartorius Stedim Lab Ltd  
Sperry Way  
Stonehouse Park  
Gloucestershire  
GL10 3UT, UK

[www.sartorius.com](http://www.sartorius.com)

Copyright by  
Sartorius Lab Instruments  
GmbH & Co. KG, Goettingen,  
Germany.

All rights reserved. No part  
of this publication may  
be reprinted or translated in  
any form or by any means  
without the prior written  
permission of Sartorius Lab  
Instruments GmbH & Co. KG.  
The status of the information,  
specifications and illustrations  
in this manual is indicated  
by the date given below.

Sartorius Lab Instruments  
GmbH & Co. KG reserves the right  
to make changes to the technology,  
features, specifications and design  
of the equipment without notice.

Status:

June 2016,  
Sartorius Lab Instruments  
GmbH & Co. KG,  
Goettingen, Germany

Printed in the EU on paper bleached  
without chlorine.  
Publication No.: SLU6097-e160609  
Ver. 06 | 2016