

IZIbio

IZI DEFROST +

Defrosting Vacuum Cleaner for Freezers

-20°C / -86°C / -152°C

Ref: IZIDEFROST+



Technical Specifications


Voltage (V)	220 -240
PIEC (W)	1200
Grid frequency (Hz)	50-60
Air flow (L/s)	62
Sound level (dBA)	61
Weight (Kg)	10

Directions for use


With IZI DEFROST +, defrosting your freezers has never been so QUICK, CLEAN and EASY.

For the first use:

- Make sure that the HEPA filter is in place
- Connect the hose to the suction outlet
- Adapt the scraper to the tip

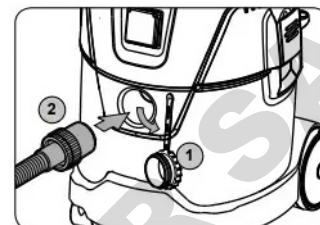

You can then turn on the device, by turning the ON/OFF button one notch to the right, and start vacuuming the frost from your freezer.

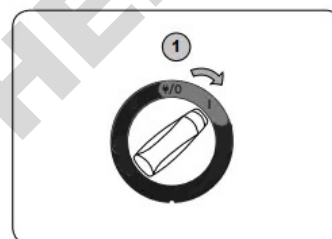

Please empty and clean the tank after each use.

If foam appears, immediately stop and empty the tank.


The device is equipped with a float system which cuts off air flow when the maximum liquid level is reached. You will hear a change in the sound of the motor and notice the suction power decreasing.

When this happens, turn off the machine and unplug it. Never collect liquid if the float is not in place.

Disconnect the hose from the inlet by pulling on it.

Release the latches by pulling them outward to free up the top of the motor. Lift the top of the motor from the tank.

You can empty the tank by tilting it backwards or sideways and pour the liquids into a floor drain or any other appropriate disposal system.

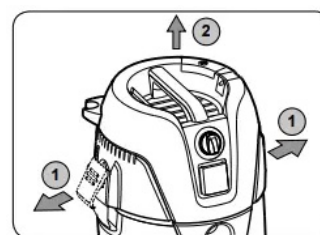

Put the upper part of the motor back in the tank and secure it with the latches.

IZI DEFROST + can be used on -20°C/-30°C, -80°C and -150°C freezers.

Warranty and Maintenance

Our IZI DEFROST + cleaner has a one-year warranty. Batteries, lamps and wear parts are not included.

If a defect occurs during the warranty period, we will either repair the device or replace it.

Please note that the warranty becomes void if the motor casing has been opened.

Changing the HEPA filter:

Turn the filter counterclockwise to remove it and dispose of it in accordance with current waste regulations.

We recommend that you change the filter every year.

(Ref: IZIHEPA.)



IZIBIO

13-15 rue Taitbout - 75 009 PARIS- FRANCE
contact@izibio.fr - 01 48 01 48 14