**AGITADOR MAGNÉTICO ANALÓGICO 20 L
ANALOGIC MAGNETIC STIRRER 20 L
AGITATEUR MAGNETIQUE ANALOGIQUE 20 L**

Serie / Series / Séries / RSLAB-1NC, RSLAB-1C

RS Lab

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This manual should be available for all users of these equipments. To get the best results and a higher duration of this equipment it is advisable to read carefully this manual and follow the processes of use.

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INSTRUCTIONS ON ENVIRONMENT PROTECTION
 At the end of its life cycle, please, does not dispose of this equipment by throwing it in the usual garbage; hand it over a collection point for the recycling of electrical and electronic appliances. It does not contain dangerous or toxic products for humans but a non-adequate disposal would damage the environment.
 The materials are recyclable as mentioned in its marking. By recycling material or by other forms of re-utilization of old appliances, you are making an important contribution to protect our environment.
 Please inquire at the community administration for the authorized disposal location.

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INDEX OF CONTENTS

1. Safety instructions	10
2. Proper use	11
3. Inspection	11
4. Trial run	11
5. Control and display	12
6. Operation	13
7. Faults	13
8. Maintenance and cleaning	13
9. Associated standards and regulation	14
10. Technical data	14
11. Accessories	15

Welcome to the Analog magnetic stirrer RSLAB-1NC (no heating) and RSL RSLAB-1C (heating) user manual. Users should read this manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

When help needed, you can always contact the service department of manufacturer or your supplier for technical support. Please provide the customer care representative with the following information: Serial number, description of problem, methods and procedures adopted to resolve the problems and your contact information.

This instrument is warranted to be free from defects in materials and workmanship under normal use and service for a period of 24 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claims under the warranty please contact your local dealer. You may also send the instrument direct to our works, enclosing the invoice copy and giving reasons for the claim.

TO GET THE BEST RESULTS AND A HIGHER DURATION OF THE EQUIPMENT IT IS ADVISABLE TO READ THOROUGHLY THIS MANUAL BEFORE OPERATING WITH THE EQUIPMENT.

1- SAFETY INSTRUCTIONS

WARNING!

Read the operating instructions carefully before use the instrument.
Ensure that only trained staff work with the instrument.

Forbid to heat the substances with low-burning point or easy-volatile.

Risk of burn!

- ◆ Caution when touching the housing parts and the heating plate. The heating plate can reach temperatures of 340 °C.
- ◆ Pay attention to the residual heat after switching off.

Protective ground contact!

- ◆ Make sure that socket is earthed (protective ground contact) before use.
- ◆ When work, wear the personal guard to avoid the risk from splashing and evaporation of liquids and release of toxic or combustible gases.
- ◆ Set up the instrument in a spacious area on a stable, clean, non-slip, dry and fireproof surface, do not operate the instrument in explosive atmospheres, with hazardous substances or under water.
- ◆ Gradually increase the speed, reduce the speed if the stirring bar breakaway because of too high speed and if the instrument is not running smoothly, or if the container moves on the base plate.
- ◆ Temperature must always be set to at least 25 °C lower than the fire point of the media used.
- ◆ Beware of hazards due to flammable materials or media with a low boiling temperature, overfilling of media and unsafe container.
- ◆ Process pathogenic materials only in closed vessels.
- ◆ The voltage started on the label must correspond to the main power supply.
- ◆ Ensure that the main power supply cable does not touch the hotplate. Do not cover the instrument.
- ◆ The instrument may only be opened by experts.
- ◆ Keep away from high magnetic field.

11- ACCESSORIES (NOT INCLUDED)

[1] 90694301

[2] 90694302


[3] 90694303

[4] 90694304

[5] 90694305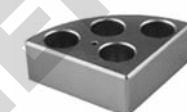

[6] 90694306

[7] 90694307

[8] 90694308

[9] 90694309

[10] 90694310

[11] 90694311

[12] 90694311

[13] 90694311

Reference

- [1] 90694301
[2] 90694302
[3] 90694303
[4] 90694304
[5] 90694305
[6] 90694306
[7] 90694307
[8] 90694308
[9] 90694309
[10] 90694310
[11] 90694311
[12] 90694312
[13] 90694313

Description

- Adapter for round bottom flask 50 ml
Adapter for round bottom flask 100 ml
Adapter for round bottom flask 250 ml
Adapter for round bottom flask 500 ml
Poly-bloc stand red, 11 holes \varnothing 12,5 mm / width 20 mm / tubes 4 ml
Poly-bloc stand purple, 4 holes \varnothing 28 mm / width 24 mm / tubos de 20 ml
Poly-bloc stand blue, 4 holes \varnothing 28 mm / width 30 mm / tubos de 30 ml
Poly-bloc stand black, 4 holes \varnothing 28 mm / width 43 mm / tubos de 40 ml
Poly-bloc stand green, 6 holes \varnothing 17,8 mm / width 26 mm / tubos de 8 ml
Poly-bloc stand gold, 4 holes \varnothing 21,6 mm / prof. 31,7 mm / tubos de 16 ml
Blue support for 4 poly-bloc stands
Blue ring for 4 poly-bloc stands
Protective cover temperature resistance 135°C

9- ASSOCIATED STANDARDS AND REGULATIONS

Construction in accordance with the following safety standards:

- EN 61010-1
- UL 3101-1
- CAN/CSA C22.2 (1010-1)
- EN61010-2-10

Construction in accordance with the following EMC standards:

- EN 61326-1

Associated EU guidelines:


- EMC guidelines: 89/336/EWG
- Machine guidelines: 73/023/EWG

10- TECHNICAL DATA

Reference	Model	Heating	Hotplate
50684010	RSLAB-1NC	NO	stainless steel
50684020	RSLAB-1NC	NO	porcelain enamel
50694010	RSLAB-1C	YES	stainless steel
50694020	RSLAB-1C	YES	porcelain enamel

Voltage:	200-240 VAC/ 100-120 VAC
Frequency:	50/60 Hz
Power:	130W / 530W (Heating model)
Stirring point position:	1
Max. stirring quantity (H₂O):	20 Liters
Max. magnetic bar:	80 mm
Motor type:	Brushless with external rotor
Max. power input of motor:	18 W
Max. power output of motor:	10 W
Speed range:	0-1500 rpm
Speed display accuracy:	1 rpm
Ø of the hotplate:	135 mm
Heating power:	500 W
Heating rage (1L water):	6 K/min
Temperature range:	RT-340 °C
Hotplate safety temperature range:	50-350 °C
Dimensions:	280x160x85 mm
Weight:	2,8 kg
Permitted ambient temperature:	5-40 °C
Permitted relative humidity:	80 %
Protection class acc. to DIN 60529:	IP42

- ◆ Observe the minimum distances between the devices, between the instruments and the wall and above the assembly (min. 100 mm).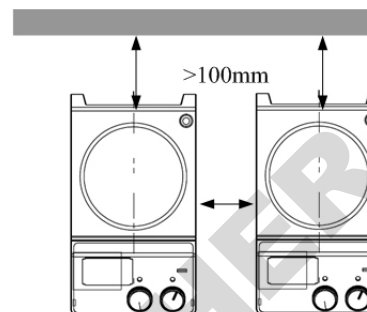

2- PROPER USE

The instrument is designed for schools, laboratories or factories. This device is not suitable for use in residential areas or other areas that may cause danger to the user or instruments as mentioned in chapter 1.

3- INSPECTION

- ◆ Unpack the equipment carefully and check for any damages which may have arisen during transport. If it happens, please contact manufacturer for technical support.
- ◆ If there is any apparent damage to the system, please do not plug it into the power line.
- ◆ The RSLAB-1NC y RSLAB-1C includes the following items:
 - Main unit: 1 unit
 - Power cable: 1 unit
 - Stirrer bar: 1 unit
 - User manual: 1 unit


4- TRIAL RUN

- ◆ Make sure the required operating voltage and power supply voltage match.
- ◆ Ensure the socket must be earthed.
- ◆ Ensure the power be off and the speed control knob and the temperature control knob to the lowest position.
- ◆ Plug in the power cable and power on the device.

- ◆ Add the medium into the vessel with a stirring bar.
- ◆ Put the vessel on the plate.
- ◆ Set the rated stirring speed and the device begins to work.
- ◆ Set the rated temperature and the device begins to work (heating model).
- ◆ Keep the temperature control knob and speed control knob slowly to the lowest position to turn off the function.
- ◆ If these operations are not normal, the device may be damaged during transportation, please contact your dealer for technical support.

5- CONTROL AND DISPLAY

1. LED heat display (heating model)
2. LED stir display
3. Heating knob (heating model)
4. Stirring knob
5. Switch ON/OFF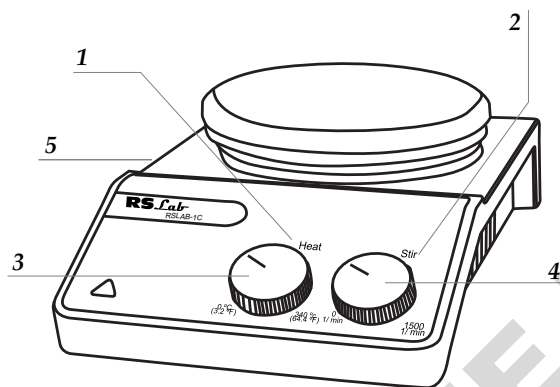

Items	Descriptions
Stirring Knob	Set the rated rotary speed in the safe stirring range from 0 to 1500 rpm. The function "stirring" is switched ON or OFF via the knob
Heating Knob (Heating model)	Set the rated temperature in the safe temperature range from room temperature to 340 °C. The function "heating" is switched ON or OFF via the knob
LED Heat (Heating model)	Lit when heating
LED Stir	Lit when stirring
Power Switch	Switch ON/OFF

6- OPERATION

- ◆ Put the device on the stable and safe place, ensure the speed control knob and the temperature control knob to the lowest position and plug in the mains power.
- ◆ Turn ON the power switch.
- ◆ Turn the speed control knob regulation to set the rated rotary speed in the safe speed limit from 0 to 1500 rpm.
- ◆ Turn the temperature control knob regulation to set the rated temperature in the safe limit from 0 to 340 °C. (Heating model)
- ◆ The instrument begins to work.

7- FAULTS

- ◆ If a unit fault happens, please power down the instrument and wait. Switch OFF the unit at the main ON/OFF switch for a few seconds.
- ◆ The stirring function will continue to operate at the speed set before the fault took place.
- ◆ The heating function will continue to operate at the set point before the fault took place.
- ◆ If the problem is not solved, take the unit to your technical support.

8- MAINTENANCE AND CLEANING

Proper maintenance can keep instruments working in a good state and lengthen its lifetime.

- ◆ Be careful not spray the cleanser into the instrument when cleaning.
- ◆ Unplug the power line when cleaning.
- ◆ Only use cleaner that we advised as below:
 - Dyes: isopropyl alcohol
 - Construction materials: water containing tenside / isopropyl alcohol
 - Cosmetics: water containing tenside / isopropyl alcohol
 - Foodstuffs: Water containing tenside.
 - Fuels: water containing tenside
- ◆ Before using other method for cleaning or decontamination, the user must ascertain with the manufacturer that this method does not harm or destroy instrument.
- ◆ Wear the proper protective gloves during cleaning of the instrument.
- ◆ The instrument must be cleaned and put it into the initial packaging carton before sending to service for repair, avoiding the contamination of hazardous.
- ◆ Use the instrument in a dry clean room and temperature stable environment.