**Bedienungsanleitung | Instructions for use |
Mode d'emploi | Instrucciones de uso | Istruzioni per l'uso**

Combisart®

Combisart® Vakuumfiltrations-Systeme zur Koloniezahl- und Partikelbestimmung |
Combisart® Vacuum Filtration Systems for Colony Count and Particle Counting |
Systèmes de filtration sous vide Combisart® pour la détermination et le
dénombrement des colonies et des particules | Sistemas de filtración al vacío Combisart®
para la determinación del número de colonias y partículas | Sistemi di filtrazione sotto
vuoto Combisart® per la conta delle colonie e l'analisi delle particelle

Deutsch	Seite 3
English	page 12
Français	page 21
Español	página 30
Italiano	pagina 39

DOMINIQUE DUTSCHER SAS

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Table of Contents

1	Use of Funnels as Combisart® System Components	13
1.1	Stainless Steel Systems 1-, 3- and 6-branch	13
1.1.1	Assembling Branch Systems	13
1.1.1.1	Biosart® 250 Funnels	13
1.1.1.2	Biosart® 100 Monitors	13
1.1.1.3	Stainless Steel Funnels	13
1.1.1.4	Glass Funnels	13
1.1.1.5	Polycarbonate Filter Holder	14
1.2	Individual Filter Holder System	14
1.2.1	Assembling Individual Filter Holder Systems	14
1.2.1.1	Biosart® 250 Funnels	14
1.2.1.2	Biosart® 100 Monitors	14
1.2.1.3	Stainless Steel Funnels	14
1.2.1.4	Glass Funnels	14
2	Assembling the Complete Vacuum Filtration System	15
2.1	Required Accessories	15
3	Microbiological Tests	16
4	Analyses	17
5	Cleaning and Care of Stainless Steel Filter Holders	17
6	Maintenance of Quick-Lock Clamps on Stainless Steel Funnels	17
7	Maintenance of Stainless Steel Three-Way Valves	18
8	Spare Parts	18
8.1	Stainless Steel Manifolds, 1-, 3- and 6-Branch	18
8.2	Individual Bases	18
8.3	Adapters	18
9	Specifications	19
9.1	Stainless Steel System Components	19
9.2	Connectors and Adapters	19
9.3	Biosart® 100 Monitors and the Biosart® 250 Funnel	19
9.4	Glass and Polycarbonate Filter Holders	19
10	Figures	
10.1	Three-Way Valve Positions	20
10.2	Multiple-Branch Systems	48
10.3	Individual Filter Holder Systems	48

Sartorius Stedim Biotech vacuum filtration systems have been specifically developed for routine residue analysis, particularly colony and particle counting.

The filter supports are designed to ensure uniform distribution of substances retained (bacteria, particles) on the surface of the membrane filter. Releasing the vacuum under sterile conditions by attaching a special air filter to the three-way valves, prevents secondary contamination on the underside of the filter. In addition, the stainless steel valves allow easy control of the vacuum supply. The special quick-lock clamp simplifies the insertion and removal of the membrane filters and guarantees secure attachment.

More information is available on the internet under www.sartorius-stedim.com/microbio in the section entitled **Stainless Steel Vacuum Systems|Combisart® Hardware Setups**.

1. Use of Funnels as Combisart® System Components

Combisart® is a modular-designed system specifically developed for residue analysis. Depending on the application, the corresponding funnel can be adapted to each base, which is connected to a vacuum source.

Multiple-branch systems with 3 or 6 filter units are employed for extensive sampling. Individual filter holder systems with just one filter unit are suitable when few filtration runs are carried out.

1.1. Stainless Steel Systems, 1-, 3- and 6-branch

Please refer to the last fold-out page for more detailed information on these systems.

1.1.1 Assembling Multiple-Branch Systems

- Insert the L-shaped glass tube into the stopper of the 5-liter filtrate flask (16672-1), then insert the stopper into the opening of the filtrate flask.
- Connect the outlet of the stainless steel manifold (hose connector) to the inlet of the filtrate flask (glass tube) using the vacuum hose.
- Attach a Minisart® SRP 25 (17575 ACK) to the side opening of the three-way valve for each filter unit.

1.1.1.1 Biosart® 250 Funnels, 16407

- Screw the single 16840 base into the filter unit thread on the manifold, turning the base until the two pins are positioned either right | left or front | back.
- Tighten the threaded nut using a 24-mm open-end wrench (spanner).
- Insert the flat silicone gasket into the single base, and place the stainless steel filter support onto the silicone gasket.
- Remove a membrane filter from the package, and center it on the stainless steel filter support. For further instructions on how to perform the analysis, please refer to Section 3 "Microbiology" and | or Section 4 "Analyses."
- Place the Biosart® 250 Funnel onto the base so that it fits precisely, and clamp down the membrane filter simply by turning the funnel clockwise.

1.1.1.2 Biosart® 100 Monitors 16401, 16402, 16403

a) Contamination-free sealing above the edge of the Monitors

- Screw the single 16840 base support into the filter unit thread on the manifold.
- Tighten the threaded nut using a 24-mm open-end wrench (spanner).
- Insert the flat silicone gasket into the single base, and place the stainless steel filter support onto the silicone gasket.
- Place the Biosart® 100 adapter (16414) onto the base support so that the pins on the support fit into the two notches on the adapter.
- You can now easily and safely attach and remove the Monitors individually. The 47-mm Monitors provide a seal via the internal groove. This unique design guarantees uniform distribution of the residue on the membrane surface when the Monitor is positioned so that it is perfectly level.

b) Sealing via the filtrate outlet of the Monitor

- Make certain that the blue silicone gasket is seated properly in the groove on the upstream side of the filter unit.
- Attach the 16835 adapter to the filter unit thread on the manifold, forming a tight seal.
- Insert the outlet of each Monitor into the stopper of the adapter. You can now easily and safely attach and remove the Monitors individually.

1.1.1.3 Stainless Steel Funnels, 40, 100 and 500 ml

- Screw the single 16840 base into the filter unit thread on the manifold until the two pins are positioned either right | left or front | back.
- Tighten the threaded nut using a 24-mm open-end wrench (spanner).
- Insert the flat silicone gasket into the base, and place the stainless steel filter support onto the silicone gasket.
- Remove a membrane filter from the package, and centre it on the stainless steel filter support. For further instructions on how to perform the analysis, please refer to Section 3 "Microbiology" and | or Section 4 "Analyses."
- Place the stainless steel funnel onto the base so that it fits precisely and lock the funnel in place with the clamp (just lightly push down on the quick-lock clamp to secure the parts; there is no need to force the clamp down to its full extent).
- If you are using a lid to prevent airborne contamination, place the silicone seal into the lid first, then close the stainless steel funnel with the lid (does not apply to the 40-ml funnel).

1.1.1.4 Glass Funnels, 16306 | 16307

- Make certain that the blue silicone gasket is seated properly in the groove on the upstream side of the filter unit.
- Attach adapter no. 16836 (for the 16306 glass funnel) or no. 16837 (for the 16307 glass funnel) to the filter unit thread on the manifold, forming a tight seal.
- Insert the base of the glass funnel into the stopper of the adapter, and attach the white filter support on top of the base.
- Remove a membrane filter from the package, and center it on the filter support. For further instructions on how to perform the analysis, please refer to Section 3 "Microbiology" and | or Section 4 "Analyses."
- You can now easily and securely connect the upper part and base of the glass funnel using the metal clamp.

1.1.1.5 Polycarbonate Filter Holder, 16511

- Make certain that the blue silicone gasket is seated properly in the groove on the upstream side of the filter unit.
- Screw the polycarbonate filter holder into the filter unit thread on the manifold, forming a tight seal.
- For further instructions on how to perform the analysis, please refer to the Directions for Use accompanying the filter holder.

1.2 Individual Filter Holder Systems

For individual systems the filter station is positioned on the top of a suction flask. Please refer to the last fold out page for more detailed information on these systems.

1.2.1 Assembling the Individual Filter Holder Systems

There are two different kinds of filtrate flasks available for the individual filter holders: a 2-liter flask, (no. 16672), with a plastic safety connector, and a 1-liter flask, (no. 16606), with a glass connector. Refer to the following funnel descriptions to obtain the order numbers for the silicone stoppers (in parentheses) which fit the 16672 2-liter flask (and the 16606 1-liter flask).

1.2.1.1 Biosart® 250 Funnels, 16407

- Attach the 16841 stainless steel individual base to the 17173 (17004) silicone stopper and insert the stopper into the opening of the filtrate flask. By turning the individual base, position the two pins so that they are either right|left or front|back.
- Insert the flat silicone gasket into the individual base, and place the stainless steel filter support onto the silicone gasket.
- Attach a Minisart® SRP 25 (17575 ACK) to the side opening of the three-way valve.
- Remove a membrane filter from the package, and center it on the stainless steel filter support. For further instructions on how to perform the analysis, please refer to Section 3 “Microbiology” and|or Section 4 “Analyses.”
- Place the Biosart® 250 funnel onto the base so that it fits precisely, and clamp down the membrane filter by simply turning the funnel clockwise.

1.2.1.2 Biosart® 100 Monitors, 16401, 16402, 16403

- Insert the 16841 stainless steel individual base into the 17173 (17004) silicone stopper and insert the stopper into the opening of the filtrate flask.
- Insert the flat silicone gasket into the individual base, and place the stainless steel filter support onto the silicone gasket.
- Attach a Minisart® SRP 25 (17575 ACK) to the side opening on the three-way valve.
- Place the Biosart® 100 Adapter (16414) onto the 16841 stainless steel individual base so that the pins on the support fit into the two notches on the adapter.
- You can now easily and safely attach and|or remove the Monitors individually. The 47-mm Monitors provide a seal via the internal groove. This unique design is to guarantee uniform distribution of the residue on the membrane surface when the Monitor is positioned absolutely level.

1.2.1.3 Stainless Steel Funnels, 40, 100 and 500 ml

- Insert the 16841 individual stainless steel base into the 17173 (17004) silicone stopper and insert the stopper into the opening of the filtrate flask. By turning the individual base, position the two pins so that they are either right|left or front|back.
- Insert the flat silicone gasket into the individual base, and place the stainless steel filter support onto the silicone gasket.
- Attach a Minisart® SRP 25 (17575 ACK) to the side opening on the three-way valve.
- Remove a membrane filter from the package, and centre it on the stainless steel filter support. For further instructions on how to perform the analysis, please refer to Section 3 “Microbiology” and|or Section 4 “Analyses.”
- Place the stainless steel funnel onto the base so that it fits precisely, and lock the funnel in place using the clamp (lightly push down on the quick-lock clamp to secure the parts; there is no need to force the clamp down to its full extent).
- If you are using a lid to prevent airborne contamination, place the silicone seal into the lid first, then close the stainless steel funnel with the lid (does not apply to the 40-ml funnel).

1.2.1.4 Glass Funnels, 16306|16307

- Insert the base of the glass funnel into the 17174 (17005) or 17175 (17006) silicone stopper and insert the stopper into the opening of the filtrate flask.
- Seat the white filter support into the base.
- Remove a membrane filter from the package, and center it on the filter support. For further instructions on how to perform the analysis, please refer to Section 3 “Microbiology” and|or Section 4 “Analyses.”
- You can now easily and safely connect the upper part and base of the glass funnel using the metal clamp.

2. Assembling the Entire Vacuum Filtration System

Sartorius Stedim Biotech vacuum filtration systems have been specially matched to one another. Problems may be encountered if components of products from other manufacturers are used with Sartorius Stedim Biotech systems. Applications technology specialists have performed numerous tests to select ideal components and materials which guarantee reliable vacuum filtration.

The manufacturer's warranty will apply only if the specified procedure is followed before, during and after vacuum filtration.

2.1 Required Accessories:

Protection against contamination:

Lid for stainless steel funnel, 100 ml	6981063
Lid seal, 100 ml	6981064
Lid for stainless steel funnel, 500 ml	6981001
Lid seal, 500 ml	6981003

Vacuum sources:

Vacuum pump for multiple-branch systems, 220 V, 50 Hz	16694-2-50-22
Vacuum pump for multiple-branch systems, 110 V, 60 Hz	16694-1-60-22
Vacuum pump for individual filter holder systems, 220 V, 50 Hz	16694-2-50-06
Vacuum pump for individual filter holder systems, 110 V, 60 Hz	16694-1-60-06
Vacuum hose	16623

Protection of vacuum pumps:

Woulff's bottle	16610
Vacusart®	17804M

Handling the membrane filters:

Stainless steel blunt-tipped forceps	16625
--------------------------------------	-------

Microbiological testing:

Laboratory incubator, 220 V, 50/60 Hz	18113
Colony counter	17649

Preparation of nutrient pads:

Dosing syringe	16685-2
Minisart®, 0.2 µm	17597K

Sterile release of vacuum:

Minisart® SRP 25, 0.2 µm	17575-ACK
--------------------------	-----------

Removal of coarse particles:

Prefilter attachment	16807
Cellulose acetate prefilter, 8 µm	12301-050-ACN

- Use the vacuum hose to connect the filtrate flask to the water trap or barrier (Woulff's bottle or Vacusart®). Before connecting the vacuum hose, attach the safety connector to the side outlet of the 5- and/or 2-liter filtrate flask (does not apply to the 16606 1-liter filtrate flask).
- Connect the vacuum pump to the water trap. When installing the Vacusart®, make sure that the filtrate overflow spills back into the suction flask and is not left to stand.
- Place the stopper onto the filtrate flask, and connect the system components as shown in the drawings and descriptions under 1.1 "Multiple-Branch Systems" and/or 1.2 "Individual Filter Holder Systems."

3. Microbiological Tests

Below is a description of the methodology to determine colony counts with reusable systems made of stainless steel, glass or polycarbonate. A detailed description of microbiological test methods used for ready-to-connect units, such as Biosart® 100 Monitors or 250 Funnels, can be found in the Instructions for use accompanying the corresponding product.

3.1 All system components that come into contact with the sample and filter membrane must be sterile. Sterility can be ensured in different ways depending on the materials utilized.

a) **Autoclaving** (for stainless steel, glass and polycarbonate): Assemble the unit so that the filter membrane is clamped securely between the base and the funnel. The filter should not be clamped too tightly; otherwise, steam will not penetrate the clamped region of the membrane.

Adjust the three-way valve to the 12 o'clock position

(does not apply to individual glass filter holders). It is not imperative to sterilize the 16844, 16842 and/or 16843 stainless steel manifolds when the adapted filter holders (Biosart® 100 or 250, stainless steel, glass, polycarbonate) are changed under sterile conditions.

Cover the inlets and outlets with aluminum foil, and autoclave according to the specifications of the system components. After sterilization, assemble the filtration funnel, ascertaining all connections are tight, and attach the funnel to other system components.

Adjust the three-way valve to the 9 o'clock position.

b) **Flaming** (only for stainless steel): Flaming prevents long waiting periods (autoclave cycles) between filtration runs and is described in detail below.

3.2 Clean the laboratory bench, and disinfect with alcohol or a commercial disinfectant.

3.3 Line up your sterile culture media, such as nutrient pads in petri dishes or agar in petri dishes. Moisten the nutrient pads with 3.0–3.5 ml of sterile distilled or demineralized water. The dosing syringe (16685-2) with a sterile Minisart® (17597-K) attached simplifies the addition of water to nutrient pads.

3.4 Place a Bunsen burner with a glass beaker containing a little alcohol and a pair of blunt-tipped forceps next to the autoclaved filter holder. If you are working with a stainless steel filter holder and will sterilize it by flaming, proceed as follows:

a) Locate the sterile packaged membrane filter within easy reach.

b) Turn on the vacuum source and remove the stainless steel funnel.

c) Using a Bunsen burner flame the filter support including the stainless steel frit by opening the valve (turn to 6 o'clock) to draw the flame through the filter support. Then close the valve (turn to 9 o'clock).

d) Hold the stainless steel funnel in one hand and flame it from below.

e) Reposition the stainless steel funnel on the base, and lightly push down on the quick-lock clamp to secure the funnel (there is no need to force the clamp down to its full extent). Using a spiral pattern, flame the funnel on the inside from the bottom to the top.

f) Flame the lid on the inside, and place it back on the stainless steel funnel.

g) To accelerate cooling of the filter holder and simplify insertion of the membrane filter, rinse the filter holder with sterile water (e.g. using a dosing syringe with a sterile Minisart® attached). To remove the water via vacuum, open the valve (turn to 6 o'clock), then close it (turn to 9 o'clock).

h) Sterilize the blunt-tipped forceps briefly in the flame and remove a sterile membrane filter (if necessary with the yellow protective paper) from the package. Using one hand, remove the stainless steel funnel with the lid from the base of the filter holder, and use your other hand to center the membrane filter (yellow protective paper or grid should face upwards) on the filter support.

i) If there is a yellow protective disc, make sure to discard it before assembling the funnel or the top part of the filter holder. Lock the funnel in place using the clamp (lightly push down on the quick-lock clamp to secure the parts; there is no need to force the clamp down to its full extent).

3.5 Pour the sample to be tested into the funnel. For samples less than 5 ml, use 10 to 20 ml of sterile water or physiological saline as diluent. Re-open the valve (turn to 6 o'clock) to suction off the liquid, then close the valve (turn to 9 o'clock).

3.6 Remove the funnel and using the flamed, blunt-tipped forceps, remove the membrane filter.

3.7 Place the membrane filter on a nutrient pad or an agar plate and incubate. Make certain no air bubbles are entrapped between the membrane and nutrient pad.

3.8 The filter holders and components should be rinsed them with 70% alcohol and air-dry for storage.

3.9 The grid printed on the membrane facilitates evaluation of the colony count. The grid optically divides the filtration surface into 130 squares, each measuring 3.1 × 3.1 mm. Therefore, one square equals 1/130 of the overall area (if you use the 6980103 stainless steel filter support with a diameter of 47 mm, available as an accessory, the filtration area is reduced to 9.6 cm²; hence, one grid square equals 1/100 of the 9.6 cm²).

Important Note: Prefiltration

Samples with a high particle load of dirt or insoluble sediment must be prefiltered for bacteriological testing. The stainless steel prefilter attachment (16807) (which also can be flamed) allows you to perform prefiltration concurrently with membrane filtration. Attach it between the stainless steel funnel and the base, and fill completely with sterile water (to prevent air from becoming entrapped and interfering with filtration), and insert a 11301-050 bacteriological prefilter (8 µm pore size). This prefilter retains any coarse particles in the sample, reducing the particulate load on the membrane filter on the stainless steel filter support. Colonies are able to grow unimpeded by interfering deposits. The prefilter attachment can also be flamed.

4. Analyses

4.1 Assemble the filtration system as described in Sections 2 and 3. A Minisart® SRP25 should be used as a particle trap during filtration to prevent exposure of the membrane underside and filter support to secondary contamination with particles. If particle contamination is irrelevant for your application, close off the Minisart® outlet with a luer lock (male) plug.

4.2 If applicable, rinse any particles off the filter holder, the membrane filter and the blunt-tipped forceps.

4.3 Remove the corresponding funnel from the base of the filter holder, and, using the blunt-tipped forceps, place the membrane filter on the filter support.

4.4 Place the funnel back onto the base of the filter holder, and lock the funnel in place with the clamp (lightly push down on the quick-lock clamp to secure the parts; there is no need to force the clamp down to its full extent). Biosart® 250 funnels can be locked in place by turning clockwise.

4.5 Pour the sample to be tested into the funnel.

4.6 Turn on the vacuum source and open the three-way valve (turn to 6 o'clock).

4.7 At completion of filtration, close the valve (turn to 9 o'clock) or turn off the vacuum pump. Remove the funnel and transfer the membrane filter for gravimetric, optical, chemical or other analysis.

4.8 The grid printed on the membrane facilitates visual evaluation of the residue. The grid optically divides the filtration surface into 130 squares, each measuring 3.1×3.1 mm. Therefore, one square equals $1/130$ of the overall area (if you use the 6980103 stainless steel filter support with a diameter of 47 mm, available as an accessory, the filtration area is reduced to 9.6 cm^2 ; hence, one grid square equals $1/100$ of the 9.6 cm^2).

5. Cleaning and Care of Stainless Steel Filter Holders

To ensure a long service life, stainless steel filter holders should be cleaned prior to initial use and after each filtration run.

Proceed as follows:

5.1 Disassemble the filter holder.

5.2 Use soft brushes, warm water and a gentle, commercially available laboratory detergent (for metal, glass, plastic) to clean all parts.

5.3 Rinse the parts with hot water and then with distilled water. Drying can be accelerated by rinsing with a 70% alcohol solution.

5.4 Dry the parts in a drying oven. Please do not use towels or cloth for drying because they may deposit lint on the filter holders.

6. Maintenance of Quick-Lock Clamps on Stainless Steel Funnels

The quick-lock clamps tend to widen after frequent use, and must be tightened to prevent leakage. Grasp the clamp with a pair of flat-nose pliers and apply slight pressure to narrow the clamp's opening.

Important Note:

Lightly push down on the quick-lock clamp to secure the filter holder parts; there is no need to force the clamp down to its full extent.

7. Maintenance of Stainless Steel Three-Way Valves

If it becomes difficult to open and close any of the valves, the valves should be removed for cleaning, drying and lubrication (e.g. generously lubricate with high-grade heavy grease). To remove a valve, proceed as follows:

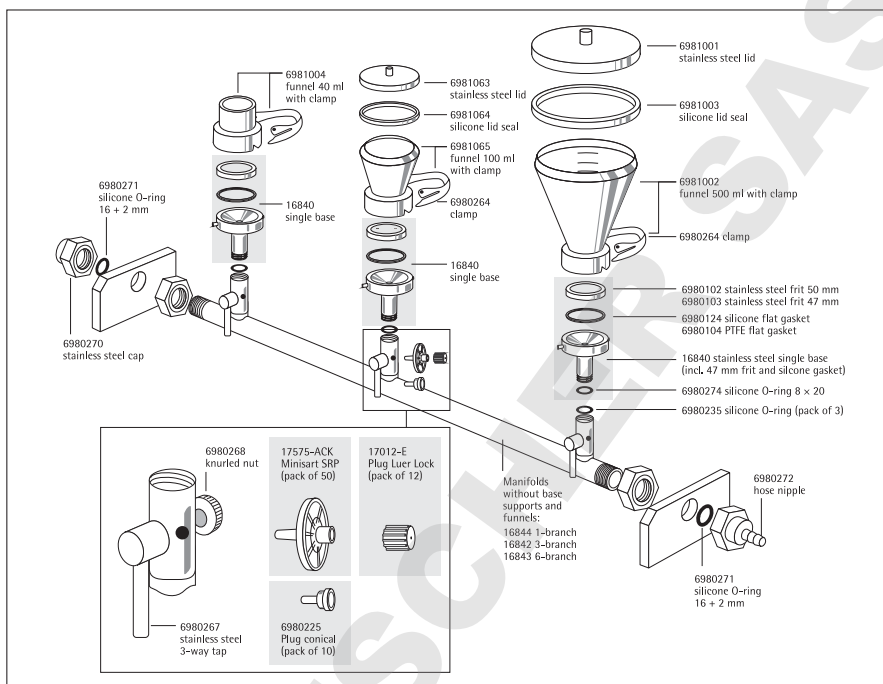
7.1 First, remove the knurled nut and the washer from the back of the valve.

7.2 Then detach the valve.


If valves from several filtration filter holders have been disassembled, then each valve must be replaced in the same tapered stem guide from which it was removed. The valve stem guide has been precisely machined to the valve's individual specifications to make a leak-tight fit.

8. Spare Parts

8.1 Spare parts for Stainless Steel Manifolds

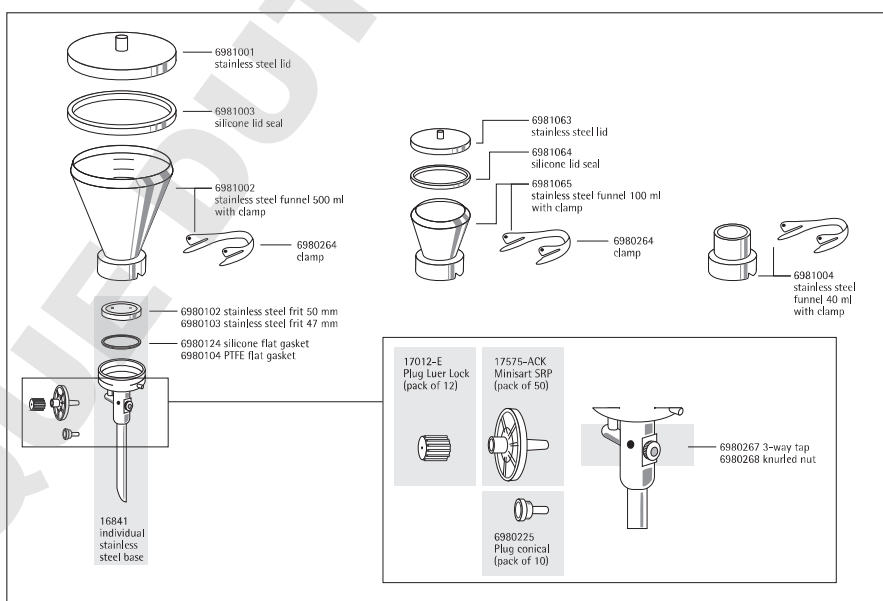8.2 Spare Parts for Individual Bases

8.3 Spare Parts for the 16835 | 16837 Adapters

Silicone stoppers for 16835	00279
Silicone stoppers for 16836	00280
Silicone stoppers for 16837	00281

Spare Parts for the 16306 | 16307 Glass Filter Holders

Please refer to the Directions for Use (FL-6001-a) accompanying the product for further information.

Spare Parts for the 16510 Polycarbonate Filter Holder

Please refer to the Directions for Use (FL-6026-n) accompanying the product for further information.

9. Specifications

9.1 Stainless Steel System Components

	Manifolds 16844 16842 16843	Funnel Lid Seal 6981004 6981065 063 064, 6981002 001 003	Base Support 16840 Single Base 16841
Material	AISI 304 stainless steel (B.S. 304S31/German std. 1.4301)	AISI 304 stainless steel (B.S. 304S31/German std. 1.4301)	AISI 304 stainless steel (B.S. 304S31/German std. 1.4301)
Gasket	Silicone, 14×2 mm (3 units)	Silicone, 77.2×85.8 (6981064) Silicone, 122×131 (6981003)	Silicone, 41×50×1
Filter diameter	–	47 mm (or 50 mm)	47 mm (or 50 mm)
Filtration area	–	12.5 cm ²	12.5 cm ²
Dimensions (L/H/D)	16842: 435/103/120 16843: 910/103/120	–	–
Maximum pressure limit	Vacuum, max. 2 bar pressure limit	Vacuum only	Vacuum only
Disinfection	–	By flaming	By flaming
Sterilization			
Autoclaving	134°C max., 30 min.	134°C max., 30 min.	134° max. C, 30 min.
Dry heat	180°C max., 30 min.	180°C max., 30 min.	180°C max., 30 min.
Inlet	TR 20×2 female thread	–	–
Outlet	Hose connector, DN 10, 10 mm	–	16840: TR 20×2 mm male thread with DN 24 hex nut 16841: 10 mm outlet connectors

9.2 Connectors and Adapters

	Filter Holder Adapters 16835 16836 16837	Biosart® 100 Adapters 16414 16415 16416
Material	AISI 304 stainless steel (B.S. 304S31/German std. 1.4301)	Silicone (16414, 16415) Polypropylene (16415)
Stoppers	Silicone	–
Max. pressure limit	Vacuum only	Vacuum only
Sterilization		
Autoclaving	134°C max., 30 min.	134°C max., 30 min.
Dry heat	180°C max., 30 min.	–
Inlet	16835: 10 mm drill hole 16836: 11 mm drill hole 16837: 14 mm drill hole	–
Outlet	TR 20×2 mm male thread	–

9.3 Biosart® 100 Monitors and the Biosart® 250 Funnel

	Biosart® 100 Monitors 16401 02 03, 47 mm	Biosart® 250 Funnel 16407
Materials:		
Housing	Polystyrene	Polypropylene
Membrane filter	Cellulose nitrate with grid	–
Pad	Cellulose	–
Plug adapter	Polyethylene	–
Filter diameter	47 mm	47 mm (or 50 mm)
Filtration area	15.5 cm ²	12.5 cm ²
Capacity	100 ml, graduated in 10-ml increments	250 ml, graduated in 50-ml increments
Max. pressure limit	Vacuum only	Vacuum only
Package quantity	50 units, sterile	50 units, sterile
Outlet	6.5×1.5 mm	–

9.4 Glass Filter Holders and the Polycarbonate Filter Holder

	Glass Filter Holder, 16306	Glass Filter Holder, 16307	Polycarbonate Filter Holder, 16511
Materials:			
Adapter base	Borosilicate glass 3.3	Borosilicate glass 3.3	Polycarbonate
Clamp	Anodized aluminum	Anodized aluminum	–
Filter support	PTFE borosilicate glass 3.3	PTFE borosilicate glass 3.3	Polypropylene
Lid	–	Silicone	Polycarbonate
Seal	Silicone, 25×3 mm	Silicone, 45×3 mm	Silicone, 40×5, 80×3, 14×2
Filter diameter	25 mm, prefilter 20 mm	47 mm, prefilter 40 mm	47 mm, prefilter 37 mm
Filtration area	3 cm ²	12.5 cm ²	11.5 cm ²
Capacity	30 ml	250 ml	250 ml
Max. pressure limit	Vacuum only	Vacuum only	Vacuum, max. 2 bar pressure limit
Sterilization			
Autoclaving	134°C max., 30 min.	134°C max., 30 min.	121°C max., 30 min.
Dry heat	180°C max., 30 min.	180°C max., 30 min.	–
Outlet	12 mm outlet connector	15 mm outlet connector	TR 20×2 mm male thread

10. Figures

10.1 Three-Way Valve Positions

The Combisart® vacuum filtration system features a high degree of flexibility and versatility.

Sartorius Stedim Biotech is the first company to offer a system where all funnel types can be adapted to all bases.

Due to the unique modular system of Combisart®, preparation and filtration time can be reduced by changing the funnel material or number of units. Special attention has been paid to the handling of multiple-branch systems (3- and | or 6-place manifolds):

- Each filter unit can be autoclaved separately from the manifold, saving space inside autoclaves.
- By removing the filter unit from the manifold, the supernatant above a blocked membrane can be easily poured off.
- The funnels can be conveniently positioned for left or right-handed use.

The Minisart® SRPs, (attached to the side opening of the three-way valves), sterile vents the air from the underside of the filter, protecting it from any contamination by microorganisms and for particles from the surrounding environment, thus ensuring reliable results.

Valve Position
OPEN
6 o'clock

Valve Position
VENTING
9 o'clock

Valve Position
CLOSE
3 o'clock

Valve Position
AUTOCLAVE
12 o'clock

For Filtration

The full vacuum draws the sample through the membrane filter. The vent filter is "off-line."

After Filtration

The vacuum between the valve and membrane filter is released under sterile conditions, without any particles being drawn into the system. Secondary contamination of the bottom of the filter is ruled out entirely.

After the Filtration Run

The residual vacuum between the pump and valve is released via the sterilizing-grade filter.

For Autoclaving

For reliable sterilization, the steam flows freely through all openings.

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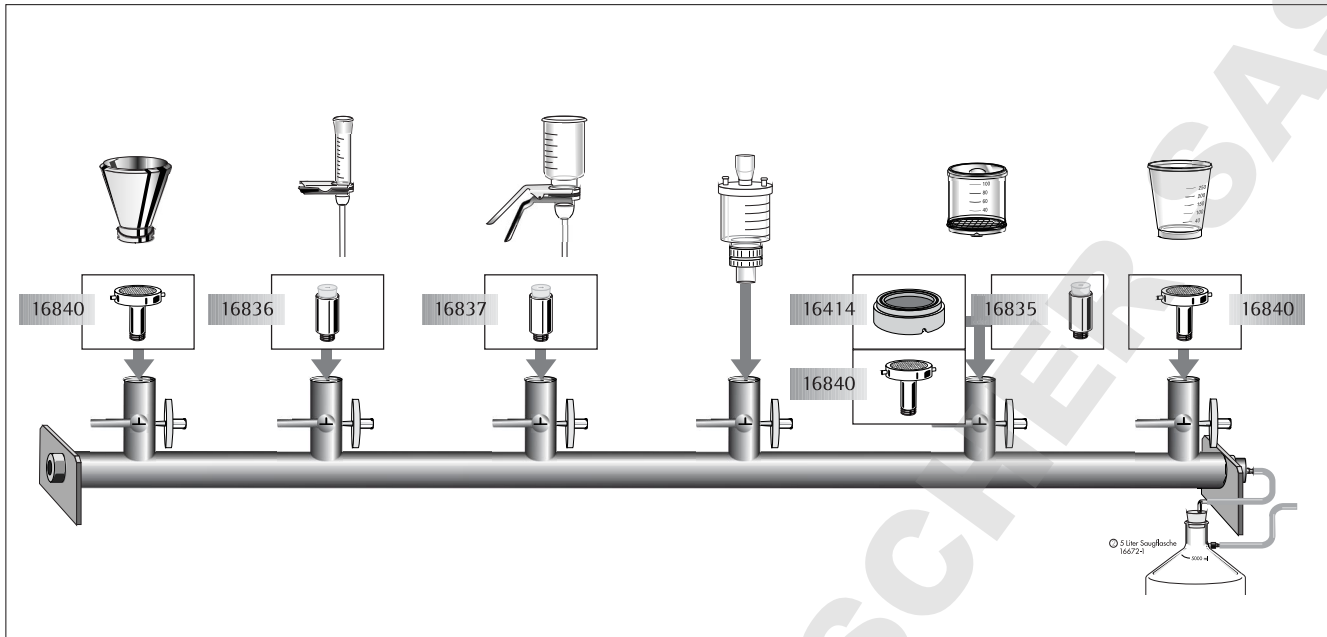
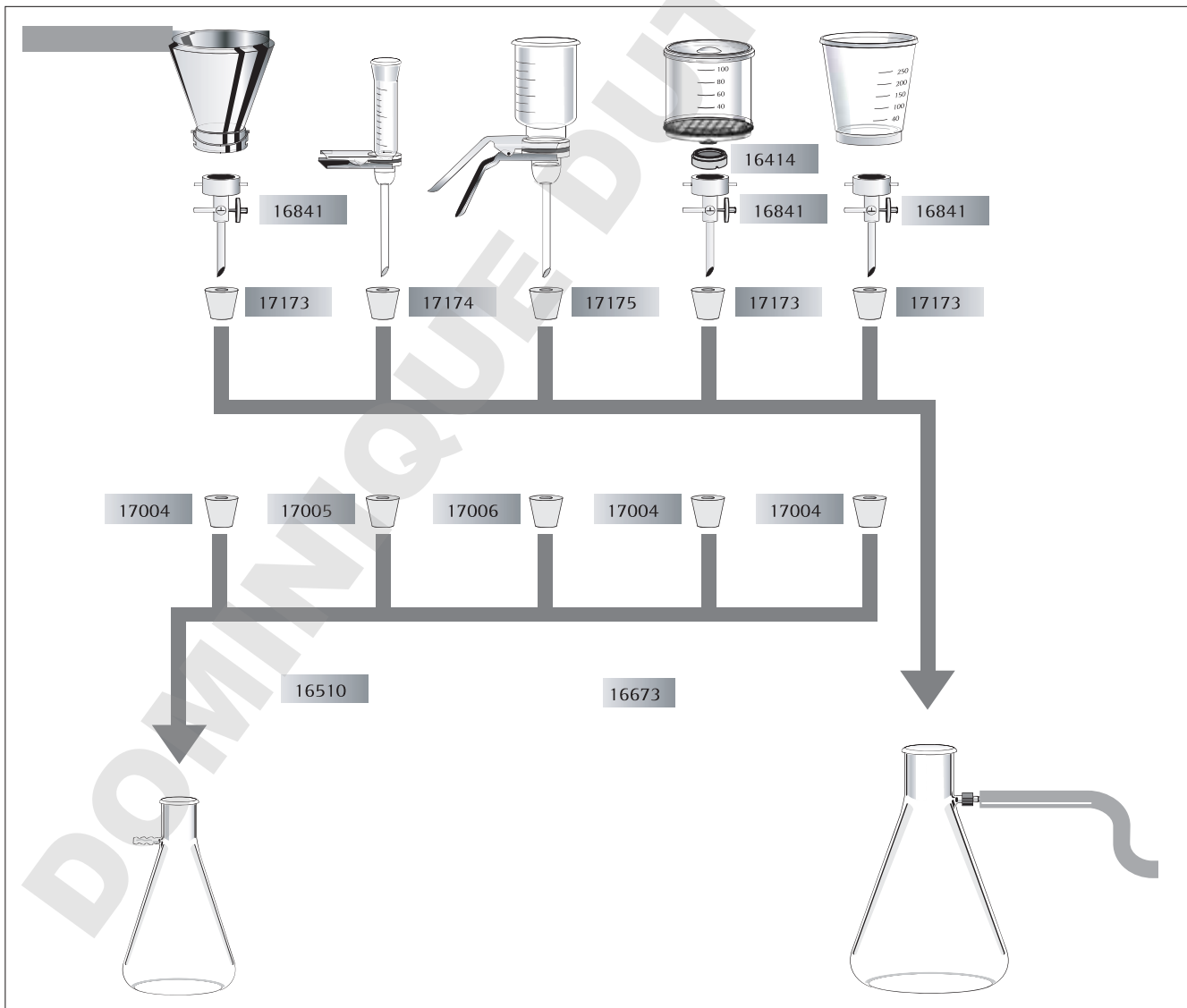
Sales and Service Contacts

For further contacts, visit www.sartorius-stedim.com

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Phone +41.44.741.05.00
Fax +41.44.741.05.09

U.K.

Sartorius Stedim UK Limited
Longmead Business Park
Blenheim Road, Epsom
Surrey KT19 9 QQ

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Asia | Pacific

China

Sartorius Stedim Beijing
Representative Office
No. 33, Yu'an Road,
Airport Industrial Zone B, Shunyi District
Beijing 101300

Phone +86.10.80426516
Fax +86.10.80426580

Sartorius Stedim Shanghai
Representative Office
Room 618, Tower 1, German Centre,
Shanghai, PRC., 201203

Phone +86.21.28986393
Fax +86.21.28986392.11

Sartorius Stedim Guangzhou Office
Room 704, Broadway Plaza,
No. 233-234 Dong Feng West Road
Guangzhou 510180

Phone +86.20.8351.7921
Fax +86.20.8351.7931

India

Sartorius Stedim India Pvt. Ltd.
10, 6th Main, 3rd Phase Peenya
KIADB Industrial Area
Bangalore – 560 058

Phone +91.80.2839.1963|0461
Fax +91.80.2839.8262

Japan

Sartorius Stedim Japan K.K.
KY Building, 8-11
Kita Shinagawa 1-chome
Shinagawa-ku
Tokyo 140-0001

Phone +81.3.3740.5407
Fax +81.3.3740.5406

Malaysia

Sartorius Stedim Malaysia Sdn. Bhd.
Lot L3-E-3B, Enterprise 4
Technology Park Malaysia
Bukit Jalil
57000 Kuala Lumpur

Phone +60.3.8996.0622
Fax +60.3.8996.0755

Singapore

Sartorius Stedim Singapore Pte. Ltd.
10, Science Park Road, The Alpha
#02-25, Singapore Science Park 2
Singapore 117684

Phone +65.6872.3966
Fax +65.6778.2494

Australia

Sartorius Stedim Australia Pty. Ltd.
Unit 17/104 Ferntree Gully Road
Waverley Business Park
East Oakleigh, Victoria 3166

Phone +61.3.9590.8800
Fax +61.3.9590.8828

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen, Germany

Phone +49.551.308.0
Fax +49.551.308.3289
www.sartorius-stedim.com

Copyright by Sartorius Stedim Biotech GmbH, Goettingen, Germany. All rights reserved. No part of this publication may be reprinted or translated in any form or by any means without the prior written permission of Sartorius Stedim Biotech GmbH. The status of the information, specifications and illustrations in this manual is indicated by the date given below. Sartorius Stedim Biotech GmbH reserves the right to make changes to the technology, features, specifications and design of the equipment without notice.

Status:
August 2008,
Sartorius Stedim Biotech GmbH,
Goettingen, Germany

Note: All product names with the syllable "sart" are either registered brand names or trade marks of Sartorius Stedim Biotech GmbH, Goettingen or for which an application for registration has already been submitted.

Combisart is a brand name of Sartorius Stedim Biotech GmbH, Goettingen, Germany and has been submitted for registration as a trade name.