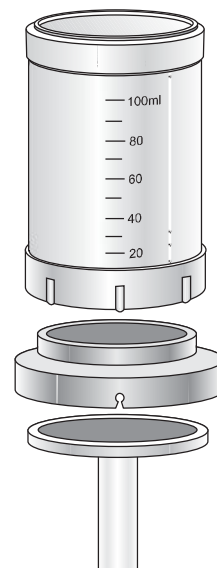
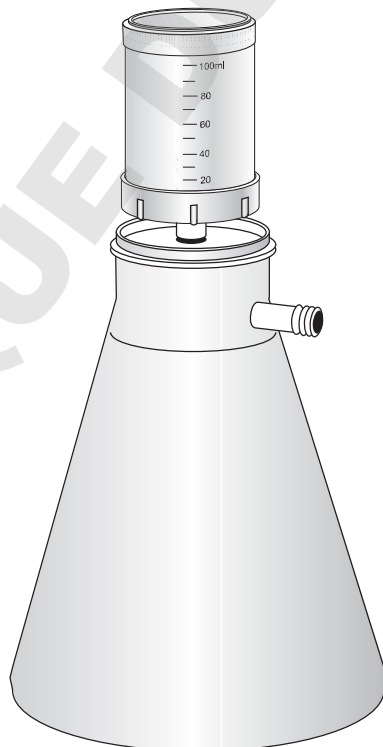
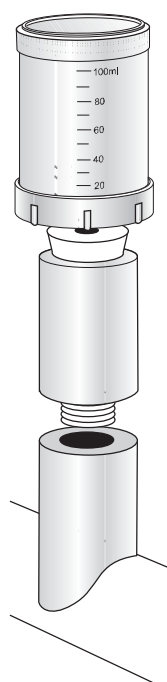
sartorius

Biosart® 100 Monitor

Instruction Manual

DOMINIQUE MUTSCHER SAS

(GB)

Biosart® 100 Monitor – Instructions for use Sartorius 100 ml Microbiological Monitor
Sterile, disposable 100 ml Microbiological Monitors with a 47 mm diameter membrane and large bottom port provides a faster flow rate and decreases testing time. Ideal for the detection of microorganisms in beverages, pharmaceuticals and water.

Materials included:

- 48 – Sterile, disposable microbiological monitors
- 48 – Bottom port plugs
- 1 – Large funnel adapter

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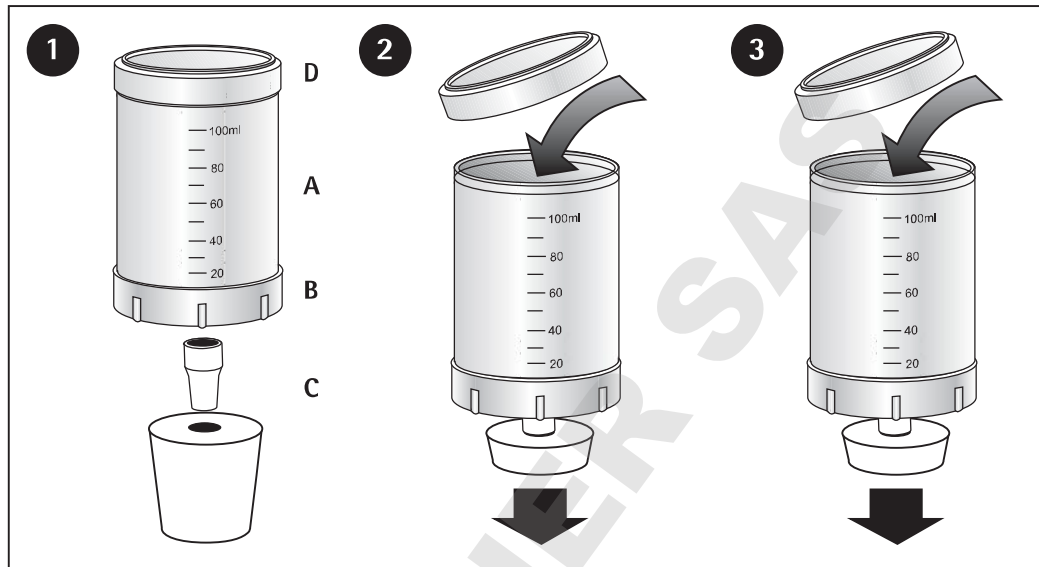
(GB)

1. Remove monitor from carton. Attach funnel adapter **C** to monitor base **B**. Securely insert the assembled unit to vacuum source.

Alternatively you can use the adapter 16424 to adapt the Biosart® 100 Monitor on the Combisart® manifold. Please leave the frit into the base and fix the adapter via bayonet on the Combisart base.

2. Remove the lid **D**, add sample and replace the lid. Attention: if there is only a small volume to filter, pour sterile buffer solution (10–20 ml) into the monitor before pouring the sample. Apply vacuum and filter the solution.

3. Turn vacuum off and rinse the inside walls of the funnel with sterile buffer solution (20–30 ml). Apply vacuum. When filtration is complete, turn vacuum off. Release vacuum from the unit with either the vacuum release cap on a suction manifold or with a three-way stopcock when utilizing sidearm vacuum flask.

4. Remove the lid of the funnel and add your selected media ampoule to the membrane and pad from the top. Be careful not to touch the tip of the media ampoule to the membrane. Replace lid. Apply a one-second vacuum to draw the media down into the pad.

5. Holding funnel **A** firmly, twist off the monitor base **B**. Discard the funnel **A**, retain the funnel lid **D** and funnel adapter **C**.

6. Push the funnel lid **D** on to the monitor base **B**. Remove the funnel adapter **C** from the monitor and replace with a port cap **F**. Invert the monitor and incubate at appropriate temperature for selected media. Ensure that the horizontal position of the petri dish is exact so there is a homogenous distribution of the media.

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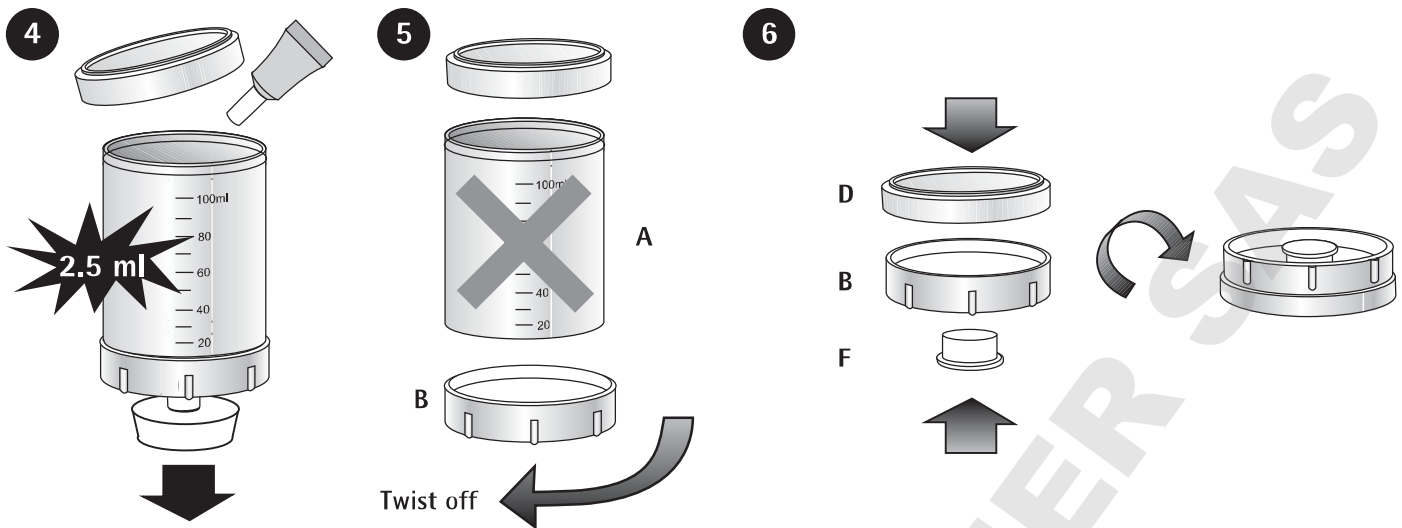
Nitrocellulose membrane filters	
Hazard statements	Precautionary statements
H228: Flammable Solid	P 210: Keep away from heat/sparks/open flames/hot surfaces. —No smoking.
	P 243: Take precautionary measures against static discharge.
DANGER	ADR 286: IATA A122

Producer: Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
D-37079 Goettingen, Germany
Phone: +49.551.308.0

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Phone +49.551.308.0
Fax +49.551.308.3289

www.sartorius.com

Specifications subject to change without notice. Printed in the EU on paper bleached without chlorine.
Publication No.: SLD6030-p150703
Ver. 07 | 2015

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Hersteller: Sartorius Stedim Biotech GmbH
Producteur: Sartorius Stedim Biotech GmbH
Produttore: Sartorius Stedim Biotech GmbH
Fabricante: Sartorius Stedim Biotech GmbH
Sartorius Lab Instruments GmbH & Co.KG