# 16306, 16307, 16315, 16316

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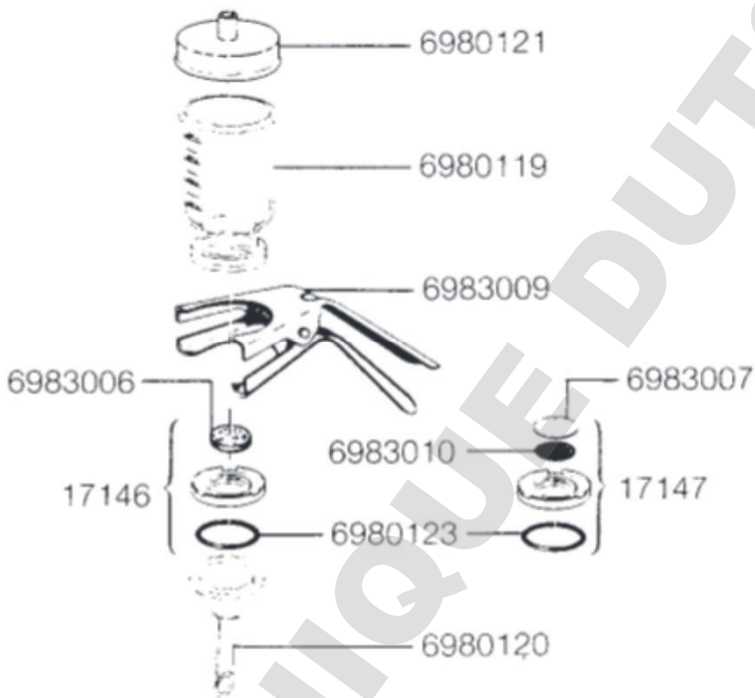
**16307**  
**16316**

## Table of Contents

1. Specifications
2. Equipment Supplied
3. Cleaning and Care of the Filter Holder
4. How to Assemble the Filter Holder
5. Clarifying Filtration
6. Sterile Filtration
7. Microbiological Testing
8. Analytical Methods
9. Recommended Accessories
10. Spare Parts

Glass vacuum filter holders are available in two types which differ only in the filter support, which is a PTFE ring. This ring contains either a glass frit or a PTFE-coated stainless steel screen and an underdrain screen.

The filter supports may only be changed as a unit. Because of the particularly even distribution of particulates of the membrane filter surface, filter supports with glass frits are preferred for use in particulate analysis.

Filter supports with a PTFE-coated screen are mainly used for clarifying or sterile filtration of liter volumes.

## 1. Specifications

Materials: Borosilicate glass (3.3 DIN-ISO 3585)

Clamp: Anodized aluminum

Filter support: a) Sintered glass frit in a PTFE ring, with a centering rim on top and silicone O-ring on the bottom, or b) PTFE-coated stainless steel screen on an underdrain screen in a PTFE ring, with a centering rim on top and a silicone O-ring on the bottom.

Filter Holder No.	16306* 16315**	16307* 16316**
Silicone O-ring	25 × 3 mm	45 × 3 mm
Filter diameter	25 mm (or 24)	50 mm (or 47)
Prefilter diameter	20 mm	40 mm
Filtration area	3 cm <sup>2</sup>	12.5 cm <sup>2</sup>
Capacity	30 ml	250 ml
Connections:		
Outer diameter of outlet	12 mm	15 mm
Operating pressure	vacuum only	vacuum only
Sterilization:	Autoclave up to 134°C maximum, or use dry heat up to 180°C maximum depending on the membrane filter used.	

\* with a glass frit


\*\* with a PTFE-coated stainless steel screen

## 2. Equipment supplied

Remove the filter holder from the packaging and check the list below to make sure you have all the necessary parts:

Equipment supplied	Shown in Fig.1 as:
1 Lid (16307 and 16316 only)	No. 1
1 Glass funnel	No. 2
1 Clamp	No. 3
1 Filter support with glass frit or PTFE-coated support screen and PTFE underdrain screen	No. 4
1 Silicone O-ring	No. 5
1 Glass base	No. 6

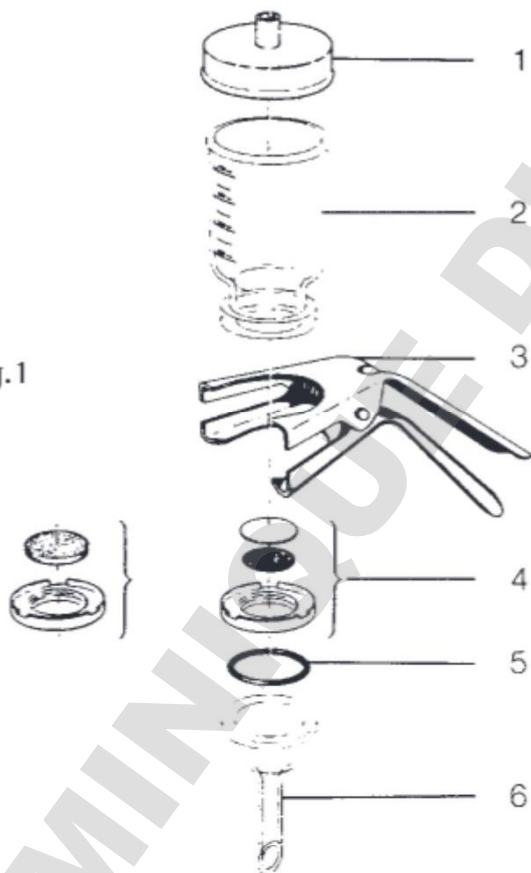Fig.1

### 3. Cleaning and Care of the Filter Holder

To ensure long service life and proper functioning of the filter holder, it must be cleaned before the first use and after every filtration run.

Proceed as follows:

- 3.1 Completely disassemble the filter holder as indicated in the list of parts (see Fig. 1).
- 3.2 Use a soft brush, warm water, and a gentle, commercially available laboratory detergent (for metal, glass and plastics) to clean all parts.
- 3.3 Rinse all parts first with hot water, then with distilled water.
- 3.4 Dry the parts with compressed air, or let them drip dry. Do not use towels or cloth as they may leave lint on the filter holder.

### 4. How to Assemble the Filter Holder and Insert the Membrane Filter

To assemble the vacuum filtration system, you will need in addition:

- 1 Stopper (see “Recommended Accessories”)
- 1 Suction flask (16672)
- 1 Vacuum source (e.g., 16612 or 16615)
- Vacuum hose (16623)
- 1 Woulff’s bottle (16610) or 1 Vacusart (17804-M)
- Pair of forceps for handling the membrane filter (see Fig. 2).

For large numbers of routine tests, we recommend our 3-place or 6-place vacuum manifold along with the corresponding number of filter holders and connectors. (See “Recommended Accessories.”)

- 4.1 Place the appropriate stopper on the glass base and insert it into the suction flask (see Fig. 3).
- 4.2 Using a vacuum hose, connect the suction flask and vacuum source, such as a water jet pump (see Fig. 4). If an electric pump is used as a vacuum source, a Woulff's bottle (see Fig. 5) or a Vacusart (17804-M) must be connected between the suction flask and the pump in order to prevent the filtrate from foaming over and entering the pump. When installing the Vacusart (17804-M), be sure that any filtrate overflow can return to the suction flask, and will not be entrapped in the hose.
- 4.3 Press-fit the silicone O-ring on the underside of the filter support and place the glass frit or support screen and underdrain screen on top. To do so, place the underdrain screen with the smooth side up in the filter support and press on the PTFE-coated support screen with the outside edge facing down so that it fits exactly into the indentation. Then put the filter support on the glass base (see Fig. 6).
- 4.4 Using forceps, remove the membrane filter from the packaging and center it on the filter support. If you need to use a prefilter, center it on the membrane filter at this point. The diameter of the prefilter may not exceed the dimensions listed in the "Specifications" (see Fig. 7).
- 4.5 Set the glass funnel on top of the base so that it fits exactly and clamp the two parts together (see Fig. 8).

## 5. Clarifying Filtration

- 5.1 Assemble the filter holder and insert the membrane filter as described in section 4.
- 5.2 Pour the liquid to be filtered into the glass funnel (see Fig. 9) and switch on the vacuum pump.

### **Important Note:**

If you are additionally using a prefilter, switch the vacuum pump on before filling the funnel with the liquid to be filtered. (Otherwise, the prefilter may float to the top of the liquid.)

- 5.3 Filtration is completed when all the liquid has passed through the membrane filter. Disconnect the vacuum source from the filter holder by pulling the hose off the hose nipple on the suction flask (if the water jet pump is used; see Fig. 10). If a Woulff's bottle and a vacuum pump have been used, switch off the pump and release vacuum via the Woulff's bottle (see Fig. 11).
- 5.4 Now remove the filter holder and stopper from the suction flask and transfer the filtrate to another container or leave it in the suction flask.

## 6. Sterile Filtration

(This is possible only with the filter support and PTFE-coated screen, and not the glass filter support.)

- 6.1 Assemble the filter holder and insert the membrane filter as described in section 4; however, do not connect the filter holder to a vacuum source or a Woulff's bottle.

**Important Note:**

Wet cellulose nitrate filters before placing them in wet filter holders. If the filter holder is dry, cellulose nitrate filter must also be dry. Always make it a point to wet cellulose acetate filters and nylon filters (Sartolon) before placing them in a filter holder.

- 6.2 Use a short length of vacuum tubing to attach a Vacusart (17804-M) to the nipple of the suction bottle (see Fig. 12).
- 6.3 Close the glass funnel with either the silicone lid and a piece of cotton or with aluminum foil (see Fig. 13).
- 6.4 Autoclave the filter holder at 121°C and 1 bar (14.5 psi) for 30 minutes. (This may be done with all Sartorius Stedim Biotech membrane filters except for gelatin and polyamide filters.)

Cellulose nitrate membrane filters (111..) and PTFE membrane filters (118..) as well as glass fiber filters (134..) may be sterilized by dry heat for either 2 hours at 180°C or for 3 hours at 160°C. (Do not wet the filters before placing them in filter holders.)

- 6.5 After sterilizing the assembled filtration unit, proceed with the filtration run as with clarifying filtration. Be sure to maintain sterile conditions.

## 7. Microbiological Testing

For microbiological testing, it is necessary to sterilize the filter holder containing the membrane filter. Either autoclave the filter at 121°C for 30 min., or use dry heat at 180°C for 2 hours. After sterilizing, perform the filtration run, then remove the membrane filter and incubate it on the required culture medium|media.

## 8. Analytical Methods

Sartorius Stedim Biotech Vacuum Filter Holders made of glass are especially suited for analysis, e.g., gravimetric determinations, particulate analysis, etc.

## 9. Recommended Accessories

Product Name	Filter Holder No.	
	16306*	16307*
	16315**	16316**
Fluoroelastomer O-ring	00118	00124
Stopper	17174	17175
Stainless steel adapter with stopper for stainless steel 3-place and 6-place manifold	16836	16837
Stainless steel 3-place manifold	16842	16842
Stainless steel 6-place manifold	16843	16843
Incubator 220 V, 50 Hz	–	18113

\* with glass frit filter support

\*\* with PTFE-coated screen for the filter support

### These additional accessories can be used with all filter holders:

16672	Suction flask, 2 l
16623	Rubber vacuum hose, 1 m
16611	Water jet pump with G 3/4" female thread
16612	Laboratory vacuum pump, 220 V, 50 Hz
16615	Laboratory vacuum pump, 110 V, 60 Hz
16692	Vacuum pump, 220 V, 50 Hz
16695	Vacuum pump, 110 V, 60 Hz
17804-M	Vacusart
16625	Stainless steel forceps

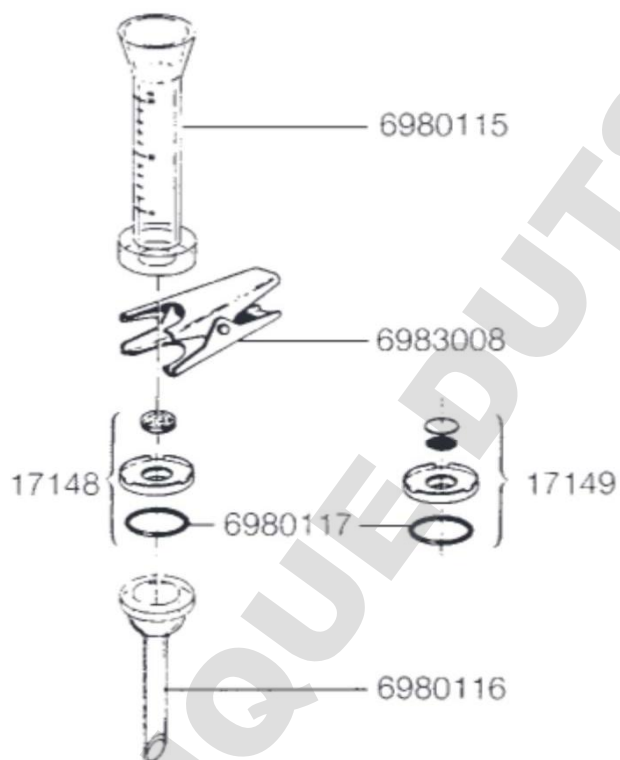
## 10. Spare parts (see diagram)

### 16306, 16315:

- 6980115 Glass funnel, 30 ml
- 6980116 Glass base, 25 mm
- 6980117 Silicone O-ring, 25 × 3 mm
- 6983008 Clamp
- 17148 PTFE ring with glass frit and silicone O-ring (for 16306)
- 17149 PTFE ring with PTFE-coated underdrain screen and support screen including a silicone O-ring (for 16315)

### 16306


### 16315

**16307, 16316:**

- 6980119 Glass funnel, 250 ml
- 6980120 Glass base, 50 mm
- 6980121 Silicone rubber lid
- 6980123 Silicone O-ring, 45 × 3 mm
- 6983006 Glass frit
- 6983007 PTFE-coated support screen
- 6983009 Clamp
- 6983010 Underdrain screen, PTFE-coated
- 17146 PTFE ring with glass frit and silicone O-ring (for 16307)
- 17147 PTFE ring with PTFE-coated underdrain screen and support screen including a silicone O-ring (for 16316)

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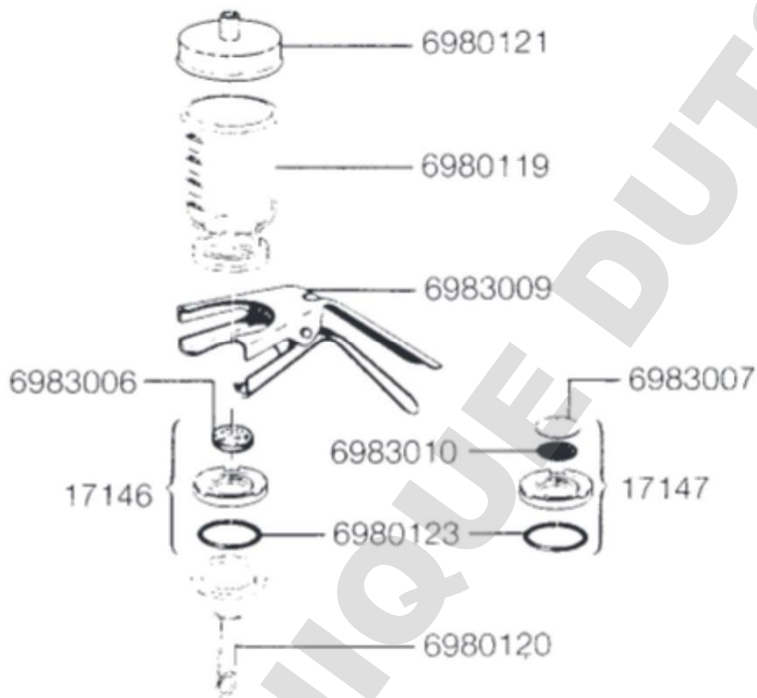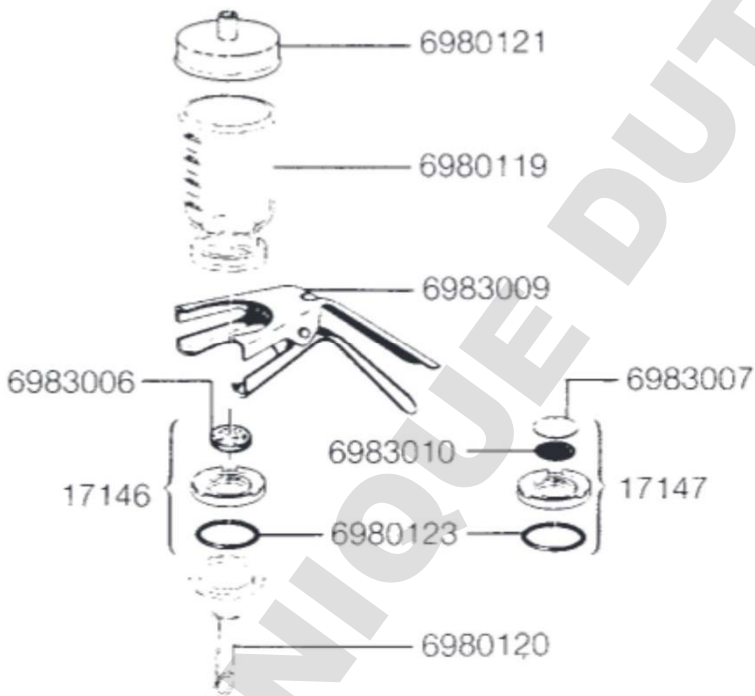
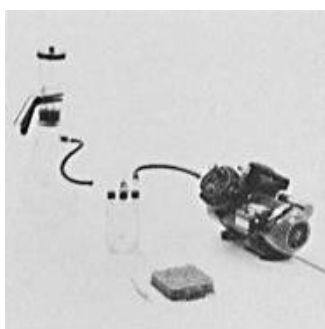
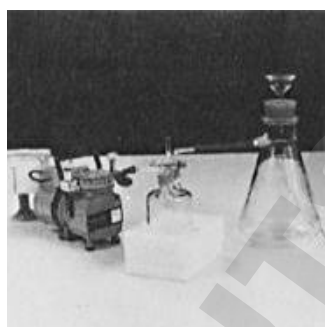
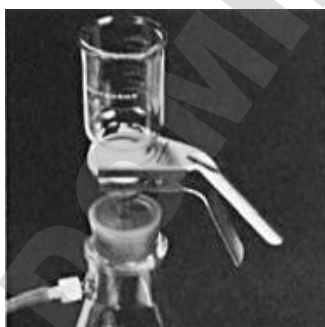
[www.sartorius-stedim.com](http://www.sartorius-stedim.com)

Copyright by Sartorius Stedim  
Biotech GmbH, Goettingen,  
Germany.

All rights reserved. No part  
of this publication may be  
reprinted or translated in any  
form or by any means without  
the prior written permission  
of Sartorius Stedim Biotech  
GmbH. The status of the  
information, specifications  
and illustrations in this  
manual is indicated by the  
date given below.

Sartorius Stedim Biotech GmbH  
reserves the right to make  
changes to the technology,  
features, specifications and  
design of the equipment  
without notice.

Status:

May 2014,

Sartorius Stedim Biotech GmbH,  
Goettingen, Germany

Printed in the EU on paper bleached  
without chlorine. | W  
Publication No.: SL-6001-p140509  
Ver. 05 | 2014