



Midisart® 2000

Ready-to-use filter for sterilizing gases and venting

Product Information

The Midisart® 2000 is the best choice for small-scale sterile air | gas filtration applications. Superior flow rates and seamless integration into single-use assemblies provide Midisart® customers with a distinct advantage that directly translates into cost savings. Individual 100% factory integrity testing verifies the quality of every filter prior to use. Moreover our new production-line with its enhanced capacity ensures a quick and reliable supply at any time.



Applications

Midisart® are excellent for

- Sterile venting of filling vessels and fermentation carboys, including culture vessels and CO₂ incubators (6 to 120 liters)
- Venting of holding tanks for sterile, distilled water and liquid culture media
- Autoclave venting
- In-line sterilization of and particulate removal from air and gases, such as sterilization of air for small fermenters

Flexibility

Midisart® 2000 were designed for maximum handling ease and safety. Tapered hose barbs ensure a simple and secure hold for 6- to 12-mm inner diameter tubing. Other connector types such as small hose barb (for tubings with 4 – 12 mm inner diameter), 8" NPT thread and TriClamp are also available. Midisart® is lightweight – only 20 g – so it will not weigh down or kink the tubing.

Reliability and safety

Midisart® 2000 is integrity testable and delivers reproducible results. The membrane is reinforced with polypropylene gauze, giving the Midisart® unit added stability and making it pressure resistant up to 3 bar | 44 psi. Midisart® 2000 entirely eliminates moisture breakthrough because of its inherently hydrophobic PTFE material. In addition, Midisart® is bio safe because all materi-

als of construction meet the requirements of the current USP Plastics Class VI testing. Midisart® 2000 units easily withstand at least 20 autoclaving cycles with no loss in performance.

Quality Control

Each Midisart® 2000 is tested by integrity prior to release. This 100% test has been defined as the final release criteria. The lot number and the individual unit number are lasered on the top part of each Midisart® 2000 housing to ensure complete traceability.

Performance

With a diameter measuring just 64 mm, Midisart® incorporates a filter area of 20 cm², which means that it is "packed" with high flow rate performance power. Midisart® multiply filtration performance in more ways than one.

Lot Number and Individual Unit Number

Lasered on the top part of each Midisart® - housing.

Integrity Testing

Midisart® 2000 filters can be integrity tested (Bubble Point Test, 60% IPA, BP_{min.} = 1.1 bar). The Midisart® Test Manifold (Order code 1Z-LB-0002) can be used to test 10 filters in parallel which simply leads to a reduction of test time by 90%.

Technical Data

Specifications

Materials

Membrane PTFE – reinforced with polypropylene gauze
Housing Polypropylene

Pore Sizes

0.2 µm
0.45 µm

Filtration Area

20 cm² | 3 square inch

Housing diameter

64 mm | 2,5"

Connectors

Hose Barb, 1/8" NPT, Triclamp, small Hose Barb

Biosafety

Materials pass current UPS Plastics, class VI
Non fiber releasing according to 21 CFR

Sterilization

Max. Temperature 134°C
Max. Autoclaving 60 Cycles
Not suitable for gamma irradiation
Gamma stable: see Midisart® BV

Technical References

Validation Guide

SLD5702-e

DOMINIQUE DUTSCHER SAS

Ordering Information

Order No.	Pore Size [µm]	Membrane	Connectors	Qty.	Sterile
17804-----E	0.45	PTFE	Hose Barb Hose Barb	12	Yes
17804-----G	0.45	PTFE	Hose Barb Hose Barb	25	Yes
17804-----NPE	0.45	PTFE	1/8" 1/8" NPT	12	Yes
17804-----NPG	0.45	PTFE	1/8" 1/8" NPT	25	Yes
17805-----E	0.2	PTFE	Hose Barb Hose Barb	12	Yes
17805-----G	0.2	PTFE	Hose Barb Hose Barb	25	Yes
17805-----NPE	0.2	PTFE	1/8" 1/8" NPT	12	Yes
17805-----NPG	0.2	PTFE	1/8" 1/8" NPT	25	Yes
17805-----UPN	0.2	PTFE	Hose Barb Hose Barb	100	No
17805-----UPQ	0.2	PTFE	Hose Barb Hose Barb	500	No
17805-----UQN	0.2	PTFE	Hose Barb Hose Barb	100*	No
17809-----UNN	0.2	PTFE	1/8" 1/8" NPT	100	No
17812-----UNN	0.2	PTFE	1/8" Hose Barb	100	No
17805-----TCN	0.2	PTFE	TriClamp TriClamp	100	No
17877-----UPN	0.2	PTFE	Small Hose Barb Small Hose Barb	100	No

* = plastic bucket – the bucket with its smooth surface can be easily transferred into clean-room areas without risking particle release from the packaging material.



Standard Hose Barb



Small Hose Barb



1/8" NPT Thread



TriClamp

Sales and Service Contacts

For further contacts, visit www.sartorius-stedim.com

Europe

Germany
Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen
Phone +49,551.308.0
Fax +49,551.308.3289

Sartorius Stedim Systems GmbH
Robert-Bosch-Strasse 5-7
34302 Guxhagen
Phone +49,5665.407.0
Fax +49,5665.407.2200

France
Sartorius Stedim FMT S.A.S.
ZI des Paluds
Avenue de Jouques - CS 91051
13781 Aubagne Cedex
Phone +33,442.845600
Fax +33,442.845619

Sartorius Stedim France SAS
ZI des Paluds
Avenue de Jouques - CS 71058
13781 Aubagne Cedex
Phone +33,442.845600
Fax +33,442.846545

Austria
Sartorius Stedim Austria GmbH
Modcenterstrasse 22
1030 Vienna
Phone +43,1.7965763.18
Fax +43,1.796576344

Belgium
Sartorius Stedim Belgium N.V.
Leuvensesteenweg, 248/B
1800 Vilvoorde
Phone +32,2.756.06.80
Fax +32,2.756.06,81

Hungary
Sartorius Stedim Hungária Kft.
Kagyló u. 5
2092 Budakeszi
Phone +36,23.457.227
Fax +36,23.457.147

Italy
Sartorius Stedim Italy S.p.A.
Via dell'Antella, 76/A
50012 Antella-Bagno a Ripoli (FI)
Phone +39,055,63,40,41
Fax +39,055,63,40,526

Netherlands
Sartorius Stedim Netherlands B.V.
Phone +31.30.60.25.080
Fax +31.30.60.25.099
filtratie.nederland@sartorius-stedim.com

Poland
Sartorius Stedim Poland Sp. z o.o.
ul. Wrzesinska 70
62-025 Kostrzyn
Phone +48,61,647,38,40
Fax +48,61,879,25,04

Russian Federation
LLC "Sartorius ICR" and LLC "Biohit"
Uralskaya str. 4, Lit. B
199155, Saint-Petersburg
Phone +7,812,327,5,327
Fax +7,812,327,5,323

Spain
Sartorius Stedim Spain, S.A.U.
Avda. de la Industria, 32
Edificio PAYMA
28108 Alcobendas (Madrid)
Phone +34,902,110,935
Fax +34,91,358,96,23

Switzerland
Sartorius Stedim Switzerland AG
Ringstrasse 24 a
8317 Tagelswangen
Phone +41,52,354,36,36
Fax +41,52,354,36,46

U.K.
Sartorius Stedim UK Ltd.
Longmead Business Centre
Blenheim Road, Epsom
Surrey KT19 9 QQ
Phone +44,1372,737159
Fax +44,1372,726171

Ukraine
LLC "Biohit"
Post Box 440 "B"
01001 Kiev, Ukraine
Phone +380,44,41,1,4918
Fax +380,50,623,3162

America

USA
Sartorius Stedim North America Inc.
5 Orville Drive, Suite 200
Bohemia, NY 11716
Toll-Free +1,800,368,7178
Fax +1,631,254,4253

Argentina
Sartorius Argentina S.A.
Int. A. Ávalos 4251
B1605ECS Munro
Buenos Aires
Phone +54,11,4721,0505
Fax +54,11,4762,2333

Brazil
Sartorius do Brasil Ltda
Avenida Senador Vergueiro 2962
São Bernardo do Campo
CEP 09600-000 - SP- Brasil
Phone +55,11,4362,8900
Fax + 55,11,4362,8901

Mexico
Sartorius de México S.A. de C.V.
Circuito Circunvalación Poniente
No. 149
Ciudad Satélite
53100, Estado de México
México
Phone +52,5555,62,1102
Fax +52,5555,62,2942

Asia | Pacific

Australia
Sartorius Stedim Australia Pty. Ltd.
Unit 5, 7-11 Rodeo Drive
Dandenong South Vic 3175
Phone +61,3,8762,1800
Fax +61,3,8762,1828

China
Sartorius Stedim Biotech (Beijing) Co. Ltd.
No. 33 Yu'an Road
Airport Industrial Park Zone B
Shunyi District, Beijing 101300
Phone +86,10,80426516
Fax +86,10,80426580

Sartorius Stedim (Shanghai)
Trading Co. Ltd
3rd Floor, North Wing, Tower 1
No. 4560 Jinke Road
Zhangjiang Hi-Tech Park
Pudong District
Shanghai 201210, China
Phone +86,21,68782300
Fax +86,21,68782332 | 68782882

Sartorius Stedim Biotech (Beijing) Co. Ltd.
Guangzhou Representative Office
Unit K, Building 23
Huihua Commerce & Trade Building
No. 80 Xianlie Middle Road
Guangzhou 510070
Phone +86,20,37618687 | 37618651
Fax +86,20,37619051

India
Sartorius Stedim India Pvt. Ltd.
#69/2-69/3, NH 48, Jakkasandra
Nelamangala Tq
562 123 Bangalore, India
Phone +91,80,4350,5250
Fax +91,80,4350,5253

Japan
Sartorius Stedim Japan K.K.
4th Fl., Daiwa Shinagawa North Bldg.
8-11, Kita-Shinagawa 1-chome
Shinagawa-ku, Tokyo, 140-0001 Japan
Phone +81,3,4331,4300
Fax +81,3,4331,4301

Malaysia
Sartorius Stedim Malaysia Sdn. Bhd.
Lot L3-E-3B, Enterprise 4
Technology Park Malaysia
Bukit Jalil
57000 Kuala Lumpur, Malaysia
Phone +60,3,8996,0622
Fax +60,3,8996,0755

Singapore
Sartorius Stedim Singapore Pte. Ltd.
1 Science Park Road,
The Capricorn, #05-08A,
Singapore Science Park II
Singapore 117528
Phone +65,6872,3966
Fax +65,6778,2494

South Korea
Sartorius Korea Biotech Co., Ltd.
8th Floor, Solid Space B/D,
PanGyoYeok-Ro 220, Bundang-Gu
SeongNam-Si, GyeongGi-Do, 463-400
Phone +82,31,622,5700
Fax +82,31,622,5799



◀ www.sartorius-stedim.com