sartorius

Operating Instructions | Betriebsanleitung | Mode d'emploi |  
Istruzioni per l'uso | Instrucciones de funcionamiento |  
操作说明 | Инструкция по эксплуатации | 操作指示 | ارشادات التشغيل

## YDP40

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## 1 Safety Precautions

### Safety

- The device complies with EU Directives and standards for electrical safety and electromagnetic compatibility. Improper use or handling, however, can result in damage and/or injury. Any improper use or operation of the device, i.e. that is not consistent with the instructions, will result in forfeiture of all claims under the manufacturer's warranty.
- Personnel need to have read and understood these installation instructions, including the safety instructions.
- In the event of use in systems and ambient conditions which have greater safety requirements, you must observe the requirements and provisions applicable in your country.

### Danger of Explosion!

Do not use this equipment in hazardous areas in which explosive materials are present.

Make sure that the voltage rating printed on the AC adapter is identical to your local mains voltage.

### Installation

Do not operate the device if the housing or AC adapter is damaged. Immediately disconnect the damaged device from the power by pulling the plug.

Do not expose the equipment to extreme temperatures, aggressive chemical vapors, moisture, shocks, vibration, or strong electromagnetic fields. Observe the conditions of operation described in the specifications!

### Installation information:

The operator shall be solely responsible for any modifications to the equipment and for connecting any cables or equipment not supplied by Sartorius. Information on operational quality is available upon request from Sartorius.

### Warning!

Only operate the printer with the supplied AC adapter HDAD36W101. Pay attention to the information on the AC adapter.

Do not compromise the IP protection of the device.

Before cleaning the AC adapter:  
Unplug from the mains power.

The device may only be opened  
by specialized personnel trained  
by Sartorius. Do not open the AC  
adapter.


Use only a slightly moistened cloth  
to clean the equipment and make  
sure that no liquid enters the printer  
housing.

Caution: The serrated edge for  
tearing off printouts is sharp;  
direct contact can result in injury.


## 2 Unpacking and Checking

Open the packaging and check the items according to the packing list.  
Please contact your local dealer if any item is missing or damaged.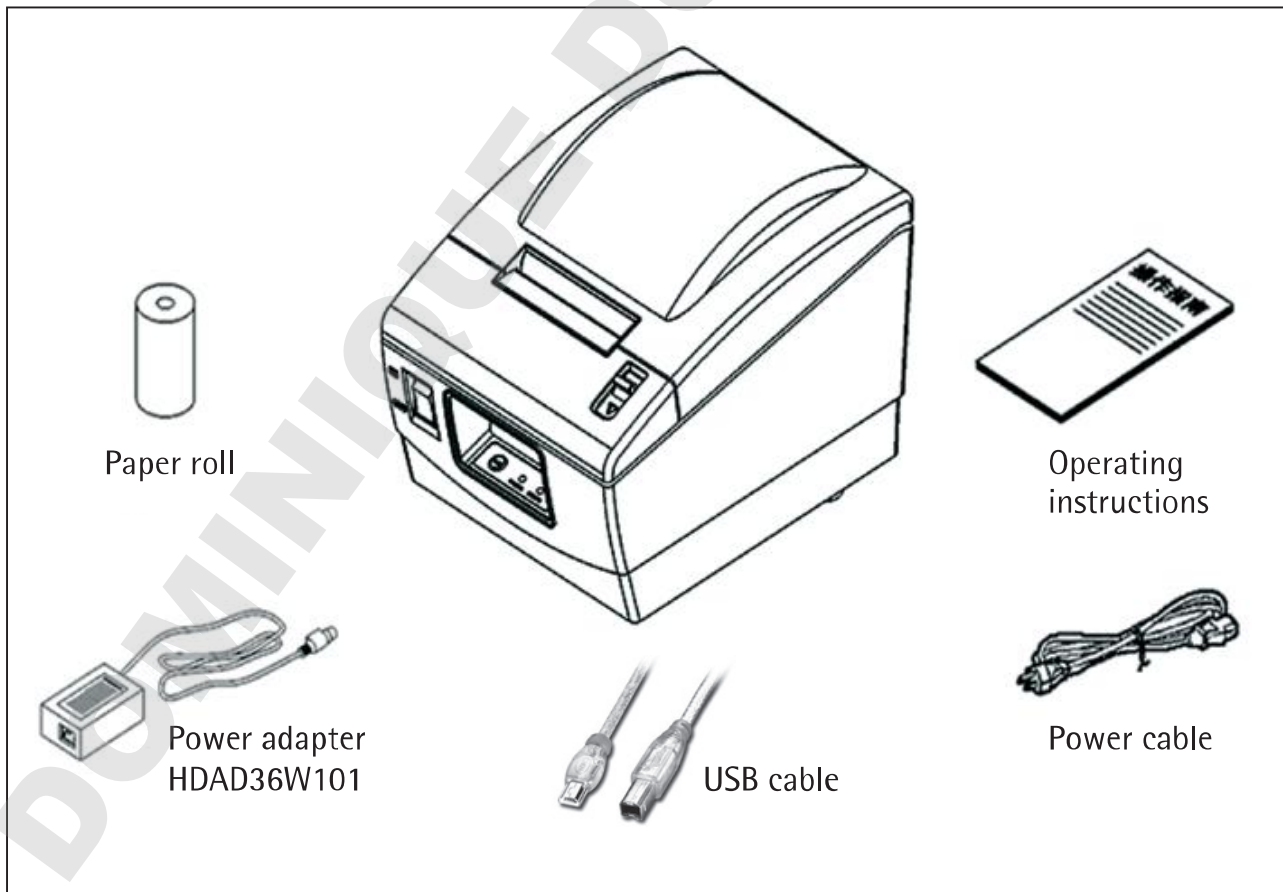

### 3 Printer Introduction

#### 1. Appearance and design (see the figure below)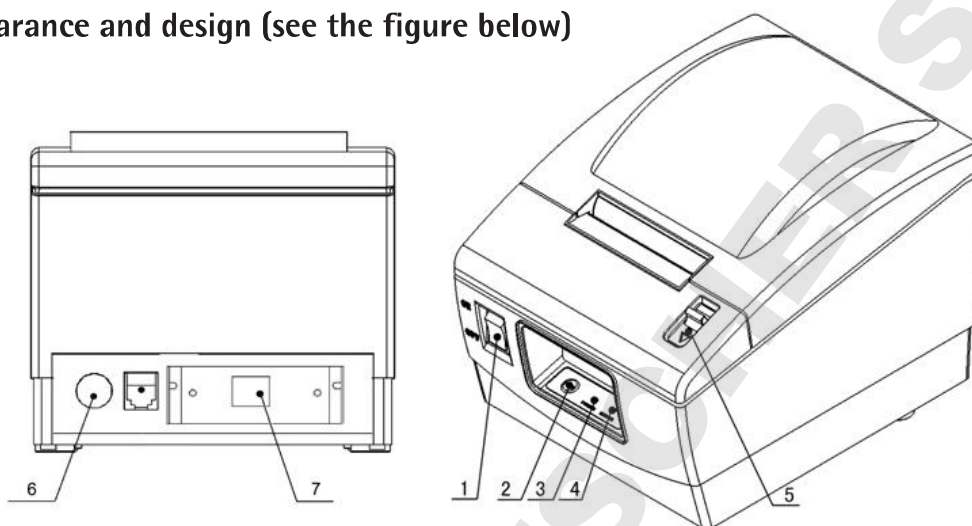

1 Power switch  
2 FEED button

3 ERROR LED  
4 POWER LED

5 Cover open lever  
6 Power interface  
7 Communication interface

#### 2. Specifications

Print method:	Direct thermal line
Print speed:	100 mm/s
Resolution:	203 × 203 dpi
Print width:	Max. 56 mm
Power supply:	DC 24 V, 1.5 A
Paper roll diameter:	Max. 80 mm
Paper type:	Thermal roll paper (external)
Operating conditions:	5...45°C, 20...90% RH (at 40°C)
Storage conditions:	-20...60°C, 20...90% RH (at 40°C)
Printer head lifetime:	50 km
Dimensions:	125 (w) × 200 (d) × 125 (h) mm
Weight:	1.2 kg
Paper thickness:	0.06 ... 0.1 mm
Paper width:	57.5 ± 0.5 mm
Paper type:	Continuous paper


#### Recommended paper:

Standard thermal paper: 69Y03287

#### Caution:


To ensure quality print and normal lifetime of print head and printer, use recommended paper supplied by Sartorius.

## 4 Printer Installation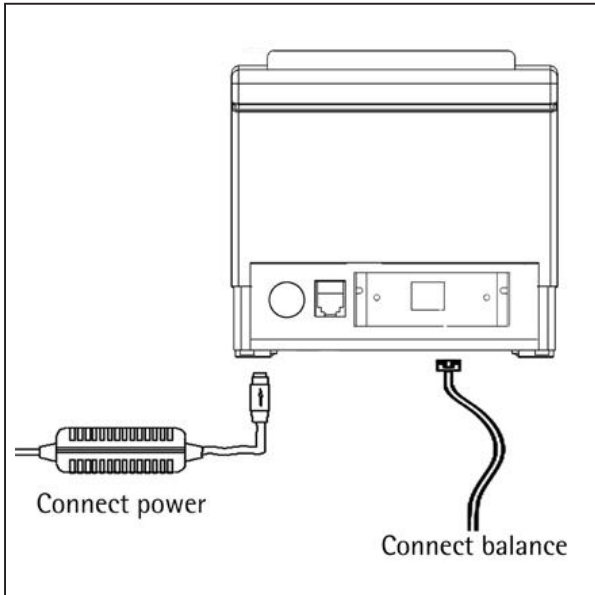


Check to make sure that the power switch of the printer is in the OFF position before connecting the printer according to the figure above.


1. Press the cover open lever and open the top cover.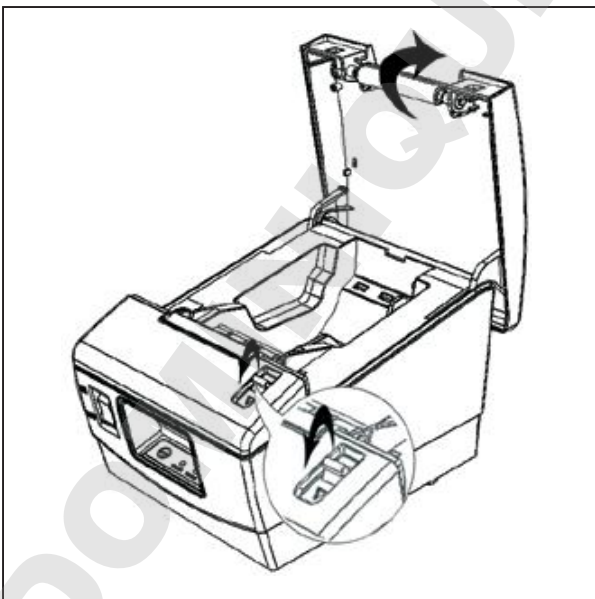

2. Load the paper roll into the unit, as shown in the figure.

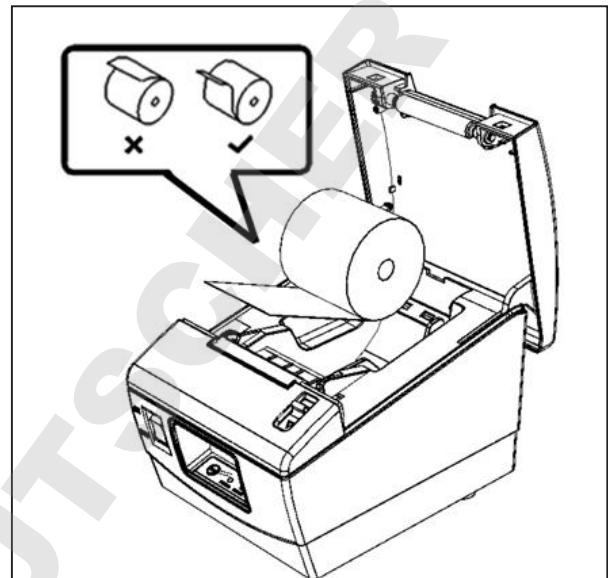
3. Switch the printer on via the power switch (1).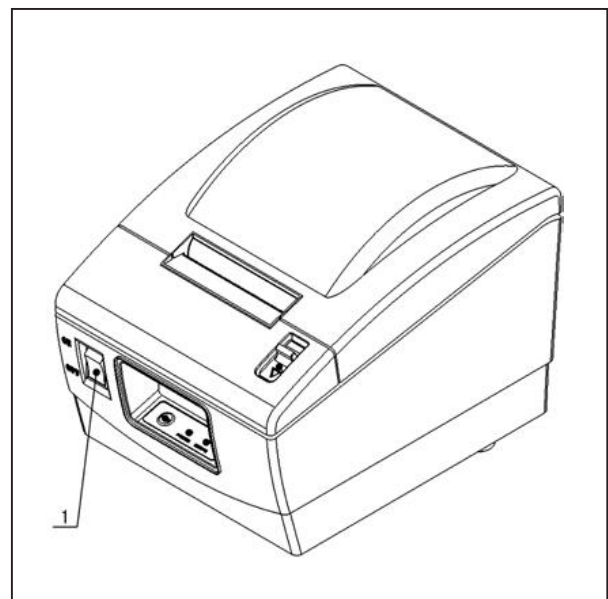

## 5 Printing a Self-test Page


4. Close the top cover and tear off the surplus paper to complete the installation.

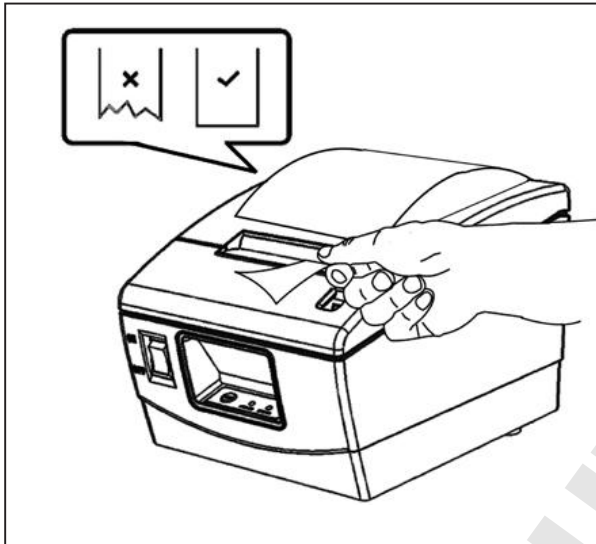Printing a self-test page is used to check whether the printer is operating in the normal mode.

1. Make sure that the paper roll has been loaded and the top cover has been closed.
2. Press and hold the feed button while turning on the printer. After hearing the printer start to feed paper, release the button; the printer will now print a self-test page. Manually tear off the paper after the printer stops feeding paper.
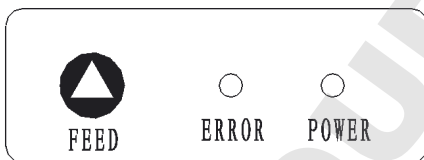
## 6 Operating

### FEED Button

If you press the FEED button for a short time, the printer will feed one line at a time. If you hold the FEED button, the printer will feed paper constantly.

Name	Status	Explanation
POWER LED (green)	Always on	The power supply of the printer is normal.
	Off	Printer is turned off.
ERROR LED (red)	Off	Printer is normal.
		Out of paper.
		Voltage is abnormal.
		Print head is overheating.

---

## 7 Routine Maintenance

### ○ Cleaning the print head and platen roller

- 1) Turn off the printer power and open the top cover.
- 2) If printing has just finished, please wait until the print head has completely cooled.
- 3) Use alcohol cotton (should be wrung out) to wipe off the stains and dust on the surface of the print head and platen roller.
- 4) Once the alcohol has completely evaporated, close the top cover.

### ⚠ **Caution:**

- Do not use hard objects (for example: forceps) on the print head and platen roller.
- To increase the lifetime of the print head, it is recommended to clean the print head and platen roller every month. If the environment is dirty, increase the cleaning times accordingly.

### ○ Cleaning the sensor

- 1) Turn off the printer power and open the top cover.
- 2) Use alcohol cotton (should be wrung out) to wipe off the stains and dust on the surface of the sensor.
- 3) Once the alcohol has completely evaporated, close the top cover.

### ○ Clearing a paper jam

- 1) Turn off the printer and press the top cover open lever to open the top cover.
- 2) Clear the paper jam and close the top cover.

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Copyright by Sartorius, Goettingen, Germany.  
All rights reserved. No part of this publication  
may be reprinted or translated in any form  
or by any means without the prior written  
permission of Sartorius.

The status of the information, specifications  
and illustrations in this manual is indicated  
by the date given below.

Sartorius reserves the right to make changes  
to the technology, features, specifications  
and design of the equipment without notice.

Status:  
April 2016,  
Sartorius Lab Instruments GmbH & Co. KG,  
Goettingen, Germany

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Ltd.No. 33 Yu An Road,