### Operation


The vacuum is controlled by the pressure sensitive turquoise control button on the top of the hand piece. By means of the plastic screw opposite the control button, the residual suction level can be set, which will prevent dripping of waste liquids from the tip or pipette. The aspirator is factory pre-set to allow a low level of constant suction; by turning the screw one full turn clockwise, the valve is shut and residual vacuum is stopped.

### Cleaning and servicing

The Hand Operator is entirely autoclavable. When working with biohazardous material, sterilise the unit before disassembling for cleaning. Please refer to the explosion drawing.

To clean, immerse the pieces in mild laboratory detergent, but do not use abrasives, as a roughened surface may affect the smooth operation of the control valve. Apply a thin coat of MOLYKOTE HP-870 (Dow Corning) lubricant as needed to the control valve to keep the action smooth (arrow).

When working with salt-rich solutions, rinse out the aspirator hand piece with distilled water after use to avoid build-up of dried deposits.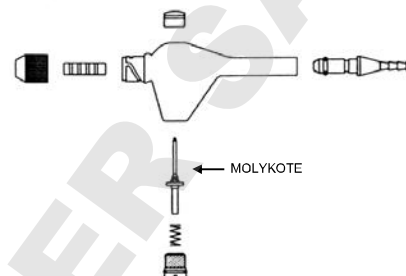

### Chemical compatibility

Avoid prolonged contact (days or weeks) with:

- concentrated inorganic acids
- sodium hypochlorite
- methylene chloride
- concentrated aldehydes
- other halogenated hydrocarbons


Resistant to:

- DMSO
- ethanol/methanol
- dimethylformamide
- acetone

Note:

- Flushing the aspirator with ethanol immediately after contact with organic solvents is the best safeguard in all cases.
- For o-ring replacement, use Viton o-rings only

### Accessories and consumables

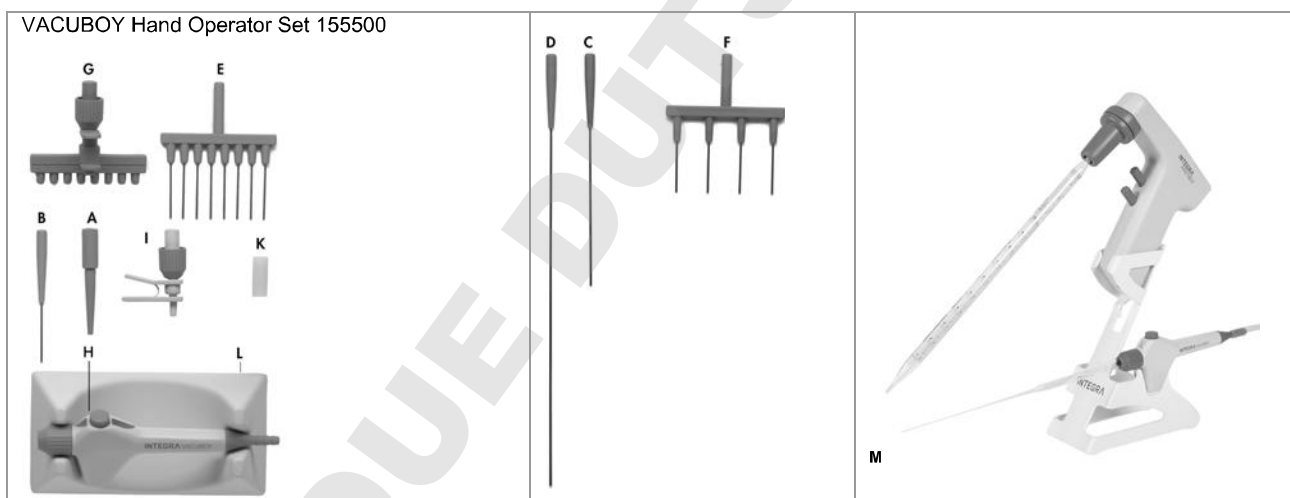<b>VACUBOY Hand Operator Set 155500</b>	
<b>VACUBOY Hand Operator:</b> hand operator (H), two aspiration adapters (B and I), rubber adapter for Pasteur pipettes (K), silicone grease	155510
<b>VACUBOY Hand Operator Set:</b> hand operator (H), aspiration adapter A, B, E, G, I, rubber adapter for Pasteur pipettes (K), stand for hand operator (L), silicone tubing 1800 mm, silicone grease	155500
<b>A</b> 1-channel plastic aspiration adapter for disposable tips	155504
<b>I</b> 1-channel plastic aspiration adapter for disposable tips with ejector	155526
<b>C</b> 1-channel stainless steel aspiration adapter 150 mm	155522
<b>D</b> 1-channel stainless steel aspiration adapter 280 mm	155525
<b>B</b> 1-channel stainless steel aspiration adapter 40 mm	155502
<b>F</b> 4-channel stainless steel aspiration adapter 40 mm	155524
<b>G</b> 8-channel plastic aspiration adapter for disposable tips with ejector	155520
<b>E</b> 8-channel stainless steel aspiration adapter 40 mm	155503
<b>K</b> Rubber adapter for Pasteur pipettes	155505
<b>L</b> Stand for VACUBOY hand operator	155501
Silicone tubing 1800 mm	158332
<b>M</b> Stand for PIPETBOY ( <i>acu</i> or <i>pro</i> ) and VACUBOY	155065

**Manufacturer:** INTEGRA Biosciences AG  
 CH-7205 Zizers, Switzerland  
[www.integra-biosciences.com](http://www.integra-biosciences.com)  
[info@integra-biosciences.com](mailto:info@integra-biosciences.com)

**Customer service:** Please contact your local INTEGRA Biosciences AG representative. You can find the name and address on [www.integra-biosciences.com](http://www.integra-biosciences.com).

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**Fabricante:** INTEGRA Biosciences AG