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### Instructions

TPP tissue culture flasks are intended for cell culture with necessary sterile handling techniques. The growth area is opto-mechanically treated for optimal cell adhesion and proliferation. The angled neck decreases the risk of media contact with the inside of the screw cap.

- Open the product packaging and inspect for intact bag and product.
- Fill medium: Introduce medium or inoculate cells with a pipette alongside the wall, or direct injection on the growth area of the flask.
- Tightly screw closed the filter screw cap with the integrated hydrophobic membrane for sterile gas exchange
- Incubate: Lay the flask on the incubator shelves for short or long term storage.
- **Open re-closable flask:** Complete access from above to flask chamber with multi-use re-closable liquid-tight lid.
- Harvest cells with large access from above to flask chamber
  - Harvest adherent cells by trypsinization or scrape with the cell scraper.
  - Harvest suspension cells by exchange of medium.

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**Advice and Recommendation**

Stable ring enables slip-free stacking of multiple flasks also in combination with the standard TPP tissue culture flasks.  
 Barrier creates a growth area of 115 cm<sup>2</sup>.  
 Both flasks are available with Filter screw cap with a hydrophobic PTFE membrane with a pore size of 0.22 µm only. If a «VENT» screw cap is used, the build-up of inside pressure of more than 0.03 bar could press open the flask.  
 A controlled level base surface and regular use of feeding medium are essential to uniform tissue growth.  
 Sources of vibrations must be eliminated to avoid circular growth pattern.

**Technical Data**

**Materials**

Flask PS  
 Screw cap PE  
 Membrane PTFE,  
 pore size 0.22 µm  
 Gasket non-toxic gasket

<b>Abmessungen</b>	<b>Measurement</b>	<b>90552</b>	<b>90652</b>
Wiederverschliessbar	Re-closable	✓	✓
Steg	Barrier	---	✓
Länge mm	Length mm	210	210
Breite mm	Width mm	122	122
Höhe mm	Heights mm	50	50
Wachstumsfläche cm <sup>2</sup>	Growth area cm <sup>2</sup>	115	115
empf Volumen ml	Recom. volume ml	15 – 45	15 – 33
Öffnung mm	Opening mm	105 x 105	105 x 105
max Volumen ml	Max volume ml	100	100