



TWIRL'EM®

TWIRL'EM®

The TWIRL'EM® provides a safe, contaminant-free and flexible container that ensures reliable test results. Its integrated secure closure system makes it a sealed and airtight solution. Proven efficient and versatile, it is used by many professionals in various industries.

The simple and economical tool that will support you in all your sampling activities.

Features and benefits

- Made of FDA compliant material
- Can accommodate solid, semi-solid and liquid samples
- Maximum bag opening to facilitate insertion of sample
- Various sizes and closure combinations available depending on the type of sample
- Suitable for use in the homogenizer
- Guaranteed sterility
- RNase, DNase and pyrogen free




Inspired by your challenges

LABPLAS


VISIT LABPLAS.COM FOR ADDITIONAL INFORMATION ABOUT OUR PRODUCTS | Tel: 1.450.649.7343



TWIRL'EM® BAG INSTRUCTION

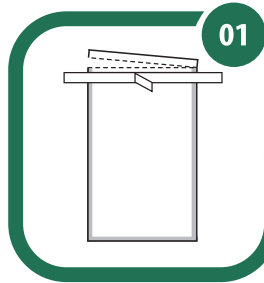

Tear off the top of the bag along the perforations

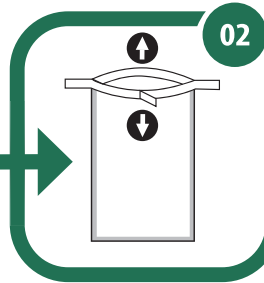

Use pull-tabs to open the bag




Aseptically insert the sample in the bag



Pull the wire-ends on each side of the bag



Twirl the bag around the wires, at least 3 times, to close the bag

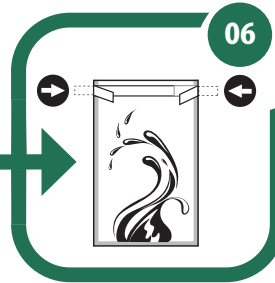

Fold both wire-ends onto the bag to provide airtight and leak-proof closure



Inspired by your challenges

LABPLAS

VISIT LABPLAS.COM FOR ADDITIONAL INFORMATION ABOUT OUR PRODUCTS | Tel: 1.450.649.7343

