R 104

<p>1. Inhalt/ Contents</p> <p>1x  1x  2x </p> <p>4x </p> <p>1x </p>	<p>2. Montage/ Assembly</p>  <p>Fig. 2.1</p>	 <p>Fig. 2.2</p>
 <p>Fig. 2.3</p>	 <p>Fig. 2.4</p>	 <p>Fig. 2.5</p>
 <p>max. 200 mm</p> <p>max. 15°</p> <p>max. 15°</p> <p>Belastung max. 0,7 kg Max. load 0,7 kg</p> <p>Drehmoment max. 0,5 Nm Max. torque 0,5 Nm</p> <p>Arbeitsbereich Working range</p> <p>370 mm</p> <p>210 mm 220 mm</p> <p>200 mm</p> <p>90 mm</p> <p>360 mm</p> <p>4x </p> <p>2x </p>		

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3. Operating and safety instructions

EN

In order to be able to use the stand **R 104** properly and safely, every user must first read the operating instructions and observe the safety instructions contained therein. Take care of these operating instructions and keep them in a place where they can be accessed by everyone.

Only staff who have been trained accordingly, know the stand **R 104**, and are authorised to carry out work in this field should use this stand.

Please follow the relevant safety instructions and guidelines, and occupational health and safety regulations for use in the laboratory. The user must select and wear his personal protective equipment according to the mixing hazard category.

Please unpack the stand **R 104** carefully and check for any damages. It is important that any damages which may have arisen during transport are ascertained when unpacking. If applicable a fact report must be set immediately (post, rail or forwarder).

Check that the contents are complete.

Assemble the stand **R 104** according to the instructions and ensure that all parts are firmly and securely connected. Assemble the stand **R 104** only on a stable, even, non-slip surface. You must also ensure that the stand **R 104** is set up in such a way that it is not liable to topple and does not start to move.

Stand **R 104** is only suitable for use in laboratories and for **IKA** laboratory instruments RW 11 basic, T 8, T 10 basic, DI 10 in conjunction with the following **IKA** accessories: cross sleeve H 44 or R 182, universal clamp R 350, clamping bracket RH 3.

Do not exceed the permitted loads, torques and distances. Always check that all connections are firm and secure before using the stand. Only use stand **R 104** with the original four rubber feet. Never use stand **R 104** if parts of the stand are damaged.

Technical Data:

Permitted ambient temperature:	°C	5 ... +40
Permitted relative humidity:	%	80
Weight:	kg	0.8

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IKA® - Werke GmbH & Co.KG
IKA® Works Inc.