Tube

ST-20/50 ...

DT-20/50 ...

BMT-20/50 ...

ST/DT/BMT - Tube aufstecken • Put on the ST/DT/BMT - Tube • Metre le ST/DT/BMT - Tube • Установка пробирки ST/DT/BMT

(RU)

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Safety instructions

- **Read the operating instructions in full before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the device.
- Follow the safety instructions, guidelines and safety and accident prevention regulations.
- **CAUTION!** For research purposes only! Not suitable for therapeutic or diagnostic use in human or veterinary medicine (in accordance with the MD Directive).
- Ensure that the cover is screwed on tightly.
- The tube must only be attached and removed while the motor is stationary.
- Ensure that the tube is firmly attached to the bayonet lock connector on the drive unit prior to operating the device.
- Ensure that the tube is able to withstand the chemical properties of the selected medium by filling the tube and storing for a suitable period of time.
- Always open the tube carefully after use as the media in the tube may heat up due to transfer of energy during operation, leading to pressurization of the container. Risk of material spraying. Protective equipment must be worn.
- Abrasion of the dispersion equipment or the rotating accessories can get into the medium you are working on.
- While working with the disperser, the user must wear his personal protective equipment, in particular safety goggles.
- The temperature of the material must not exceed 40 °C.
- Note the operating instructions of the disperser unit.

Product information

ST20/50..., comprise mixing vessel and stirring tool, DT20/50..., comprise mixing vessel and dispersion tool, BM17/20/50..., comprise mixing vessel, stirring tool and ball.
 ST20, ST 20-M, ST50, ST50-M, DT 20-M, DT50, DT50-M and BM17/20, BM17/20-M, BM17/50, BM17/50-M are not sterile and not free of DNA or RNA.
 ST20-gamma, ST20-M-gamma, DT20-gamma, DT20-M-gamma and BM17/20-gamma, BM17/20-M-gamma are sterilized.
 DT20-M and DT50-M are tube covers with a permeable membrane for needs or pipette points.
 The tubes are ideal for applications, with which no cross contamination is permissible.
 The tubes ST20, DT20 and BM17/20 are appropriate only for a single-use. Multiple-use can destroy the tube and the disperser unit.

Application instruction

Store the tubes in a dry place, in their packaging and at room temperature. Note the expiry date of the tubes. Protect the tubes against ultraviolet light.
Attention! Never run the DT20/50... dry.
 For ST 20..., DT20..., and BM17/20..., the fill level of the medium must be between the 2, respectively 5 ml and max. 15 ml marks. For ST 20-M..., DT20-M..., and BM17/20-M..., the fill level of the medium must be between the 15 ml and max. 50 ml marks.
 Do not turn the tube after filling the medium when the kit's tightness is ensured no longer.
 Always use the tube after filling the medium at room temperature. The tightness may be affected if the solid bodies in the medium are frozen or too hard. Dispensing causes the medium to heat up. Likewise the optimal dispersion duration and rotating frequency must be determined by attempts. Usually a few seconds are sufficient. maximum duration lines bring no improvements, increase however the sample temperature substantially.
 Longer application times may be necessary according to the regulations. Observe the laboratory regulations. Dispose of the tubes after use according to the regulations. Observe the laboratory regulations.

Materials and chemical resistance

The containers are manufactured from polypropylene (PP), polyetheretherketone (PEEK), rotor/stator are manufactured from polyetheretherketone (PEEK) and polyetheretherketone with teflon (PEEK+PTFE), at DT20/50 eco are rotor/stator of polyethersulfone (PES) and polyetheretherketone with teflon (PEEK+PTFE).
 The balls are manufactured from either stainless steel A19 304 or soda-lime glass.
 The used plastic materials conform with the FDA (Food and Drug Administration Approval) rules.
 The parts have a good stability against weak acids, chlorides, hypochlorites and many other chemicals.

Technical data

Designation	For disperser unit	Working range (ml)	Stator (Ø) (mm)	Rotor (Ø) (mm)	Circumferential speed at 6000 rpm (m/s)
ST20...	Tube Drive 2-15	2-15	-	-	-
DT20...	Tube Drive 5-15	5-15	26	20	6,5
BM17/20...	Tube Drive 2-15	2-15	-	-	-
ST50...	Tube Drive 15-50	15-50	-	-	-
DT50...	Tube Drive 15-50	15-50	26	20	6,5
BM17/50...	Tube Drive 15-50	15-50	-	-	-

Warranty

In accordance with IKA's warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs. The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

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